







BINDER Scientific  
Product catalog

# Table of contents

	03	BINDER	
	11	Growth	1
	13	Cell and tissue cultures – C   CB	
	23	Microbiology – BD   BF   KB	
	37	In vivo and In vitro cultures – KBW   KBWF	
	47	Storage	2
	49	Sample storage – UF V	
	55	Drying and tempering	3
	57	Drying and Heat treatment – ED   FD   FED	
	71	Material testing – FP   M	
	81	Safety drying – FDL   MDL	
	91	Vacuum drying – VD   VDL	
	103	Environmental simulation	4
	105	Constant climate testing – KBF	
	111	Constant climate testing with illumination – KBF P   KBF LQC	
	121	Dynamic constant climate testing – KMF	
	127	Process documentation APT-COM™	5
	131	Service and BINDER INDIVIDUAL	6
	140	Product selection chart	



# BINDER – The Mission

## | Technology for a better world

BINDER's core purpose is to make a significant technical contribution to improving the health and safety of mankind. Our aim is to provide our customers with products that best support their daily laboratory work and allow them to continuously improve results.

We strive unremittingly for perfection and to differentiate ourselves from others through continuous innovation. To this end, the percentage of our revenues that we invest in research and development annually is well above the industry average.

Consequently, our customers are at the center of all our reflections and activities.

They not only work with our products, they also inspire and drive us to improve every day.

Everyone at BINDER is dedicated to meeting our customers' requirement beyond their expectations.

To ensure the highest quality standards, all our products are solely manufactured at our state-of-the-art plant in Germany. Our exclusive focus on simulation chambers makes us the world's largest specialist in the market. We intend to maintain our step ahead and in that regard, we will work with the same responsibility, passion and desire as we have since the foundation of BINDER.

Peter Michael Binder

*President & CEO*





## BINDER – The company

### | Best conditions for your success

BINDER is a family business dedicated to simulation chambers. We are the world's largest specialist in simulation chambers for the scientific and industrial laboratory.

More than 23,000 units leave our plant in Tuttlingen annually.

Proven cutting-edge technologies, pioneering innovations and absolute precision characterize the BINDER brand image. Our focus lies in the perfect simulation of biological, chemical and physical environmental conditions for a number of industries. Our logo's three red triangles stand for: superior products, best service package and professional consultation. Our tagline embodies these values by providing: *"best conditions for your success"*.

In addition to our range of products, we offer our customers a comprehensive selection of options and accessories, making it possible to optimize standard units to your requirements profile. BINDER INDIVIDUAL offers flexible solutions adapted to the customer's particular requirements.

Whatever requirements may be, we can support and guide you locally from our headquarters in Tuttlingen and from our four BINDER offices in New York, Moscow, Hong Kong and Shanghai, as well as through our sales partners in more than 135 countries.





## BINDER – Environmental responsibility

### | Good for you. Good for the environment.

For nearly three decades, we have been offering more to our customers: more innovation, safety and local support. This is also the case for the environment. We take into account ecology and sustainable environmental protection not only in the development and production of our chambers, but in everything we do as a company.

With this in mind, BINDER is committed to use material combinations that can be separated to enable recycling. This is why asbestos or mercury are not used in any BINDER chamber today or will ever be in the future. We consciously choose synthetic materials made of recycled resin in our manufacturing process without any compromise in quality.

- ▶ **Less energy consumption** – The geothermal heating system at the Research & Development Center has a transmission heat requirement of 40 % below that required by the German Energy Saving Ordinance (EnEV). Daylight harvesting system: large windows allow optimum use of daylight supported by daylight-dependent automatic blinds; green roof for heat insulation; energy savings by the use of state-of-the-art equipment and systems and adaptive lighting systems. We are dedicated to reducing energy consumption which is why a combined heating and power station is in planning.
- ▶ **Stringent selection of suppliers in accordance with the European Ecodesign Directive** – Reusable packaging with upstream suppliers is used whenever possible. Materials such as steel, plastics, cardboard and paper are consistently recycled. Renewable and recycled materials are given preference wherever possible.
- ▶ **Less environmental impact** – The extensive oil-free steel processing performed with automatic machines allows BINDER to dispense with cleaning procedures detrimental to the environment before surface coating. Only solvent-free powder coating systems are used at BINDER.
- ▶ **Less resource consumption** – Use of up to 100 % recyclable packaging; 98 % powder reprocessing thanks to closed circuit coating system; 70 % less paper consumption thanks to digital workflow.

URKUNDE

an  
Deutsche Gesellschaft  
für Züchtungslehre

am 20.01.2011  
BINDER Innovations-Preis 2011

Herrn Univ.-Prof. Dr. rer. nat. Christof Bär

Er hat sich um die Entwicklung  
„Agly-Tissue Culture“  
im Rahmen der Projekt „Innovations-Preis 2011“ verdient gemacht.

Datum: 20.01.2011

BINDER





# BINDER – Research and progress

## The BINDER Innovation Prize: An excellent idea.

Since 1998, the BINDER Innovation Prize has been awarded annually by the German Society for Cell Biology (DGZ). The prize is endowed by BINDER. It is awarded to candidates for their outstanding work in the field of basic cell biology research.

With this award, BINDER commits to the continuous advancement of life sciences and deepens its close cooperation with members of the science and research community. Especially for projects involving cell culture, BINDER is an ideal partner for the realization of sophisticated endeavors.

- |  |  |
|--|--|
| <p><b>1998</b>   Dr. Ronald Frank, <i>Society for Biotechnological Research Braunschweig</i>, for his work in the field of SPOT analysis</p> <p><b>1999</b>   Dr. Ludger Fink, <i>Institute for Pathology at the Justus Liebig University of Giessen</i>, for his work on mRNA analysis according to laser-assisted cell-picking from complex tissues</p> <p><b>2000</b>   Dr. Maria Wartenberg and Dr. Heinrich Sauter, <i>Institute for Neurophysiology in Cologne</i>, for their work: "Thalidomide inhibits angiogenesis in embryoid bodies by the generation of hydroxyl radicals"</p> <p><b>2001</b>   Dr. Nicole Maas-Szabowski and Dr. Axel Szabowski, <i>German Cancer Research Center in Heidelberg</i>, for their in-vitro research on the complex interactions between the human subcutis and cutis</p> <p><b>2003</b>   Dr. Thomas Korff and Prof. Hellmut Augustin, <i>Cancer Biology Clinic in Freiburg</i>, for fundamental research on an in-vitro model of the endothelium differentiation and blood vessel formation (angiogenesis)</p> <p><b>2004</b>   Dr. M. Christina Cardoso, <i>Max-Delbrück Center for Molecular Medicine (MDC) Franz-Volhard-Klinik, Berlin</i>, for her work: "A look at genome duplication in living cells and the development of cell cycle markers"</p> | <p><b>2005</b>   Prof. Frauke Melchior, <i>Georg-August-University, Göttingen</i>, for her work in the field of regulating cellular processes through ubiquitin-related proteins of the SUMO family</p> <p><b>2006</b>   Prof. Dr. Ivan Dikic, <i>University Frankfurt/Main</i>, for his work: "Ubiquitin and Ub-like modifiers in the heart of cell signalling"</p> <p><b>2007</b>   Dr. Jan-Michael Peters, <i>Research Institute of Molecular Pathology, Vienna</i>, for "Regulation of sister chromatid cohesion in mammalian cells"</p> <p><b>2008</b>   Prof. Dr. Heinrich Leonhardt and Dr. Ulrich Rothbauer, <i>Ludwig Maximilians, University Munich BioCenter, Department of Biology</i> for their work: "Targeting and tracing antigens in living cells with fluorescent nanobodies"</p> <p><b>2009</b>   Prof. Dr. Ludger Hengst, <i>Medical University of Innsbruck</i>, for his work in protein research, with the topic: CDK-Inhibitors – connecting signal transduction and cell cycle control</p> <p><b>2010</b>   Prof. Anne Spang, <i>University of Basel</i> for her work: Intracellular transport in yeast</p> <p><b>2011</b>   Prof. Dr. Christian Behl, <i>Institute for Pathobiochemistry, Johannes Gutenberg University Mainz</i> for the study of the biochemistry of aged cells in conjunction with the course of neurodegenerative processes</p> |
|--|--|



# Growth

## | Always the right conditions

Precise adherence to the pre-selected values for temperature, humidity, CO<sub>2</sub>, O<sub>2</sub> and light as well as reliable reproducibility of growth conditions are the two most important requirements for simulation chambers. On the other hand, they should make documenting your work with living organisms as easy as possible.

BINDER is aware of these requirements and works continuously to provide you with the perfect CO<sub>2</sub> incubators, microbiological incubators and growth chambers. Our product range offers innovative technical solutions and comprehensive equipment options. We can also meet special technical requirements through BINDER INDIVIDUAL.







# Cell and tissue cultures

## C | CB series

### Absolutely superior: BINDER CO<sub>2</sub> incubators

- ▶ Absolute stability
  - ▶ Reproducibility of the best culture conditions
  - ▶ High temperature uniformity provided by the APT.line™ preheating chamber
  - ▶ Saturated relative humidity up to 95 % r. H.
  - ▶ Stable pH values thanks to drift-free CO<sub>2</sub> / O<sub>2</sub> IR sensor technology
- ▶ Absolute reliability
  - ▶ Reliable decontamination concept
  - ▶ Validated and recognized 180 °C hot air sterilization routine
  - ▶ Condensation-free interior due to Permadyr™
- ▶ Absolutely clean
  - ▶ Reduction of interior contaminable surface to a minimum
  - ▶ Easy-to-clean interior

The BINDER CO<sub>2</sub> incubators ensure optimal and reproducible growth conditions for your cultures. The proven APT.line™ preheating chamber and Permadyr™ eliminate the risk of condensation. You can depend on hot air sterilization at 180 °C, to eliminate bacteria and spores. The unique seamless, deep-drawn chamber without obstructions makes cleaning the chamber simple. The CO<sub>2</sub> incubators of the C and CB series offer you all these benefits standard for successful cell and tissue cultivation standard.



Cell culture



Tissue engineering



In vitro fertilization



# The standard for CO<sub>2</sub> incubators

## C series

The C 150 series ensures optimal growth conditions up to 50 °C. With a highly effective decontamination concept, it meets all basic requirements for cell and tissue cultures. A choice of options allows for an incubator configuration based on the application requirements.



► C 150 model



150

Available size (liters)

### ► EQUIPMENT


- Temperature range 7 °C above ambient temperature to 50 °C
- Hot-air sterilization at 180 °C
- Microprocessor with LED display for temperature and CO<sub>2</sub> concentration
- Various alarm and status displays
- Automatic diagnostic system with visual and audible alarm, as well as zero-voltage relay contact for central monitoring
- Drift-free infrared CO<sub>2</sub> measurement system
- Seamless, deep-drawn inner chamber made of stainless steel with integrated shelf support system
- Water pan
- Independent temperature safety device class 3.1 (DIN 12880) with visual and audible temperature alarm
- Gas mixing head
- Tightly-fitted inner glass door
- Door lock
- Door hinged right or left
- Units are stackable
- Three perforated shelves, stainless steel

## ► C SERIES | BEST GROWTH CONDITIONS:




### Uniform growth conditions

- APT.line™ preheating chamber
  - Homogeneous temperature distribution
  - Exact temperatures at all levels
- High humidity at 95 % r. H.
  - Minimizes the risk of evaporation

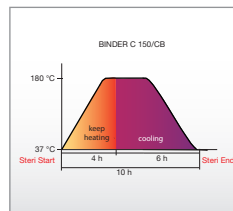

### Reliable decontamination

- 180 °C hot-air sterilization
  - Complete elimination of contaminants
  - Automatic sterilization cycle
  - Globally recognized method

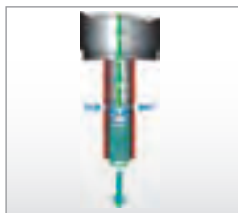

### Stable pH value

- Drift-free CO<sub>2</sub> sensor technology
  - Gas mixing head ensuring an homogeneous CO<sub>2</sub> distribution
  - Homogeneous conditions quickly attained



### Easy cleaning


- Seamless, deep-drawn inner chamber
  - High-quality stainless steel materials
  - Minimization of contaminable surface
  - Time-saving manual cleaning

## ► OPTIONS

- Access ports of various diameters with silicone plug
- Perforated rack, stainless steel
- Gas cylinder connection set
- Analog outputs 4 – 20 mA for temperature and CO<sub>2</sub> measurements
- BINDER Gas Supply Service (external gas tank changer)
- Stacking adapter
- Base on casters
- Calibration certificate for temperature and CO<sub>2</sub>
- Temperature measurement according to DIN 12880



Base on casters

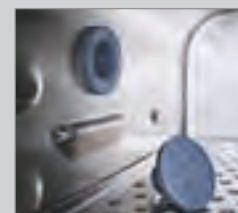

Silicone access port with two silicone plugs



Gas Supply Service



CO<sub>2</sub> incubator with integrated turning gear

► BINDER  
INDIVIDUAL  
More information  
on page 133

# Technical data for your planning and installation

## C series



	C 150
<b>Exterior dimensions</b>	
Width (mm/inch)	680 / 26.77
Height (including feet) (mm/inch)	820 / 32.28
Depth (mm/inch)	815 / 32.09
plus i-triangle, door handle, connections (mm/inch)	105 / 4.13
<b>Interior dimensions</b>	
Width (mm/inch)	500 / 19.69
Height (mm/inch)	600 / 23.62
Depth (mm/inch)	500 / 19.69
Interior volume (l/cu.ft.)	150 / 5.4
Perforated shelves (no. standard/max.)	3 / 6
Weight (empty) (kg/lbs.)	95 / 210
<b>Temperature data</b>	
Temperature range, 7 °C / 12.6 °F above ambient up to °C/°F	50
Temperature uniformity at 37 °C / 98.6 °F (±K)	0.4
Temperature fluctuation (±K)	0.1
Recovery time after door was opened for 30 sec <sup>1)</sup> at 37 °C / 98.6 °F (min)	5 / 122
<b>CO<sub>2</sub> data</b>	
CO <sub>2</sub> range (Vol.-% CO <sub>2</sub> )	0 – 20
Setting accuracy (Vol.-% CO <sub>2</sub> )	0.1
Recovery time after door was opened for 30 sec <sup>1)</sup> up to 5 vol. % (min)	7
CO <sub>2</sub> measurement	IR
<b>Humidity data</b>	
Humidity average value (% r. H.)	95 ±2
<b>Electrical data</b>	
IP system of housing protection acc. to EN 60529	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230
Nominal power (kW)	1.4
Energy consumption <sup>2)</sup> at 37 °C / 98.6 °F (Wh/h)	110
<b>Model no.</b>	
	9040-0078

<sup>1)</sup> to 98 % of the set value. The recovery times of the gas concentrations inside the chamber following door opening refer to a connection pressure of 2.0 bar. Decreasing supply pressure leads to longer recovery times. // <sup>2)</sup> Use this value for dimensioning air conditioning systems. /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.




Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



► DIMENSIONS

C 150

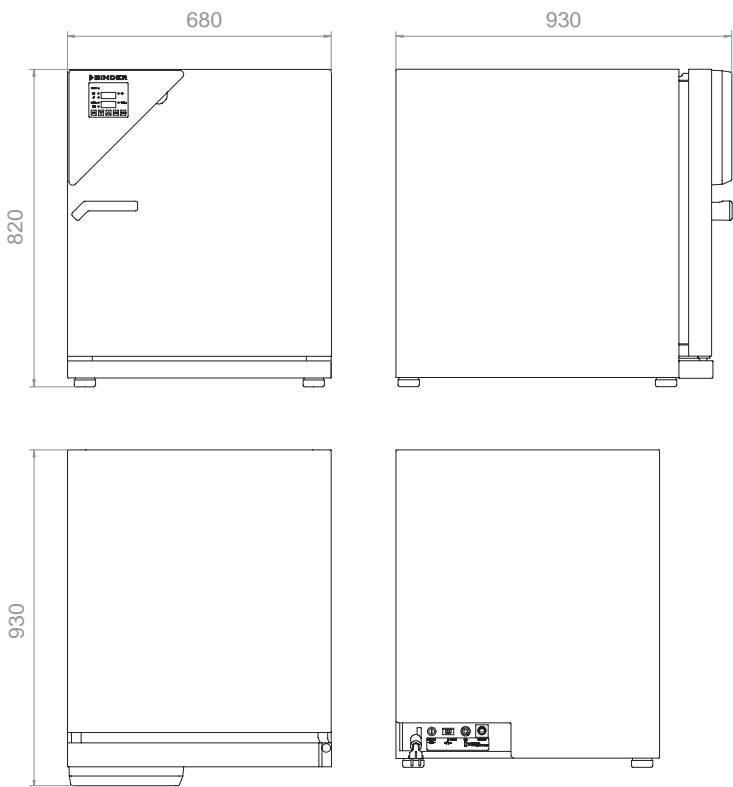

► INSTALLATION REQUIREMENTS

	C 150
Wall clearance rear (mm/inch)	100 / 3.94
Wall clearance side (mm/inch)	50 / 1.97
Connection hose nozzle DN6 for CO <sub>2</sub> for hosewith internal diameter (mm/inch)	6 / 0.24
Nominal voltage (±10 %) 50/60 Hz (V)	230
Nominal power (kW)	1.4

# CO<sub>2</sub> incubator with additional process control CB series | CB series with O<sub>2</sub> control

The CB series offers optimal growth conditions up to 60 °C combined with a highly effective 180 °C sterilization routine. A wide range of options such as O<sub>2</sub> control for hypoxic conditions, glass door or roller system are available for a variety of application-specific solutions. Our premium class incubator meets the highest demands on accuracy, safety and data communication.



► CB 150 model  
with divided inner glass door option



Available sizes (liters)

## ► EQUIPMENT

- Temperature range from 7 °C above ambient temperature to 60 °C
- MCS controller for temperature and CO<sub>2</sub> concentration
  - User-friendly LCD screen
  - Easy-to-read menu guidance
  - Integrated electronic chart recorder
  - Variety of options for the graphic display of process parameters
  - Real-time clock
- Hot air sterilization at 180 °C (DIN 58947)
- VENTAIR™ air jacket system
- Drift-free infrared CO<sub>2</sub> measurement system
- Patented gas mixing nozzle
- Permadyr™ System; condensation-free double-pan humidification system, maintains dry interior walls
- Seamless deep-drawn inner chamber made of stainless steel with integrated shelf support system
- Electronic error self-diagnostic system with visual and audible alarms, as well as potential free contact for central monitoring
- Independent temperature safety device class 3.1 (DIN 12880) with visual and audible temperature alarm
- Tightly-fitted inner glass door
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Three perforated shelves, stainless steel
- Door lock



## ► CB SERIES | BEST GROWTH CONDITIONS:




### Uniform growth conditions

- VENTAIR™ air jacket system
  - Homogeneous temperature distribution at all levels
  - Extremely fast recovery times



### Easy cleaning

- Seamless, deep-drawn inner chamber
  - High-quality stainless steel materials (1.4301)
  - No permanent fixtures
  - Removable water pan
  - Integrated shelf support system

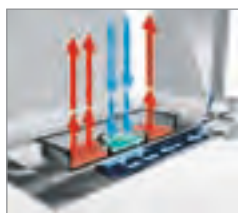


### High humidity

- Permady™ double-pan humidification system
  - Saturated relative humidity up to 95 % r. H.
  - Rapid humidity recovery times
  - Optimal evaporation protection for cultures




### Stable pH values

- Infrared measurement system
  - Drift-free CO<sub>2</sub>/O<sub>2</sub> sensor technology
  - Rapid response time
  - Precise measurement results

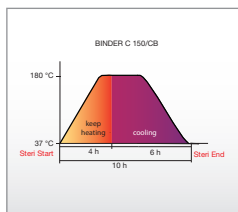

### Reliable decontamination

- 180 °C hot-air sterilization
  - Complete elimination of contaminants
  - Automatic sterilization cycle
  - Globally recognized method

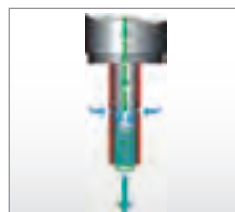

- Gas mixing head with Venturi effect
  - Homogeneous CO<sub>2</sub>/O<sub>2</sub> distribution
  - Fast effective aeration

## ► OPTIONS

- Silicone access port with two silicone plugs
- Perforated shelf made of stainless steel
- Divided shelf made of stainless steel for gas-tight divided inner glass door
- Lockable controller keyboard
- Gas tank connection kit CO<sub>2</sub>, O<sub>2</sub>, N<sub>2</sub>
- BINDER Gas Supply Service
- Unit conversion from RS 422 to Ethernet interface
- 4 – 20 mA analog outputs for temperature and CO<sub>2</sub> measurements
- Switched LEMO inner chamber socket
- Independent Intelligent Fail Safe monitoring function
- Calibration certificate for temperature and CO<sub>2</sub> or O<sub>2</sub>
- Temperature measurement according to DIN 12880
- Current feedthrough (8-pin) for low voltage with LEMO socket
- Internal CO<sub>2</sub>, O<sub>2</sub> or N<sub>2</sub> tank changer for connecting two gas tanks
- CELLROLL Set
- Stacking adapter:
  - Base on casters
  - Stable vibration-free stacking frame for two CO<sub>2</sub> incubators



Stacking adapter for thermal decoupled stacking of two CO<sub>2</sub> incubators (C and CB)



Gas Supply Service



CO<sub>2</sub> incubator with 9-compartment glass door

► BINDER  
INDIVIDUAL  
More information  
on page 133

# Technical data for your planning and installation

## CB series | CB series with O<sub>2</sub> control



	CB 53	CB 150	CB 210
<b>Exterior dimensions</b>			
Width (mm/inch)	580 / 22.83	680 / 26.77	740 / 29.13
Height (including feet) (mm/inch)	720 / 28.35	920 / 36.22	1070 / 42.13
Depth (mm/inch)	550 / 21.65	715 / 28.15	715 / 28.15
Depth plus instrument triangle, door handle (mm/inch)	75 / 2.95	75 / 2.95	75 / 2.95
<b>Interior dimensions</b>			
Width (mm/inch)	400 / 15.75	500 / 19.69	560 / 22.05
Height (mm/inch)	400 / 15.75	600 / 23.62	750 / 29.53
Depth (mm/inch)	330 / 12.99	500 / 19.69	500 / 19.69
Interior volume (l/cu.ft.)	53 / 1.9	150 / 5.4	210 / 7.5
Perforated shelves (no. standard/max.)	2 / 3	3 / 6	3 / 8
Load per shelf (kg/lbs.)	10 / 22	10 / 22	10 / 22
Permitted total load (kg/lbs.)	30 / 66	30 / 66	30 / 66
Weight (empty) (kg/lbs.)	60 / 132	107 / 236	121 / 267
<b>Temperature data</b>			
Temperature range, above ambient (°C/°F)	5 – 60 / 9 – 140	7 – 60 / 12.6 – 140	7 – 60 / 12.6 – 140
Temperature uniformity at 37 °C / 98.6 °F (±K)	0.3	0.3	0.4
Temperature fluctuation at 37 °C / 98.6 °F (±K)	0.1	0.1	0.1
Recovery time after door was opened for 30 sec <sup>1)</sup> at 37 °C / 98.6 °F (min)	3	3	3
<b>CO<sub>2</sub> data</b>			
CO <sub>2</sub> range (Vol.-% CO <sub>2</sub> )	0 – 20	0 – 20	0 – 20
Setting accuracy (Vol.-% CO <sub>2</sub> )	0.1	0.1	0.1
Recovery time after door was opened for 30 sec <sup>1)</sup> at 5 vol. % (min)	6	7	9
<b>Humidity data</b>			
Humidity average value (% r. H.)	95 ±2	95 ±2	95 ±2
<b>Model no. standard equipment</b>	<b>9040-0069</b>	<b>9040-0038</b>	<b>9040-0039</b>
<b>O<sub>2</sub> data</b>			
O <sub>2</sub> range (vol.-% O <sub>2</sub> )	0.2 – 95	0.2 – 95	0.2 – 95
Setting accuracy (vol.-% O <sub>2</sub> )	0.1	0.1	0.1
Recovery time <sup>1)</sup> after door was opened for 30 sec at 1 vol.-% O <sub>2</sub> (min)	25	33	54
Recovery time <sup>1)</sup> after door was opened for 30 sec at 5 vol.-% O <sub>2</sub> (min)	9	10	11
<b>Model no. with O<sub>2</sub> control</b>	<b>9040-0071</b>	<b>9040-0044</b>	<b>9040-0045</b>


<sup>1)</sup> to 98 % of the set value. The recovery times of the gas concentrations inside the chamber following door opening refer to a connection pressure of 2.0 bar. Decreasing supply pressure leads to longer recovery times. // <sup>2)</sup> Use this value for dimensioning air conditioning systems. /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.




Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)

## ► DIMENSIONS


### CB 53

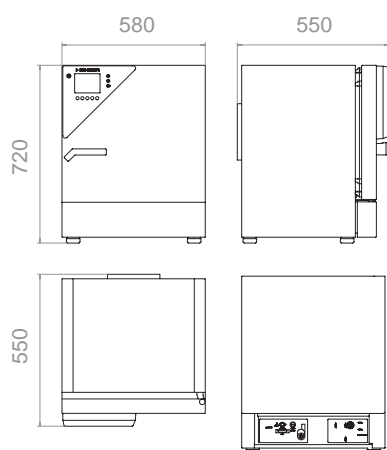

### CB 150

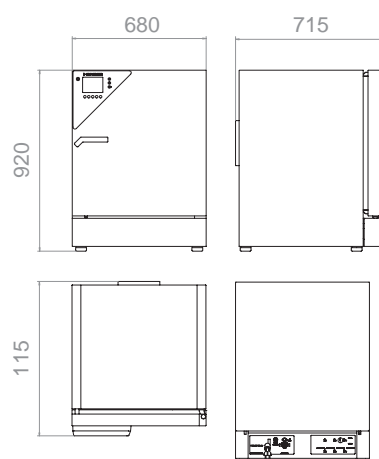

### CB 210

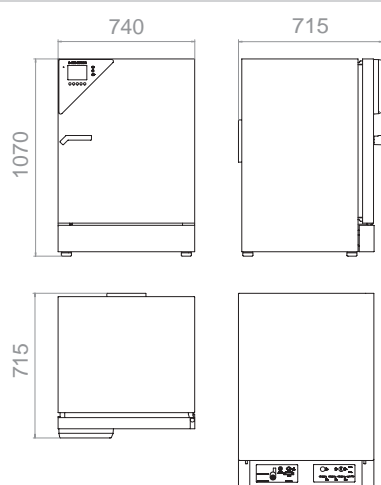

## ► INSTALLATION REQUIREMENTS

	CB 53	CB 150	CB 210
Connection hose nozzle DN6 for CO <sub>2</sub> for hose with internal diameter (mm/inch)	6 / 0.24	6 / 0.24	6 / 0.24
Connection hose nozzle DN6 for O <sub>2</sub> /N <sub>2</sub> for hose with internal diameter (mm/inch)	6 / 0.24	6 / 0.24	6 / 0.24
IP system of housing protection acc. to EN 60529	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230	230	230
Nominal power (kW)	1	1.3	1.5
Energy consumption <sup>2)</sup> at 37 °C / 98.6 °F (Wh/h)	70	120	130
Wall clearance side (mm/inch)	50 / 1.97	50 / 1.97	50 / 1.97
Wall clearance rear (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94



# Microbiology

## BD | BF | KB series

### Outstanding in detail: BINDER Microbiological incubators

- ▶ Outstanding precision
  - ▶ Reliably best culture conditions
  - ▶ High temperature uniformity thanks to APT.line™ preheating chamber
  - ▶ Tenths of a degree accuracy, precise temperature setting
  - ▶ Adjustable fan speed
- ▶ Outstandingly simple
  - ▶ Reliable decontamination concept
  - ▶ Disinfection routine at 100 °C
  - ▶ Inner chamber made of stainless steel with autoclavable fixtures
  - ▶ Inner glass door: “First look, then open”

When working with microorganisms, nothing is more crucial than precision incubation which remains stable in the long term. With the APT.line™ preheating chamber and the cooling system, we have developed two components that stand for unique precision – and results that are reproducible in each routine test. Whether for food microbiology or hygiene control, BINDER is your supplier of choice in the area of microbiological incubators. We offer a range of incubators that meet any application need.



Qualitative proof



Quantitative proof



In vitro tests



# Incubators with gravity convection

## BD series

BD series incubators from BINDER with gravity convection are specially designed for long-term and stable continuous operation, ideal for gentle incubation of organisms, as well as conditioning of heat-sensitive media.




► BD 23 model



Available sizes (liters)

### ► EQUIPMENT

- Temperature range 5 °C above ambient temperature to 100 °C
- DS control with integrated timer 0 to 99 hrs
- Digital temperature setting with an accuracy of a tenth of a degree
- Independent temperature safety device class 3.1 (DIN 12880) with visual alarm
- Adjustable ventilation flap slide and rear exhaust Ø 50 mm
- Inner glass door
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Units up to 115 liters are stackable
- Two chrome-plated racks included

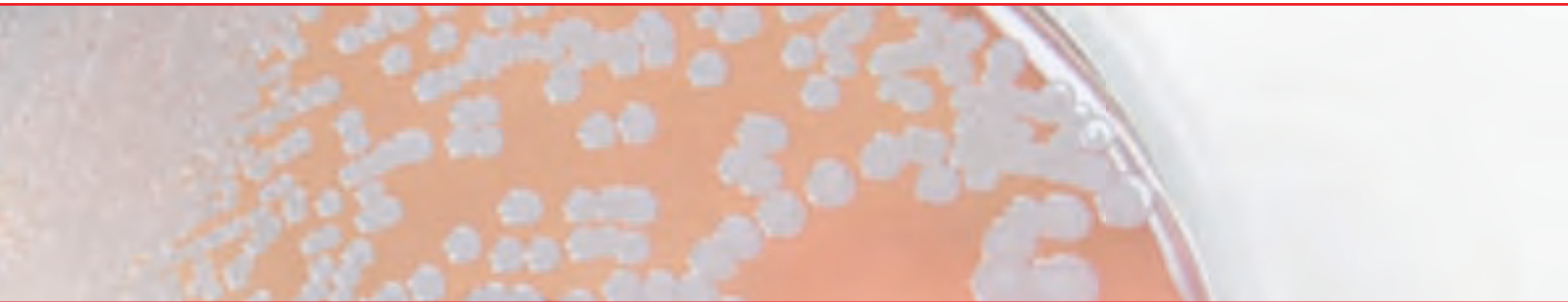

## ► BD SERIES | BEST CULTURE CONDITIONS:



### Reproducible growth conditions

- BINDER APT.line™ preheating chamber
  - Gentle temperature distribution with gravity convection
  - Homogeneous temperature at all levels



### Minimal operating costs

- Energy saving
  - Hermetic door with 2-points door closure
  - Low heat dissipation due to 60 mm insulation
- Stable atmosphere
  - Inner glass door



### Reliable decontamination

- Disinfection routine at 100 °C
  - Reduced risk of cross contamination
- Simple and fast cleaning
  - High-quality stainless steel inner chamber
  - No permanent fixtures

## ► OPTIONS

- Access ports of various diameters with silicone plug
- Racks, chrome-plated or stainless steel
- Perforated shelf
- Door lock
- Switchable waterproof interior socket 230 V AC
- Switchable audio alarm for overheating
- Additional PT 100 temperature sensor
- Temperature measurement according to DIN 12880
- Calibration certificate
- Extension to factory calibration certificate (additional measuring point)
- Data Logger Kits and Logger Software



Data Logger Kits



Additional PT 100 temperature sensor



Access ports with silicone plug



BD 53 incubator with left-hinged door

► **BINDER  
INDIVIDUAL**  
More information  
on page 133

# Technical data for your planning and installation

## BD series




	BD 23	BD 53	BD 115	BD 240	BD 400	BD 720
<b>Exterior dimensions</b>						
Width (mm/inch)	435 / 17.13	635 / 25.0	835 / 32.87	1035 / 40.75	1235 / 48.62	1235 / 48.62
Height (including feet/casters) (mm/inch)	495 / 19.49	620 / 24.41	705 / 27.76	825 / 32.48	1025 / 40.35	1530 / 60.24
Depth (mm/inch)	520 / 20.47	575 / 22.64	645 / 25.39	745 / 29.33	765 / 30.12	865 / 34.06
plus door handle, I-panel and exhaust duct (mm/inch)	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13
Quantity of doors	1	1	1	2	2	2
<b>Interior dimensions</b>						
Width (mm/inch)	222 / 8.74	400 / 15.75	600 / 23.62	800 / 31.50	1000 / 39.37	1000 / 39.37
Height (mm/inch)	330 / 12.99	400 / 15.75	480 / 18.90	600 / 23.62	800 / 31.50	1200 / 47.24
Depth (mm/inch)	277 / 10.91	330 / 12.99	400 / 15.75	500 / 19.69	500 / 19.69	600 / 23.62
Interior volume (l/cu.ft.)	20 / 0.7	53 / 1.9	115 / 4.1	240 / 8.6	400 / 14.3	720 / 25.7
Quantity of shelves (standard/max.)	2 / 3	2 / 4	2 / 5	2 / 7	2 / 10	2 / 15
Load per shelf (kg/lbs.)	12 / 26	15 / 33	20 / 44	30 / 66	35 / 77	45 / 99
Permitted total load (kg/lbs.)	25 / 55	40 / 88	50 / 110	70 / 155	90 / 199	120 / 265
Weight (empty) (kg/lbs.)	26 / 57	43 / 95	61 / 135	93 / 205	135 / 298	191 / 422
<b>Temperature data</b>						
Temperature range, 5 °C / 9 °F above ambient up to °C/°F	100 / 211.8	100 / 211.8	100 / 211.8	100 / 211.8	100 / 211.8	100 / 211.8
Temperature uniformity at 37 °C / 98.6 °F (±K)	0.5	0.5	0.4	0.5	0.5	0.5
Temperature fluctuation at 37 °C / 98.6 °F (±K)	0.2	0.1	0.1	0.1	0.1	0.1
Heating up time <sup>1)</sup> to 37 °C / 98.6 °F (min)	49	38	62	70	105	84
Recovery time after door was opened 30 sec <sup>1)</sup> at 37 °C / 98.6 °F (min)	3	5	5	5	6	4
<b>Electrical data</b>						
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230	230	230	230	230	230
Nominal power (kW)	0.2	0.4	0.4	0.68	0.85	1.25
Energy consumption at 37 °C / 98.6 °F (Wh/h)	11	11	20	33	56	80
<b>Model no.</b>						
	9010-0187	9010-0081	9010-0088	9010-0095	9010-0073	9010-0074

<sup>1)</sup> to 98 % of the set value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)

## ► DIMENSIONS

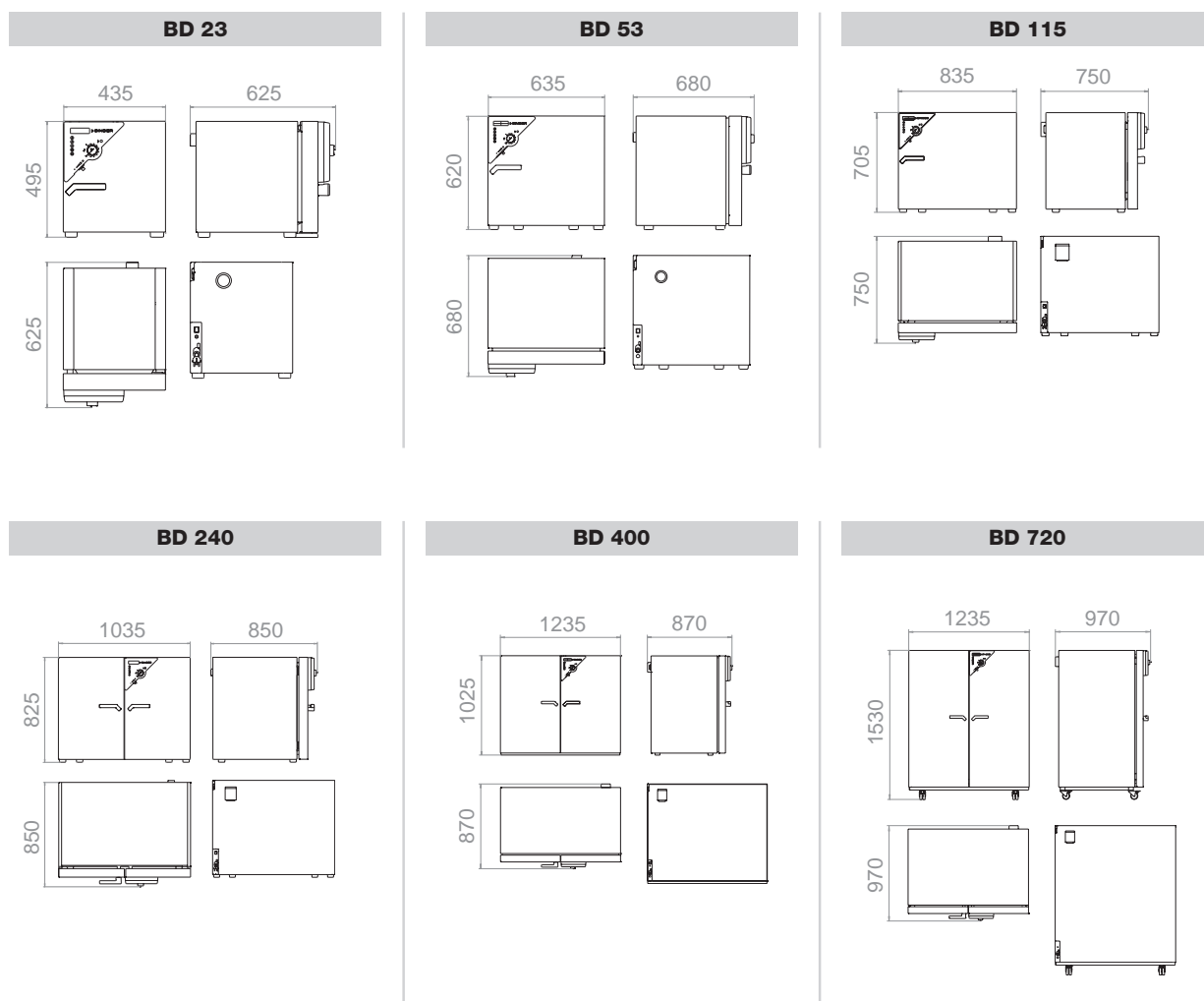

## ► INSTALLATION REQUIREMENTS

	BD 23	BD 53	BD 115	BD 240	BD 400	BD 720
Wall clearance rear (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94	100 / 3.94	100 / 3.94	100 / 3.94
Wall clearance side (mm/inch)	100 / 3.94	160 / 6.30	160 / 6.30	160 / 6.30	160 / 6.30	160 / 6.30
Nominal voltage (±10 %) 50/60 Hz (V)	230	230	230	230	230	230
Nominal power (kW)	0.2	0.4	0.4	0.68	0.85	1.25


# Incubators with mechanical convection

## BF series

The BF series is suitable for demanding incubation tasks, particularly for full loads and large batch throughputs. The incubator with mechanical convection provides optimal homogeneous conditions and has quick recovery times.



► BF 240 model

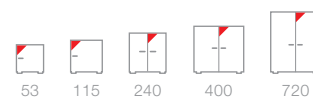

Available sizes (liters)

### ► EQUIPMENT

- Temperature range 5 °C above ambient temperature to 100 °C
- MS controller with several timer functions
- Timer functions: delayed ON, delayed OFF, temperature dependent delayed OFF
- Digital temperature setting with an accuracy of a tenth of a degree
- Adjustable fan speed
- Adjustable ventilation flap slide and rear exhaust Ø 50 mm
- Inner glass door
- Independent temperature safety device class 3.1 (DIN 12880) with visual alarm
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Adjustable intervals for printer
- Units up to 115 liters are stackable
- Two chrome-plated racks

## ► BF SERIES | BEST CULTURE CONDITIONS:



### Reproducible growth conditions

- BINDER APT.line™ preheating chamber
  - Homogeneous temperature conditions even under full load
  - Effective and gentle incubation
  - Rapid heating up and recovery times



### Reliable decontamination

- Disinfection routine at 100 °C
  - Reduced risk of cross contamination
- Simple and fast cleaning
  - High-quality stainless steel inner chamber
  - No permanent fixtures



### Minimal operating costs

- Energy saving
  - Hermetic door with 2-points door closure
  - Low heat dissipation due to 60 mm insulation
- Stable atmosphere
  - Inner glass door



### Standard-compliant documentation

- FDA-compliant APT-COM™ software
- Years of proven and recognized validation and documentation materials
- Standard RS 422 interface for network connection

## ► OPTIONS

- Access ports of various diameters with silicone plug
- Racks, chrome-plated or stainless steel
- Perforated shelf
- Door lock
- Switchable waterproof interior socket 230 V AC
- Switchable audio alarm for overheating
- Additional PT 100 temperature sensor
- Temperature measurement according to DIN 12880
- Calibration certificate
- Extension to calibration certificate
- Data Logger Kits and Logger Software



Data Logger Kits



Additional PT 100 temperature sensor



Calibration certificate



Incubator with keyboard lock

► **BINDER  
INDIVIDUAL**  
More information  
on page 133

# Technical data for your planning and installation

## BF series



	BF 53	BF 115	BF 240	BF 400	BF 720
<b>Exterior dimensions</b>					
Width (mm/inch)	635 / 25.0	835 / 32.87	1035 / 40.75	1235 / 48.62	1235 / 48.62
Height (including feet/castors) (mm/inch)	620 / 24.41	705 / 27.76	825 / 32.48	1025 / 40.35	1530 / 60.24
Depth (mm/inch)	575 / 22.64	645 / 25.39	745 / 29.33	765 / 30.12	865 / 34.06
plus door handle and exhaust duct (mm/inch)	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13
Quantity of doors	1	1	2	2	2
<b>Interior dimensions</b>					
Width (mm/inch)	400 / 15.75	600 / 23.62	800 / 31.50	1000 / 39.37	1000 / 39.37
Height (mm/inch)	400 / 15.75	480 / 18.90	600 / 23.62	800 / 31.50	1200 / 47.24
Depth (mm/inch)	330 / 12.99	400 / 15.75	500 / 19.69	500 / 19.69	600 / 23.62
Interior volume (l/cu.ft.)	53 / 1.9	115 / 4.1	240 / 8.6	400 / 14.3	720 / 25.7
Quantity of shelves (standard/max.)	2 / 5	2 / 5	2 / 7	2 / 10	2 / 15
Load per shelf (kg/lbs.)	15 / 33	20 / 44	30 / 66	35 / 77	45 / 99
Permitted total load (kg/lbs.)	40 / 88	50 / 110	70 / 155	90 / 199	120 / 265
Weight of the unit (empty) (kg/lbs.)	43 / 95	64 / 141	104 / 230	145 / 320	180 / 397
<b>Temperature data</b>					
Temperature range, 5 °C / 9 °F above ambient up to °C/°F	100 / 212	100 / 212	100 / 212	100 / 212	100 / 212
Temperature uniformity <sup>1)</sup> at 37 °C / 98.6 °F (±K)	0.4	0.3	0.3	0.4	0.4
Temperature fluctuation at 37 °C / 98.6 °F (±K)	0.2	0.2	0.2	0.2	0.1
Heating-up time <sup>2)</sup> to 37 °C / 98.6 °F (min)	12	22	12	18	21
Recov. time after door was opened for 30 sec. <sup>2)</sup> at 37 °C / 98.6 °F (min)	1	1	1	2	1
Air change at 70 °C / 158 °F (x/h)	59	29	19	17	11
<b>Electrical data</b>					
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230	230	230	230	230
Nominal power (kW)	0.4	0.4	0.68	0.85	1.25
Energy consumption at 37 °C / 98.6 °F (Wh/h)	11	20	33	56	80
<b>Model no.</b>					
	9010-0235	9010-0237	9010-0239	9010-0241	9010-0243

<sup>1)</sup> value without window // <sup>2)</sup> to 98 % of the set value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %.






The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



## ► DIMENSIONS

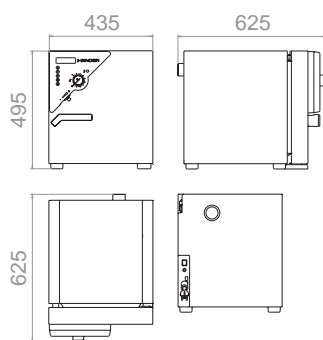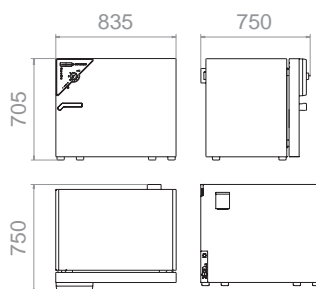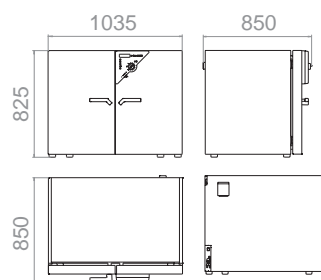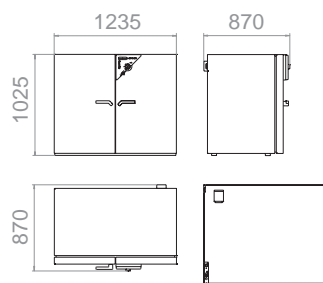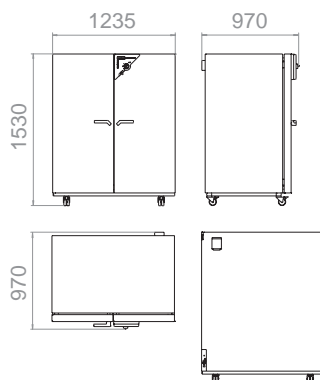**BF 53****BF 115****BF 240****BF 400****BF 720**

## ► INSTALLATION REQUIREMENTS

	BF 53	BF 115	BF 240	BF 400	BF 720
Wall clearance rear (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94	100 / 3.94	100 / 3.94
Wall clearance side (mm/inch)	160 / 6.30	160 / 6.30	160 / 6.30	160 / 6.30	160 / 6.30
Nominal voltage ( $\pm 10\%$ ) 50/60 Hz (V)	230	230	230	230	230
Nominal power (kW)	0.4	0.4	0.68	0.85	1.25


# Refrigerated incubator with mechanical convection KB series



The BINDER refrigerated incubators are all-rounders for the incubation of microorganisms. The KB series controls a temperature range of -5 °C to 100 °C. With their comprehensive program functions, they offer a wide range of capabilities and deliver reproducible test results.




► KB 240 model

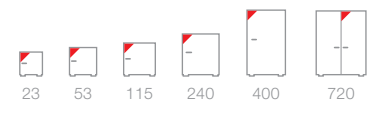

Available sizes (liters)

## ► EQUIPMENT

- Temperature range -5 °C to 100 °C
- MP controller with two programs with 10 sections each, alternatively switchable to one program with 20 sections
- Integrated weekly program timer with real-time function
- Adjustable ramp function via program editor
- Digital temperature setting with an accuracy of a tenth of a degree
- Adjustable fan speed
- Elapsed time indicator
- Independent temperature safety device class 3.1 (DIN 12880) with visual and audible alarm
- Inner glass door
- RS 422 interface for communication software APT-COM™ DataControlSystem
- Adjustable intervals for printer
- Units up to 115 liters are stackable
- Two shelves, stainless steel

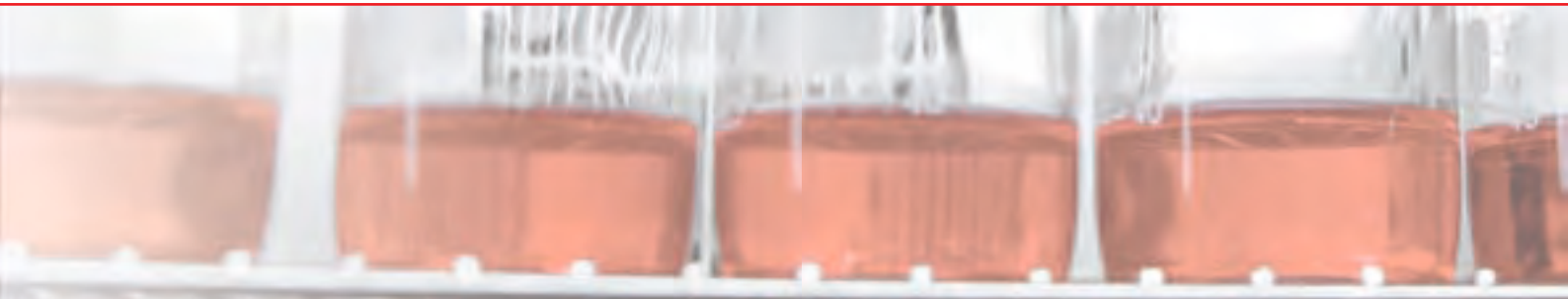

## ► KB SERIES | BEST CULTURE CONDITIONS:



### Wide range of programming options

- Multifunctional MP controller
  - 2 programs with 10 sections each
  - Weekly program function
  - Tenths of a degree accuracy, precise temperature setting
  - Standard RS 422 interface



### Reliable decontamination

- Disinfection routine at 100 °C
  - Reduced risk of cross contamination
- Simple and fast cleaning
  - High-quality stainless steel inner chamber
  - No permanent fixtures



### Reproducible incubation conditions

- APT.line™ preheating chamber
  - Homogeneous temperature conditions
  - Effective and gentle incubation
  - Rapid heating up and recovery times thanks to mechanical convection
- DCT™ cooling system
  - Minimum risk of evaporation
  - Incubation at 37 °C with integrated heat sources
  - Incubation even at high environmental conditions



### Minimal operating costs

- Energy saving
  - Hermetic door with 2-points door latch
  - Low heat dissipation due to 60 mm insulation
- Stable atmosphere
  - Inner glass door

## ► OPTIONS


- Access ports of various diameters with silicone plug
- Racks, chrome-plated or stainless steel
- Perforated shelf
- Reinforced shelf with additional mounting
- Door lock
- Waterproof interior socket 230 V AC
- Switchable audio alarm for over temperature
- Additional PT 100 temperature sensor
- Analog output for temperature 4 – 20 mA with 6-pin DIN socket
- Zero-voltage relay outputs accessible via 6-pin DIN socket
- Temperature measurement according to DIN 12880
- Independent temperature safety device class 3.3 (DIN 12880)
- Calibration certificate and extension to calibration certificate
- Data Logger Kits and Logger Software



Data Logger Kits



Watertight inner chamber socket

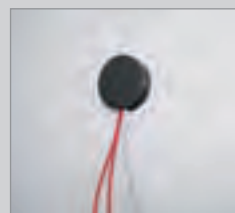

Various access ports



Refrigerated incubator with full-view glass doors

► **BINDER  
INDIVIDUAL**  
More information  
on page 133

# Technical data for your planning and installation

## KB series




	KB 23	KB 53	KB 115	KB 240	KB 400	KB 720
<b>Exterior dimensions</b>						
Width (mm/inch)	435 / 17.13	635 / 25.0	835 / 32.87	930 / 36.61	930 / 36.61	1255 / 49.41
Height (incl. feet/casters) (mm/inch)	620 / 24.41	840 / 33.07	1025 / 40.35	1460 / 57.48	1945 / 76.57	1925 / 75.79
Depth including door handle, I-panel, connection, cable (mm/inch)	600 / 23.62	665 / 26.18	730 / 28.74	800 / 31.50	800 / 31.50	890 / 35.04
Quantity of doors	1	1	1	2	2	2
<b>Interior dimensions</b>						
Width (mm/inch)	222 / 8.74	400 / 15.75	600 / 23.62	650 / 25.59	650 / 25.59	970 / 38.19
Height (mm/inch)	330 / 12.99	400 / 15.75	480 / 18.90	785 / 30.91	1270 / 50.0	1250 / 49.21
Depth (mm/inch)	277 / 10.91	330 / 12.99	400 / 15.75	485 / 19.09	485 / 19.09	576 / 22.68
Interior volume (l/cu.ft.)	20 / 0.71	53 / 1.87	115 / 4.06	247 / 8.73	400 / 14.13	698 / 24.66
Quantity of racks (standard/max.)	2 / 3	2 / 4	2 / 5	2 / 9	2 / 15	2 / 15
Load per rack (kg/lbs.)	10 / 22	15 / 33	20 / 44	30 / 66	30 / 66	45 / 99
Permitted total load (kg/lbs.)	25 / 55	40 / 88	50 / 110	100 / 221	120 / 265	150 / 331
Weight (empty) (kg/lbs.)	44 / 97	72 / 159	105 / 232	170 / 375	220 / 486	309 / 682
<b>Temperature data</b>						
Temperature range (°C/°F) <sup>1)</sup>	0 – 100 / 32 – 212	-5 – 100 / 23 – 212	-5 – 100 / 23 – 212	-5 – 100 / 23 – 212	-5 – 100 / 23 – 212	-5 – 100 / 23 – 212
Temperature uniformity at 25 °C / 77 °F (±K)	0.3	0.2	0.1	0.2	0.2	0.2
Temperature fluctuation max. (±K)	0.2	0.1	0.1	0.1	0.1	0.1
Recovery time <sup>2)</sup> after door was open for 30 sec at 5 °C / 41 °F (min)	5	4	5	16	14	12
<b>Electrical data</b>						
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230 1N~	230 1N~	230 1N~	200 – 240 1N~	200 – 240 1N~	200 – 240 1N~
Nominal power (kW)	0.34	0.46	0.46	1.20	1.4	2.3
Energy consumption <sup>3)</sup> at 40 °C (Wh/h)	60	215	115	260	420	510
<b>Model no.</b>						
	9020-0112	9020-0114	9020-0116	9020-0162	9020-0178	9020-0111

<sup>1)</sup> Lower values are valid up to an ambient temperature of max. 25 °C / 77 °F // <sup>2)</sup> to 98 % of the set value // <sup>3)</sup> These energy consumption values can be used upon calculation of air conditioning systems. The useable interior height depends on the position of racks. /// All technical data is specified for units with standard equipment at an ambient temperature of 25 °C / 77 °F and a mains voltage fluctuation of ±10 %. The temperature data is determined in accordance with factory standard following DIN 12880, observing the recommended wall clearances of 10 % of the height, width, and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



► DIMENSIONS

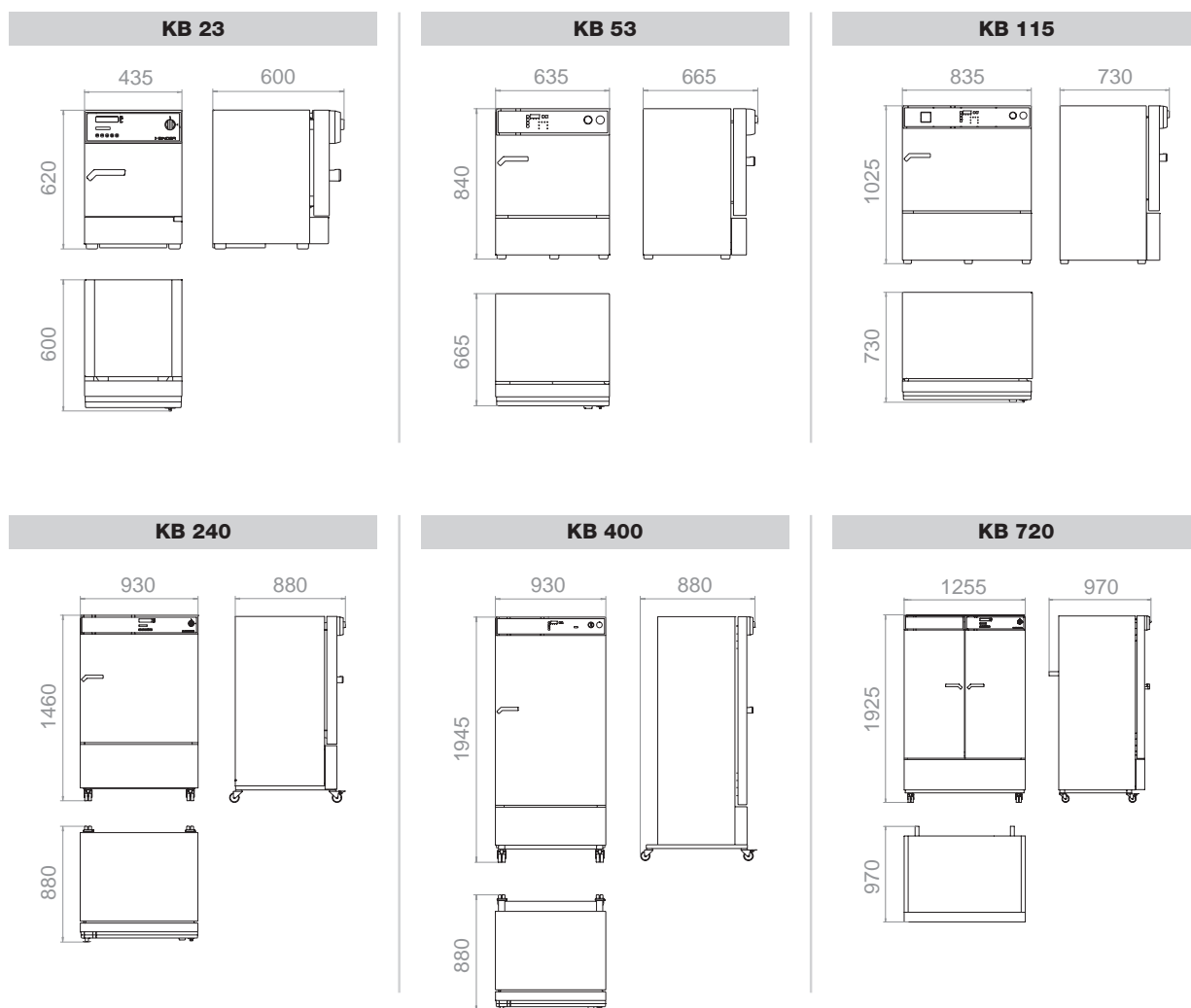

► INSTALLATION REQUIREMENTS

	KB 23	KB 53	KB 115	KB 240	KB 400	KB 720
Wall clearance rear (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94	100 / 3.94	100 / 3.94	100 / 3.94
Wall clearance side (mm/inch)	100 / 3.94	160 / 6.30	160 / 6.30	100 / 3.94	100 / 3.94	100 / 3.94
Nominal voltage ( $\pm 10\%$ ) 50/60 Hz (V)	230 1N~	230 1N~	230 1N~	200 – 240 1N~	200 – 240 1N~	200 – 240 1N~
Nominal power (kW)	0.34	0.46	0.46	1.2	1.4	2.3





# In vivo and In vitro cultures

## KBW | KBWF series

### Perfect conditions: BINDER Growth chambers

- ▶ Perfect climate
  - ▶ Responsive vapor pressure humidification ensure extremely low climatic gradients
  - ▶ Gentle air circulation provided by double-sided horizontal air flow
- ▶ Perfect light
  - ▶ Patented lighting system ensures homogeneous light distribution
  - ▶ Height-adjustable lighting system, three lamp types to choose from
- ▶ Perfect handling
  - ▶ Reliable decontamination concept
  - ▶ Easy-to-clean interior volume made of stainless steel 1.4301
  - ▶ Integrated weekly program in real-time

Homogeneous lighting conditions combined to constant temperature and humidity conditions are the outstanding features of the BINDER growth chambers. The APT.line™ preheating chamber and the responsive humidification system ensure powerful cooling and uniform light distribution for the best culture conditions. The KBW and KBWF series complements our comprehensive product portfolio and is a reliable partner for diverse applications.



In vitro cultures



In vivo cultures



Quality tests


# Growth chambers with illumination

## KBW series

The BINDER KBW series growth chambers achieve homogeneous light distribution and ensure consistent temperature conditions with its natural illumination. Minimized response times guarantee that all test and growth parameters remain balanced. The weekly program timer integrated in the controller is a real plus in handling.



► KBW 240 model

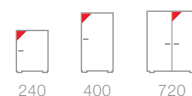

Available sizes (liters)

### ► EQUIPMENT

- Temperature range 0 °C to 70 °C (without illumination)
- Temperature range 5 °C to 60 °C (with illumination)
- Variable positionable illumination cassettes, each with 5 daylight fluorescent illumination tubes
- MP controller with two programs with 10 sections each, alternatively switchable to one program with 20 sections
- Integrated weekly program timer with real-time function
- Digital temperature setting with an accuracy of a tenth of a degree
- Adjustable ramp function via program editor
- Adjustable fan speed
- Elapsed time indicator
- Access port Ø 30 mm, left side
- Inner glass door
- Independent temperature safety device class 3.1 (DIN 12880) with visual and audible alarm
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Adjustable intervals for printer
- Two racks, stainless steel



## ► KBW SERIES | BEST GROWTH CONDITIONS:



### Precise temperature conditions

- APT.line™ preheating chamber
  - Homogeneous temperature conditions
  - Constant and gentle air circulation
  - Small temperature gradients of  $\pm 0.1$  K
- DCT™ cooling system
  - Minimized risk of evaporation
  - Large capacity reserves for lighting applications



### Wide range of programming options

Multifunctional MP controller

- Two program with 10 sections each
  - Variable positionable illumination cassettes
  - Tenths of a degree accuracy, precise temperature setting
  - Standard RS 422 interface




### Homogeneous light distribution

- Flexible positioning of illumination cassettes
- Fully removable
- Three different light spectra, individually selectable



### Comprehensive additional services

- Data Logger Kit
- Years of proven and recognized validation and documentation materials
- Customer-specific modifications

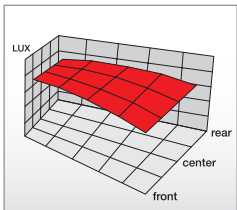

- Special reflector geometries guarantee a uniform illumination of the interior

## ► OPTIONS

- Access ports of various diameters with silicone plug
- Shelves, stainless steel
- Perforated shelf
- Reinforced rack with additional mounting
- Door lock
- FLUORA® growth fluorescent tube set
- Arabidopsis fluorescent tubes set
- Waterproof interior socket 230 V AC
- Ethernet interface
- Additional PT 100 temperature sensor
- Analog output for temperature 4 – 20 mA with 6-pin DIN socket
- Zero-voltage relay outputs accessible via 6-pin DIN socket
- Temperature measurement according to DIN 12880
- Independent temperature safety device class 3.3 (DIN 12880)
- Calibration certificate and extension to calibration certificate
- Data Logger Kits and Logger Software



FLUORA® or Arabidopsis illumination options



Data Logger Kits



Inner chamber socket



Growth chamber with viewing window

► **BINDER  
INDIVIDUAL**  
More information  
on page 133

# Technical data for your planning and installation

## KBW series



	KBW 240	KBW 400	KBW 720
<b>Exterior dimensions</b>			
Width (including access port) (mm/inch)	930 / 36.61	930 / 36.61	1255 / 49.41
Height (including castors) (mm/inch)	1460 / 57.48	1945 / 76.57	1925 / 75.79
Depth (mm/inch)	800 / 31.50	800 / 31.50	890 / 35.04
including door handle, I-panel, connection (mm/inch)	880 / 34.65	880 / 34.65	970 / 38.19
Quantity of doors	1	1	2
<b>Interior dimensions</b>			
Width (mm/inch)	650 / 25.59	650 / 25.59	970 / 38.19
Height (mm/inch)	785 / 30.91	1270 / 50.00	1250 / 49.21
Depth (mm/inch)	485 / 19.09	485 / 19.09	576 / 22.68
Interior volume (l/cu.ft.)	247 / 8.73	400 / 14.13	698 / 24.66
Quantity of racks (standard/max.)	2 / 9	3 / 12	3 / 12
Quantity of light cassettes (standard/max.)	2	3	3
Load per rack (kg/lbs.)	30 / 66	30 / 66	45 / 99
Permitted total load (kg/lbs.)	100 / 221	120 / 265	150 / 331
Weight (empty) (kg/lbs.)	202 / 446	267 / 589	377 / 832
<b>Temperature data</b>			
Temperature range <sup>1)</sup> (°C/°F)	0 – 70 / 32 – 158	0 – 70 / 32 – 158	0 – 70 / 32 – 158
Temperature uniformity (±K)	0.5	0.5	0.5
Temperature fluctuation (±K)	0.1	0.1	0.1
<b>Temperature data with 100 % light intensity</b>			
Temperature range <sup>1)</sup> (°C/°F)	5 – 60 / 41 – 140	5 – 60 / 41 – 140	5 – 60 / 41 – 140
Temperature uniformity (±K)	0.5	1.5	1.0
Temperature fluctuation (±K)	0.1	0.1	0.1
<b>Electrical data</b>			
Housing protection acc. to EN 60529	20	20	20
Nominal voltage (±10 %) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~	200 – 240 1N~
Nominal power (kW)	1.4	1.6	2.7
Energy consumption with 100 % light intensity <sup>2)</sup> at 25 °C / 77 °F (Wh/h)	495	660	820
<b>Model no.</b>			
	9020-0164	9020-0180	9020-0172

<sup>1)</sup> Lower values are valid up to an ambient temperature of max. 25 °C / 77 °F // <sup>2)</sup> These values can be used upon calculation of air conditioning systems. // <sup>3)</sup> Average value, measured with a spherical sensor 12 cm / 4.7 in below the light cassette. The values given in W/m² refer to global radiation // All technical data is specified for units with standard equipment at an ambient temperature of 25 °C / 77 °F and a mains voltage fluctuation of +/-10 %. The temperature data is determined in accordance with factory standard following DIN 12880, respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.




Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)






## ► DIMENSIONS

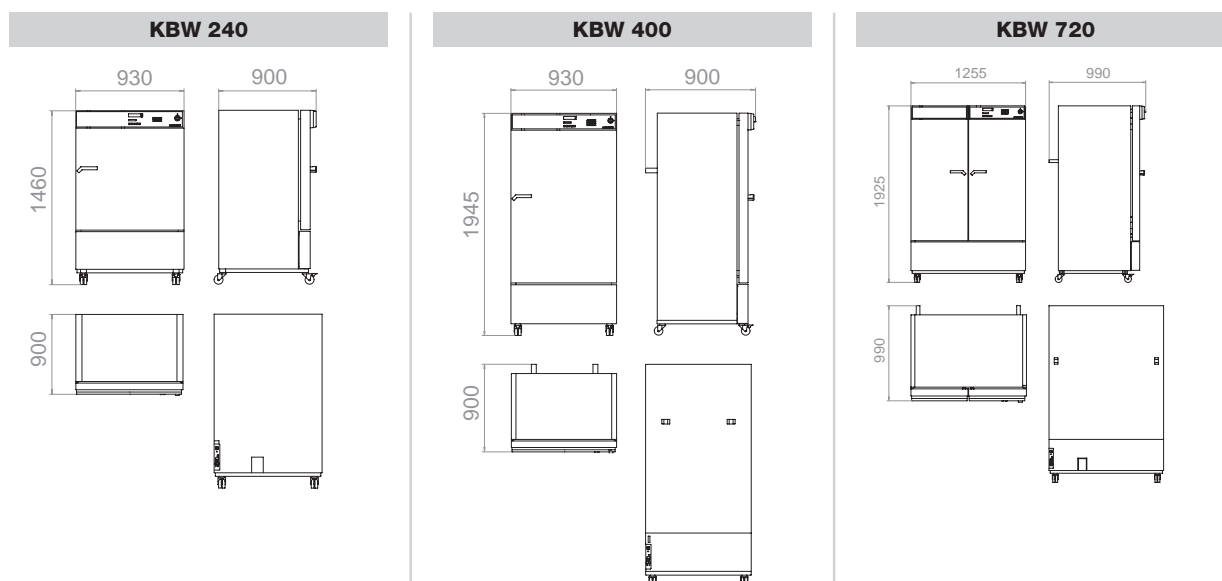

## ► INSTALLATION REQUIREMENTS

	KBW 240	KBW 400	KBW 720
Wall clearance rear (minimum) (spacers) (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94
Wall clearance side (minimum) (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94
Nominal voltage ( $\pm 10\%$ ) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~	200 – 240 1N~
Nominal power (kW)	1.4	1.6	2.7

## ► LIGHT SPECTRA

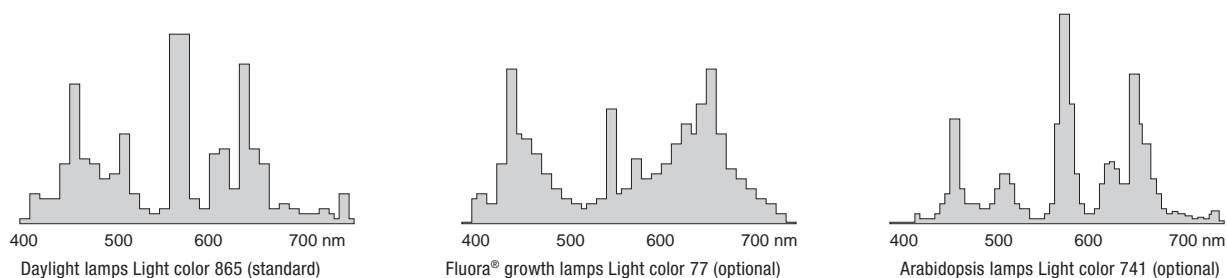

Illumination data per light cassette	KBW 240	KBW 400	KBW 720
Daylight tubes <sup>3</sup> (Lux / W/m <sup>2</sup> )	10.000 / 36	10.000 / 31	14.200 / 38
Fluora® growth lamps <sup>3</sup> (Lux / W/m <sup>2</sup> )	6.500 / 34	7.500 / 28	10.500 / 31
Arabidopsis lamps <sup>3</sup> (Lux / W/m <sup>2</sup> )	10.000 / 40	11.000 / 35	14.400 / 38

# Growth chambers with illumination and humidity KBWF series

Through its versatile programming options, KBWF series achieves the perfect combination of temperature, humidity and light. The wide climate range enables any climate simulation precisely and consistently, over long periods of time, including programming for day/night cycles.



► KBWF 720 model



240

720

Available sizes (liters)

## ► EQUIPMENT

- Temperature range: 0 °C to 70 °C (without illumination and humidity)
- Temperature range: 10 °C to 60 °C (with illumination and humidity)
- Humidity range without illumination: 10 % to 80 % r. H.
- Humidity range with illumination: 10 % to 75 % r. H.
- Two variable positionable illumination cassettes, each with 5 daylight fluorescent illumination tubes
- MCS controller for temperature, humidity, and lighting control which can store 25 programs of 100 sections each, for a maximum of 500 program segments (for programming of day/night cycles)
  - Integrated electronic chart recorder
  - Variety of options for the graphic display of process parameters
  - Real-time clock
- Controlled humidification and dehumidification system with capacitive humidity sensor
- Independent temperature safety device class 3.1 (DIN 12880) with visual and audible temperature alarm
- Access port Ø 30 mm, left side
- Inner glass door
- Complete safety connection kit for water supply and drainage, including water hose, (total length 6 m)
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Two racks, stainless steel

## ► KBWF SERIES | BEST GROWTH CONDITIONS:



### Precise temperature conditions

- APT.line™ preheating chamber
  - Homogeneous temperature conditions
  - Constant and gentle air circulation
  - Small temperature gradients of  $\pm 0.1K$
- DCT™ cooling system
  - Minimized risk of evaporation
  - Large capacity reserves for lighting applications



### Individual program control

MSC controller color screen

- Programmable day and night cycles
- Variety of options for the graphic display of process parameters
- Standard RS 422 interface




### Homogeneous light distribution

- Flexible positioning of illumination cassettes
- Fully removable
- Three different light spectra, individually selectable



### Comprehensive additional services

- Data Logger Kit
- Years of proven and recognized validation and documentation materials
- Customer-specific modifications

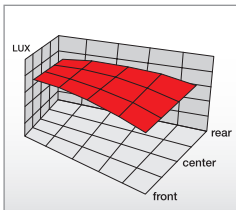

- Special reflector geometries guarantee a uniform illumination of the interior



### Fast and precise humidification

- Vapor pressure humidification with fast response times
- Drift-free, capacitive humidity sensor
- Short recovery times after door opening
- Finely adjustable humidity control

## ► OPTIONS

- Access ports of various diameters with silicone plug
- Shelves, stainless steel
- Perforated shelf
- Reinforced rack with additional mounting
- Door lock
- FLUORA® growth fluorescent tube set
- Arabidopsis fluorescent tubes set
- Ethernet interface
- Additional PT 100 temperature sensor
- Analog output for temperature 4 – 20 mA with 6-pin DIN socket
- Zero-voltage relay outputs accessible via 6-pin DIN socket
- Lockable controller keyboard
- Ethernet interface
- External water supply set
- BINDER PURE AQUA SERVICE with single-use cartridge
- Measurement of temperature accuracy according to DIN 12880
- Calibration certificate and extension to calibration certificate
- Data Logger Kits and Logger Software




FLUORA® or Arabidopsis illumination options



BINDER PURE AQUA SERVICE



External water supply set

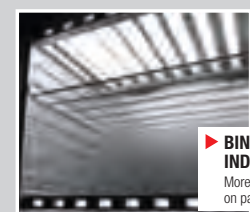

KBWF 720 with increased illumination

► **BINDER  
INDIVIDUAL**  
More information  
on page 133


# Technical data for your planning and installation

## KBWF series




	KBWF 240	KBWF 720
<b>Exterior dimensions</b>		
Width (including access port) (mm/inch)	930 / 36.61	1225 / 49.41
Height (including castors) (mm/inch)	1460 / 57.48	1925 / 75.79
Depth (mm/inch)	800 / 31.50	890 / 35.04
Depth including door handle, I-triangle, connection, cable (mm/inch)	880 / 34.65	970 / 38.19
Number of doors	1	2
<b>Interior dimensions</b>		
Width (mm/inch)	650 / 25.59	973 / 38.31
Height (mm/inch)	785 / 30.91	1250 / 49.21
Depth (mm/inch)	485 / 19.09	576 / 22.68
Interior volume (l/cu.ft.)	247 / 8.73	700 / 24.73
Number of racks (standard/max.)	2 / 7	3 / 12
Number of light cassettes	2	3
Load per rack (kg/lbs.)	30 / 66	45 / 99
Permitted total load (kg/lbs.)	100 / 221	150 / 331
Weight (empty) (kg/lbs.)	214 / 472	374 / 826
<b>Temperature data (without humidity)</b>		
Temperature range without light cassettes <sup>1)</sup> (°C/°F)	0 – 70 / 32 – 158	0 – 70 / 32 – 158
Temperature range with light cassettes, with illumination <sup>1)</sup> (°C/°F)	10 – 60 / 50 – 140	10 – 60 / 50 – 140
<b>Climatic data</b>		
Temperature range without light cassettes <sup>1)</sup> (°C/°F)	0 – 70 / 32 – 158	0 – 70 / 32 – 158
Temperature range with light cassettes, with illumination <sup>1)</sup> (°C/°F)	10 – 60 / 50 – 140	20 – 60 / 68 – 140
Temperature uniformity with illumination (±K)	0.5 – 1.0	1.0 – 1.5
Temperature fluctuation with illumination (±K)	0.1 – 0.5	0.1 – 0.5
Humidity range without light cassettes (% r. H.)	10 – 80	10 – 80
Humidity range with light cassettes, with illumination <sup>1)</sup> (°C/°F)	10 – 75	10 – 75
Humidity fluctuation with illumination (± % r. H.)	2	≤ 2.5
<b>Electrical data</b>		
Housing protection acc. to EN 60529	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~
Nominal power at 240 V (kW)	2.4	3.5
Energy consumption <sup>3)</sup> at 37 °C / 98.6 °F / 75 % r. H. (Wh/h)	600	2350
<b>Model no.</b>		
	9020-0166	9020-0173

► **TEMPERATURE-HUMIDITY CHART  
KBWF 240 WITH LIGHT**

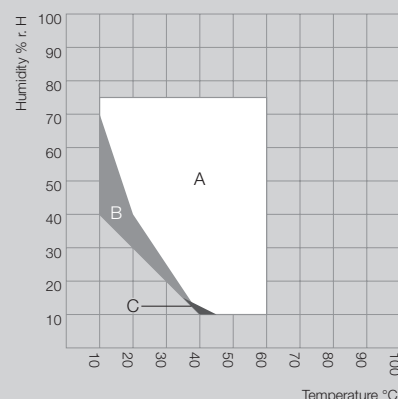

► **TEMPERATURE-HUMIDITY CHART  
KBWF 720 WITH LIGHT**

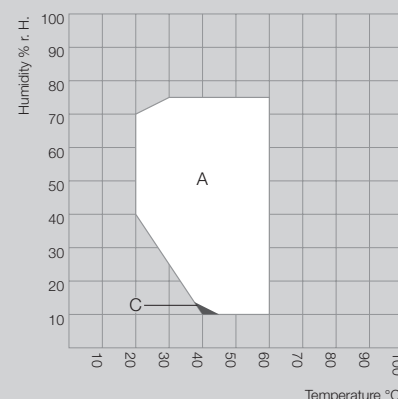


A: Guaranteed condensation free range  
B: Temporary usable operation (up to 24 h)  
C: Deviations of technical data may be possible

<sup>1)</sup> Lower values are valid up to an ambient temperature of max. 25 °C / 77 °F // <sup>2)</sup> Average value, measured with a spherical sensor (±10 %) by 12 cm / 4.7 in below the light cassette. The values given in W/m<sup>2</sup> refer to global radiation. // <sup>3)</sup> Use this value for dimensioning air conditioning systems /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with BINDER factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)


## ► DIMENSIONS

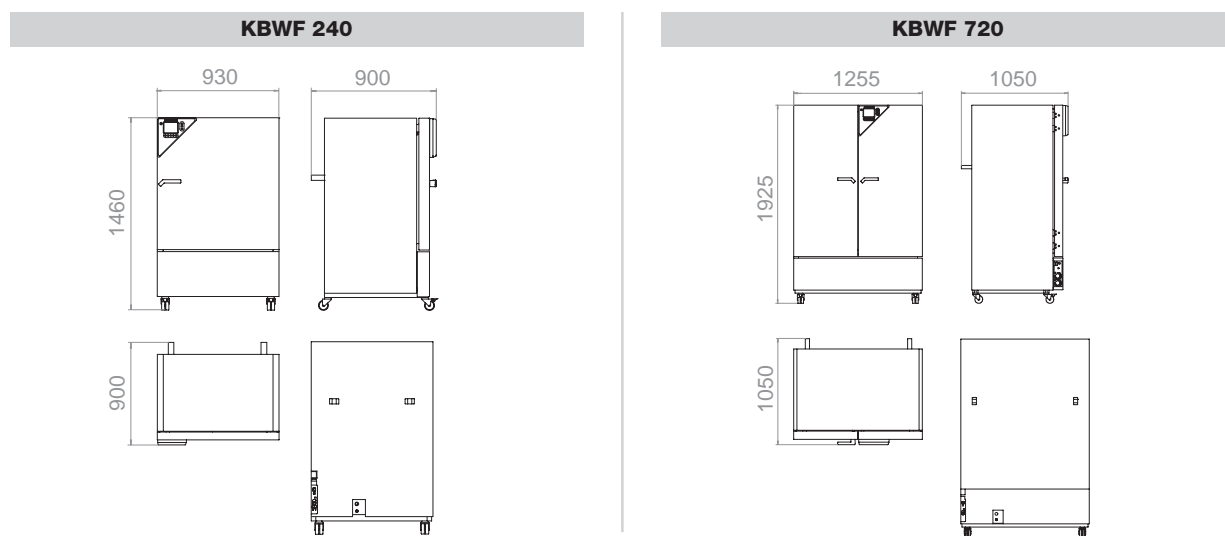

## ► INSTALLATION REQUIREMENTS

	KBWF 240	KBWF 720
Wall clearance rear (minimum) (spacer) (mm/inch)	100 / 3.94	100 / 3.94
Wall clearance side (minimum) (mm/inch)	100 / 3.94	200 / 7.87
Nominal voltage ( $\pm 10\%$ ) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~
Nominal power at 240 V (kW)	2.4	3.5

## ► LIGHT SPECTRA

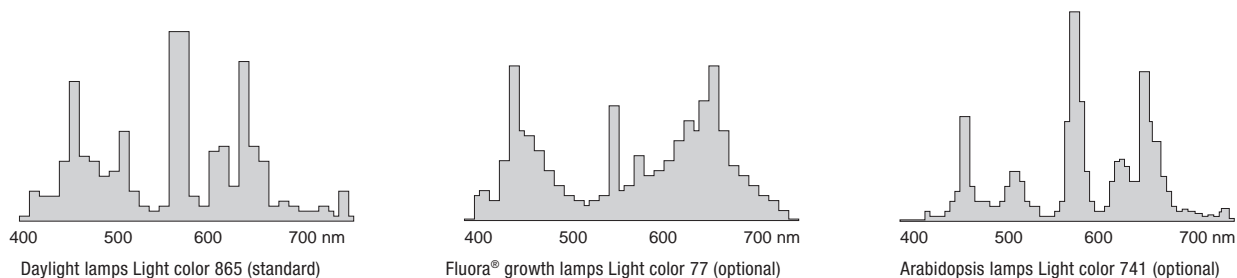

Illumination data per light cassette	KBWF 240	KBWF 720
Daylight tubes <sup>2)</sup> (Lux / W/m <sup>2</sup> )	10.000 / 36	14.200 / 38
Fluora® growth lamps <sup>2)</sup> (Lux / W/m <sup>2</sup> )	6.500 / 34	10.500 / 31
Arabidopsis lamps <sup>2)</sup> (Lux / W/m <sup>2</sup> )	10.000 / 40	14.400 / 38





# Storage

## | Always the right concept

An ultra low temperature freezer must reliably cool to maintain the value of the samples. This crucial laboratory requirement is supplemented by the necessity of a practical and user-oriented safety and control concept, e. g. by the effortless traceability of the temperature profile throughout the storage period or opening of the chamber easily.

BINDER rose to this challenge and developed a revolutionary ULT freezer that sets new security standards and fits seamlessly in BINDER's products range dedicated to the scientific laboratory.



**Sample storage**  
UF V





# Sample storage

## UF V series

### Ultra safe: BINDER Ultra low temperature freezer

- ▶ Ultra innovative
  - ▶ Personalized access control with RFID technology
  - ▶ Convenient RFID read-out of all door opening events
  - ▶ 24-hour exchange service in case of emergency
- ▶ Ultra simple
  - ▶ Door opens easily at the push of a button
  - ▶ Toolless filter change
  - ▶ Simple data read-out through a data logger with USB port
- ▶ Ultra efficient
  - ▶ Vacuum insulation panels
  - ▶ Optimal sample capacity for up to 52.800 samples

Safety without sacrificing reliability: true to this maxim, BINDER ultra low temperature freezers offer maximum process reliability and ease of use through innovative solutions. Our product includes everything you could wish for in sample storage: from reliable operation to user-friendly and system integration, to emergency exchange service.



Biological samples



Pharmaceutical agents




Blood banks

# -86 °C Ultra low temperature freezers

## UF V series

The new BINDER ultra low temperature freezers ensure safe long-term storage of samples at -86 °C. The UF V series is based on a multi-layer security concept, meeting any laboratory staff's needs beyond expectations.

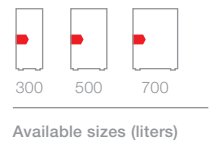

### ► EQUIPMENT

- Temperature range: -40 °C to -86 °C
- GUARD.CONTROL™ automatic door mechanism for easy opening and closing
- Personalized access control through RFID technology
- VIP.COMPLETE™ internal large-surface vacuum insulation panels (VIP) on all 6 sides
- Two-stage cascade refrigeration with hermetically-sealed compressors
- Non-flammable refrigerant, CFC-free
- Heated duplex door gasket prevents icing
- User-friendly microprocessor controller with LED display
- Alarm test system
- Integrated visual and audible alarm system: in case of power failure, temperature alarm or battery failure
- Battery back-up system: Provides full functionality of alarms in case of power failure for 72 hours
- RS 422 interface
- Zero-voltage relay alarm output
- Analog output 4 – 20 mA
- Inner chamber made of high-quality stainless steel
- Robust, lockable casters
- Easy front access for air filter cleaning
- SERVICE.PRO™ 2-year warranty and 4-year extended warranty

## ► UF V SERIES | MAXIMUM PROCESS SAFETY:



### Safe and convenient operation

#### GUARD.CONTROL™

- Personalized access control with RFID key card
- Complete user log
- Simple data read-out through RFID reader
- Door mechanism for easy opening and closing at the push of a button



### Reliable back-up service

#### SERVICE.PRO™

- 24-hour emergency exchange service (currently valid in G, F and GB)
- 24/7 Hotline service after online registration
- 2+4 (2 year warranty and 4-year extended warranty)



### Simple data management

#### DATA.SECURE™

- Temperature recording via data logger with USB connection
- Ease of integration into the security / alarm infrastructure
- Integrated PT 100 for independent data recording



### Reliable refrigeration

#### VIP.COMPLETE

- Large-surface insulation with Vacuum Insulation Panels
- Maximum interior space utilization thanks to the narrow design (up to 52.800 samples)



### Easy maintenance and cleaning

- Easy front access for air filter cleaning
- Defrosting kit facilitates the de-icing process

## ► OPTIONS

- Access port 30 mm, rear panel
- Sample divider
- Cardboard cryo boxes, 50 mm, 75 mm without dividers
- Cardboard divider for cryo boxes, 49/64/81/100 cells
- Polypropylene cryo boxes, 50 mm / 75 mm with 81-cells and codification
- Various side access racks, for 50 and 75 mm boxes
- Various sliding drawer racks UF V, for 50 and 75 mm boxes
- CO<sub>2</sub> back-up system
- Ethernet interface
- Additional PT 100 sensor with LEMO socket
- Remote alarm GSM box
- Advanced voltage booster
- RFID reader
- PDF Data Logger with USB port
- Circular chart recorder
- Defrosting kit including de-icing tray, wiper and scraper
- Water cooling
- Deep temperature gloves
- Magnetic pouch



Temperature recording with PDF Data Logger via USB



Wide selection of racks and boxes



Defrosting kit (de-icing tray, wiper and scraper)



**BINDER  
INDIVIDUAL**  
More information  
on page 133

Solutions customized to your requirements




# Technical data for your planning and installation

## UF V series




	UF V 300	UF V 500	UF V 700
<b>Exterior dimensions</b>			
Width (mm/inch)	755 / 29.72	900 / 35.43	1200 / 47.24
Height (mm/inch)	1970 / 77.56	1970 / 77.56	1970 / 77.56
Depth including controller, connection (mm/inch)	935 / 36.81	935 / 36.81	935 / 36.81
Wall clearance rear (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94
Wall clearance side (side without hinges) (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94
Wall clearance side (side with hinges) (mm/inch)	245 / 9.65	245 / 9.65	245 / 9.65
<b>Interior dimensions</b>			
Width of interior (mm/inch)	473 / 18.62	619 / 24.37	911 / 35.87
Width of shelf (mm/inch)	442 / 17.40	588 / 23.15	880 / 34.65
Height of interior (mm/inch)	1300 / 51.18	1300 / 51.18	1300 / 51.18
Depth of interior / Depth of shelf (mm/inch)	600 / 23.62	600 / 23.62	600 / 23.62
Interior volume, total (l/cu.ft.)	345 / 12.2	460 / 16.25	685 / 24.20
Number of shelves	3	3	3
Number of compartments / compartment doors	4 / 4	4 / 4	4 / 4
Permitted load per shelf (kg/lbs.)	40 / 88	50 / 110	65 / 143
Permitted total load (kg/lbs.)	160 / 353	200 / 442	260 / 574
Weight of the unit (empty) (kg/lbs.)	300 / 662	320 / 706	360 / 795
<b>Temperature data</b>			
Setting range °C/°F	-90/-40/-130/-40	-90/-40/-130/-40	-90/-40/-130/-40
Measuring range, ambient up to °C/°F	-90 / -130	-90 / -130	-90 / -130
Average temperature uniformity (±K)	2.5	2.5	2.5
Pull-down time from +25 °C / 77 °F to -80 °C / -112 °F <sup>1)</sup> (hours)	9	9	9
Pull-up time in case of power failure from -80 °C / -122 °F to -60 °C / -76 °F <sup>2)</sup> (hours)	3	3	3
<b>Electrical data UF V</b>			
Housing protection acc. to EN 60529	20	20	20
Nominal voltage (+10 %) 50 Hz (V)	230	230	230
Energy consumption at -80 °C / -112 °F with an ambient temperature of +25 °C / 77 °F (kWh/day)	12	13.9	16.2
Average heat dissipation at set-point -80 °C / -112 °F with an ambient temperature of +25 °C / 77 °F <sup>2)</sup> (W)	500	580	680
<b>Model no. UFV with right hinged door</b>			
	9020-0165	9020-0232	9020-0233
<b>Model no. UFV with left hinged door</b>			
	9020-0214	9020-0213	9020-0138

### ► COOLING DOWN TIME / DEFROST TIME AFTER POWER FAILURE

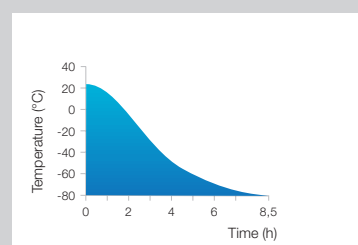

Cooling down time without loading at 25 °C ambient temperature

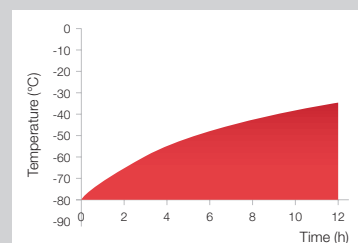

Defrost time without loading at 25 °C ambient temperature

<sup>1)</sup> to 98 % of the set value // <sup>2)</sup> Use this value for sizing air conditioning systems // <sup>3)</sup> pay attention to wall clearance /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.




Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)




► LOADING VARIANTS (DRAWER RACKS AND BOXES)


UF V 300

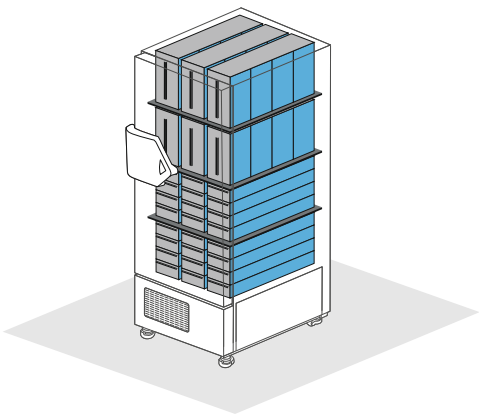

UF V 500

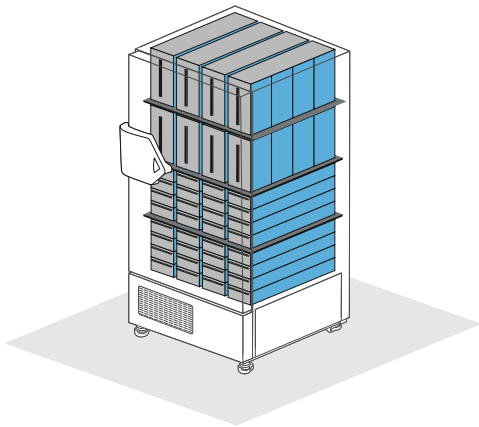

UF V 700

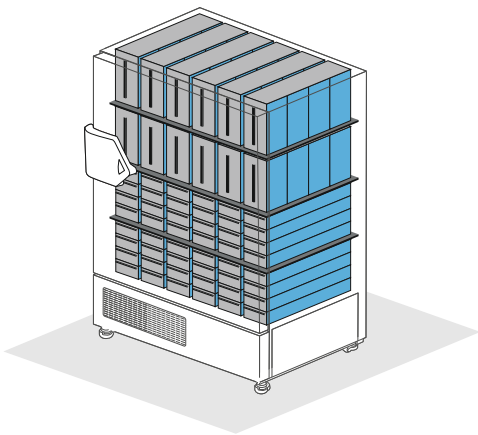

	UF V 300	UF V 500	UF V 700
Racks per compartment	3	4	6
Racks per chamber	12	16	24
Cryo boxes 50 mm (max. no.)	264	352	528
Cryo boxes 75 mm (max. no.)	168	224	336






# Drying and tempering

## | Always the right temperature

Test chambers used in scientific laboratories, must achieve temperature stress testing requirements far beyond temperature accuracy, heat distribution and residue-free drying. This is the case with a chamber used to dry flammable materials that must meet safety standards.

BINDER masters these requirements like no other and offers you a wide range of drying and heating ovens, as well as material test chambers. Choose from ovens with gravity (natural) or mechanical (forced) convection for safety or vacuum drying. BINDER INDIVIDUAL meets your special requirements for these products as well.

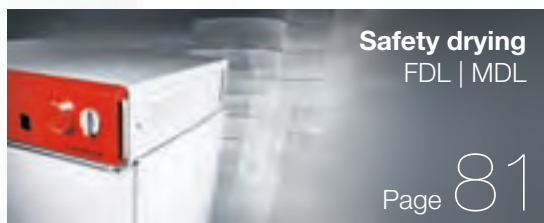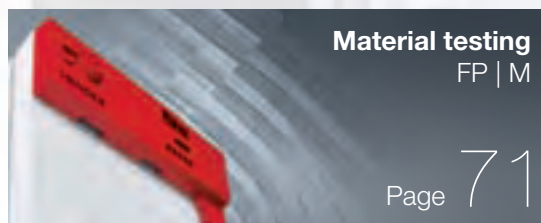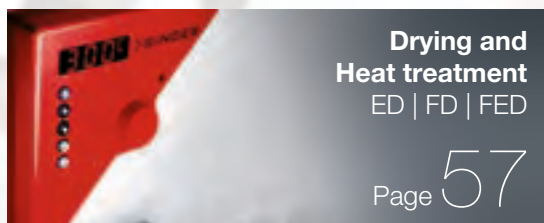





# Drying and Heating ovens

## ED | FD | FED series

### Unsurpassed efficiency: BINDER Drying and Heating ovens

- ▶ Unsurpassed precision
  - ▶ Wide temperature range of 5 °C above ambient temperature to 300 °C
  - ▶ Fast, uniform tempering
  - ▶ Identical test conditions throughout the chamber interior
  - ▶ High standard according to DIN 12880 (27-point measurement)
- ▶ Unsurpassed reliability
  - ▶ Durable construction based on corrosion-resistant materials
  - ▶ Large capacity reserves
  - ▶ Short heating up and recovery times
  - ▶ Material temperature protection through an independent temperature safety device Class 2
- ▶ Unsurpassed versatility
  - ▶ Available with gravity (natural) and mechanical (forced) convection
  - ▶ Digital multifunctional controller with advanced timer functions
  - ▶ Communication interface
  - ▶ Adjustable front ventilation flap slide

Whether by gravity (natural) or mechanical (forced) convection, our drying and heating ovens provide high standards of quality and process stability. The wide temperature range of 5 °C above ambient temperature to 300 °C allows short heating up times and provide large capacity reserves. The excellent thermal insulation also reduces operating costs.



Glass drying



Sterilization



Plastics tempering



# Drying ovens with gravity convection

## ED series

Routine drying and sterilization applications up to 300 °C and precision heat storage are the strength of ED drying ovens. Drying processes run much more efficiently with gravity (natural) convection and the high rate of air exchange.



► ED 53 model



Available sizes (liters)

### ► EQUIPMENT


- Temperature range from 5 °C above ambient temperature to 300 °C
- DS control with integrated timer 0 to 99 hrs
- Digital temperature setting with an accuracy of one degree
- Independent adjustable temperature safety device class 2 (DIN 12880), with visual temperature alarm
- Adjustable ventilation by means of front ventilation flap slide and rear exhaust Ø 50 mm
- Optional RS 422 Interface for APT-COM™ Data Control System communication software
- Units up to 115 liters are stackable
- Two chrome-plated racks included

## ► ED SERIES | BEST DRYING RESULTS:




### Uniform drying conditions

- APT.line™ preheating chamber
  - Homogeneous temperature control with gravity (natural) convection
  - Identical test conditions throughout the chamber interior independent of sample size and quantity

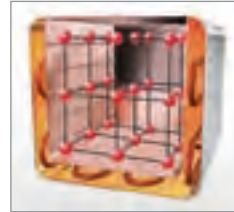

### Best quality and precision guaranteed

- High standard according to DIN 12880 (27-point measurement according to factory standard)
- Short delivery times
- Minimal maintenance and operating costs

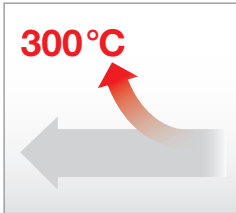

### Broad range of applications

- Standard temperature range up to 300 °C
- Large capacity reserves
- Short heating up times



### Convenient work environment

- Hermetic door closure with 2-points door closure
- Low heat dissipation due to 60 mm insulation
- Rack with tilt protection for easy loading and unloading
- Complete stainless steel inner chamber without permanent fixtures
- Easy cleaning

## ► OPTIONS

- Access ports of various diameters with silicone plugs
- Racks, chrome-plated or stainless steel
- Perforated shelf, stainless steel
- Independent temperature safety device class 3.1 according to DIN 12880
- Door with window and interior lighting
- Door lock
- Door gasket made of FKM (Viton)
- Switchable audio alarm for overheating
- Analog output for temperature 4 – 20 mA with 6-pin DIN socket
- Temperature measurement according to DIN 12880
- Calibration certificate
- Extension to factory calibration certificate (additional measuring point)
- Data Logger Kits and Logger Software



Access ports with silicone plugs



Calibrations and validations



Door with window and interior lighting



Drying oven with special door access port for loading test materials

► **BINDER  
INDIVIDUAL**  
More information  
on page 133

# Technical data for your planning and installation

## ED series



	ED 23	ED 53	ED 115	ED 240	ED 400	ED 720
<b>Exterior dimensions</b>						
Width (mm/inch)	435 / 17.13	635 / 25.0	835 / 32.87	1035 / 40.75	1235 / 48.62	1235 / 48.62
Height (including feet/castors) (mm/inch)	495 / 19.49	620 / 24.41	705 / 27.76	825 / 32.48	1025 / 40.35	1530 / 60.24
Depth (mm/inch)	520 / 20.47	575 / 22.64	645 / 25.39	745 / 29.33	765 / 30.12	865 / 34.06
plus door handle, I-panel and exhaust duct (mm/inch)	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13
Quantity of doors	1	1	1	2	2	2
<b>Interior dimensions</b>						
Width (mm/inch)	222 / 8.74	400 / 15.75	600 / 23.62	800 / 31.50	1000 / 39.37	1000 / 39.37
Height (mm/inch)	330 / 12.99	400 / 15.75	480 / 18.90	600 / 23.62	800 / 31.50	1200 / 47.24
Depth (mm/inch)	300 / 11.81	350 / 13.78	420 / 16.54	520 / 20.47	520 / 20.47	620 / 24.41
Interior volume (l/cu.ft.)	20 / 0.7	53 / 1.9	115 / 4.1	240 / 8.6	400 / 14.3	720 / 25.7
Quantity of racks (standard/max.)	2 / 3	2 / 5	2 / 6	2 / 7	2 / 10	2 / 15
Load per rack (kg/lbs.)	12 / 26	15 / 33	20 / 44	30 / 66	35 / 77	45 / 99
Permitted total load (kg/lbs.)	25 / 55	40 / 88	50 / 110	70 / 155	90 / 199	120 / 265
Weight of the unit (empty) (kg/lbs.)	22 / 49	42 / 93	57 / 126	86 / 190	125 / 276	174 / 384
<b>Temperature data</b>						
Temperature range, by 5 °C/9 °F above ambient up to °C/°F	300 / 572	300 / 572	300 / 572	300 / 572	300 / 572	300 / 572
Temperature uniformity <sup>1)</sup> at 150 °C/302 °F (±K)	2.5	3.2	2.5	2.5	3	2.8
Temperature fluctuation at 70 °C/158 °F (±K)	0.3	0.3	0.3	0.3	0.3	0.3
Heating-up time <sup>2)</sup> to 150 °C/302 °F (min)	24	27	29	48	62	69
Recov. time after door was opened for 30 sec. <sup>2)</sup> at 150 °C/302 °F (min)	5	6	9	13	20	14
Air change at 150 °C/302 °F (x/h)	13	19	10	10	10	9
<b>Electrical data</b>						
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230 1 N ~	230 1 N ~	230 1 N ~	230 1 N ~	400 3 N ~	400 3 N ~
Nominal power (kW)	0.8	1.2	1.6	2.7	3.4	5.0
Energy consumption at 150 °C/302 °F (Wh/h)	148	210	300	447	672	750
Model no. without RS 422	9010-0190	9010-0078	9010-0096	-	-	-
Model no. with RS 422	9010-0191	9010-0079	9010-0097	9010-0101	9010-0075	9010-0076

<sup>1)</sup> value without window // <sup>2)</sup> to 98 % of the set value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %.


The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber.

All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)

► DIMENSIONS

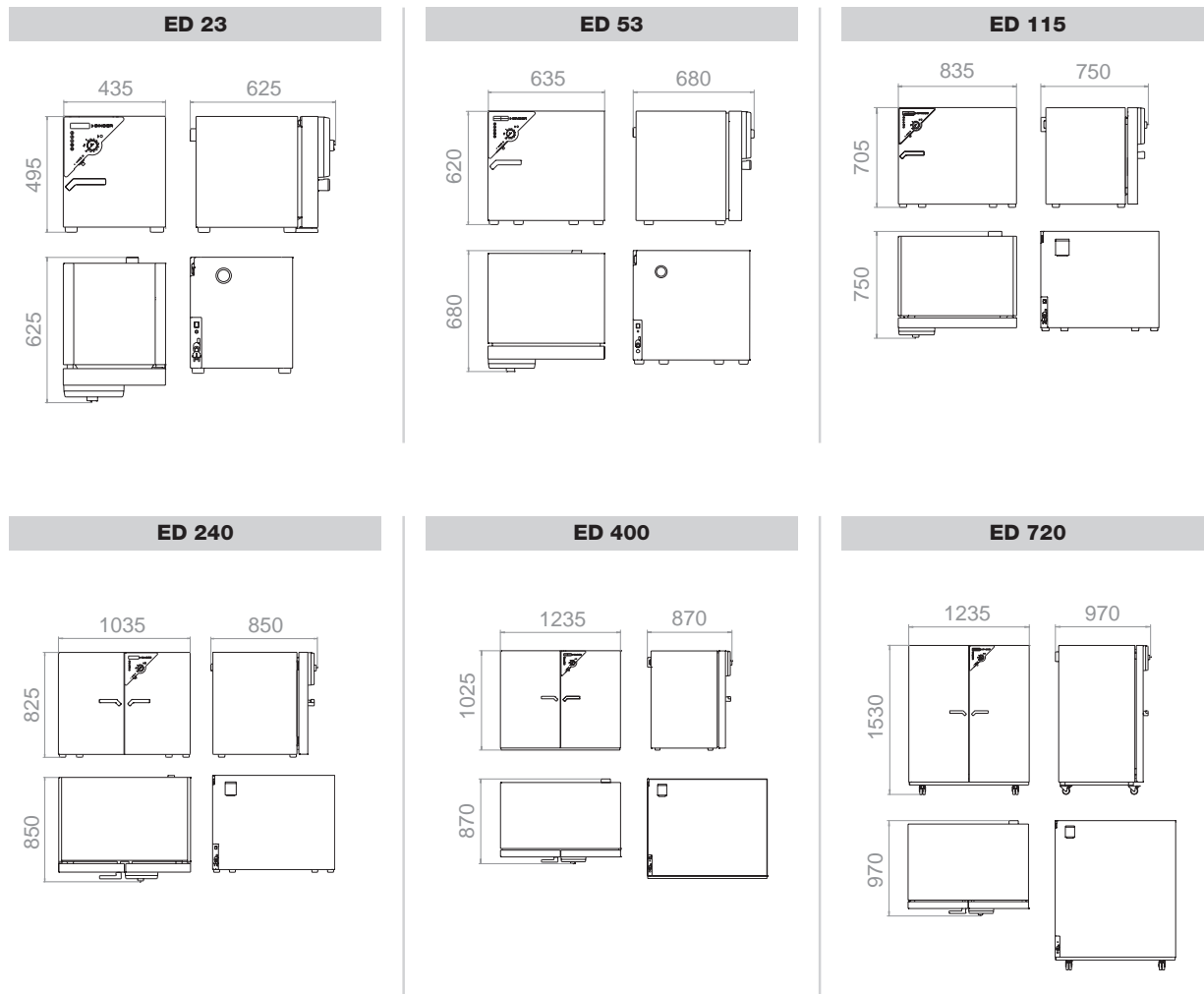

► INSTALLATION REQUIREMENTS

	ED 23	ED 53	ED 115	ED 240	ED 400	ED 720
Nominal voltage ( $\pm 10\%$ ) 50/60 Hz (V)	230 1 N ~	230 1 N ~	230 1 N ~	230 1 N ~	400 3 N ~	400 3 N ~
Nominal power (kW)	0.8	1.2	1.6	2.7	3.4	5

# Drying ovens with mechanical convection

## FD series

The FD series is the optimal choice when fast drying and sterilization are required. With fully homogeneous temperature distribution, quick dynamics and powerful fan, the FD saves valuable time.



► FD 115 model  
with option window



### ► EQUIPMENT


- Temperature range from 5 °C above ambient temperature to 300 °C
- DS control with integrated timer 0 to 99 hrs
- Digital temperature setting with an accuracy of one degree
- Independent adjustable temperature safety device class 2 (DIN 12880), with visual temperature alarm
- Adjustable ventilation by means of front ventilation flap slide and rear exhaust Ø 50 mm
- Units up to 115 liters are stackable
- Two chrome-plated racks included

## ► FD SERIES | BEST DRYING RESULTS:




### Uniform drying conditions

- APT.line™ preheating chamber
  - Homogeneous temperature control with mechanical (forced) convection
  - Identical test conditions throughout the chamber interior independent of sample size and quantity

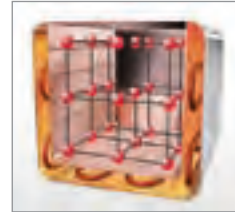

### Best quality and precision guaranteed

- High standard according to DIN 12880 (27-point measurement)
- Short delivery times
- Minimal maintenance and operating costs

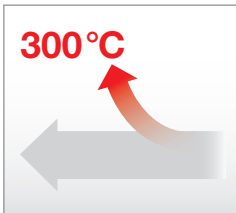

### Broad range of applications

- Standard temperature range up to 300 °C
- Large capacity reserves
- Short heating up times




### Convenient work environment

- User-friendly microprocessor control
- Hermetic door closure with 2-points door closure
- Low heat dissipation due to 60 mm insulation
- Rack with tilt protection for easy loading and unloading
- Complete stainless steel inner chamber without permanent fixtures
- Easy cleaning

## ► OPTIONS

- Access ports of various diameters with silicone plugs
- Racks, chrome-plated or stainless steel
- Perforated shelf, stainless steel
- Independent temperature safety device class 3.1 according to DIN 12880
- Door with window and interior lighting
- Door lock
- Door gasket made of FKM (Viton)
- Switchable audio alarm for overheating
- Analog output for temperature 4 – 20 mA with 6-pin DIN socket
- Temperature measurement according to DIN 12880
- Calibration certificate
- Extension to factory calibration certificate (additional measuring point)
- Data Logger Kits and Logger Software

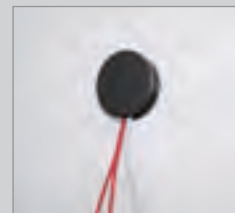

Access ports with silicone plugs



Reinforced shelves



Door with window and interior lighting



Drying oven with partitioned inner chamber and suspensions

► **BINDER  
INDIVIDUAL**  
More information  
on page 133



# Technical data for your planning and installation

## FD series



	FD 23	FD 53	FD 115	FD 240
<b>Exterior dimensions</b>				
Width (mm/inch)	435 / 17.13	635 / 25.0	835 / 32.87	1035 / 40.75
Height (including feet/castors) (mm/inch)	495 / 19.49	620 / 24.41	705 / 27.76	825 / 32.48
Depth (mm/inch)	520 / 20.47	575 / 22.64	645 / 25.39	745 / 29.33
plus door handle, I-panel and exhaust duct (mm/inch)	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13
Quantity of doors	1	1	1	2
<b>Interior dimensions</b>				
Width (mm/inch)	222 / 8.74	400 / 15.75	600 / 23.62	800 / 31.50
Height (mm/inch)	330 / 12.99	400 / 15.75	480 / 18.90	600 / 23.62
Depth (mm/inch)	300 / 11.81	340 / 13.39	410 / 16.14	510 / 20.08
Interior volume (l/cu.ft.)	20 / 0.7	53 / 1.9	115 / 4.1	240 / 8.6
Quantity of racks (standard/max.)	2 / 3	2 / 5	2 / 6	2 / 7
Load per rack (kg/lbs.)	12 / 26	15 / 33	20 / 44	30 / 66
Permitted total load (kg/lbs.)	25 / 55	40 / 88	50 / 110	70 / 155
Weight of the unit (empty) (kg/lbs.)	33 / 73	44 / 97	62 / 137	96 / 212
<b>Temperature data</b>				
Temperature range, 5 °C/9 °F above ambient up to °C/°F	300 / 572	300 / 572	300 / 572	300 / 572
Temperature uniformity <sup>1)</sup> at 150 °C/302 °F (±K)	2.2	2	1.8	2.0
Temperature fluctuation (<= ±K)	0.3	0.3	0.3	0.3
Heating-up time <sup>2)</sup> to 150 °C/302 °F (min)	22	24	28	24
Recov. time after door was opened for 30 sec. <sup>1) 2)</sup> at 150 °C/302 °F (min)	4	5	5	6
Air change at 150 °C/302 °F (x/h)	64	64	32	20
<b>Electrical data</b>				
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230 1 N ~	230 1 N ~	230 1 N ~	230 1 N ~
Nominal power (kW)	0.8	1.2	1.6	2.7
Energy consumption at 150 °C/302 °F (Wh/h)	300	429	544	850
<b>Model no.</b>				
	9010-0194	9010-0082	9010-0102	9010-0104

<sup>1)</sup> value without window // <sup>2)</sup> to 98 % of the set value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %.

The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber.





All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



► DIMENSIONS

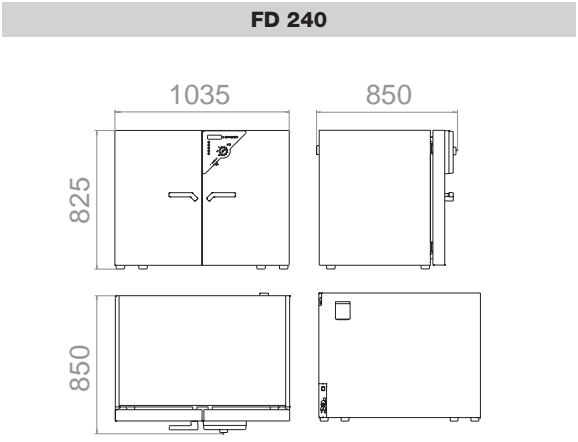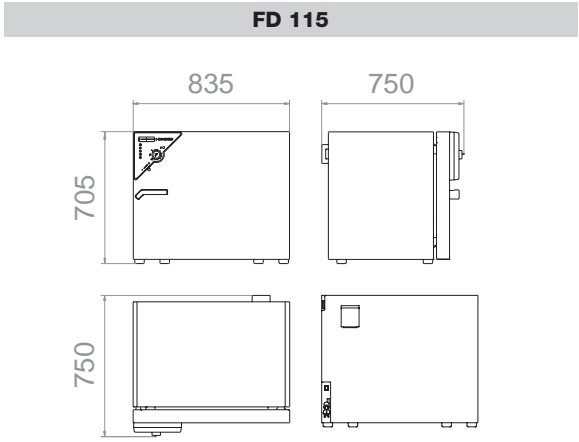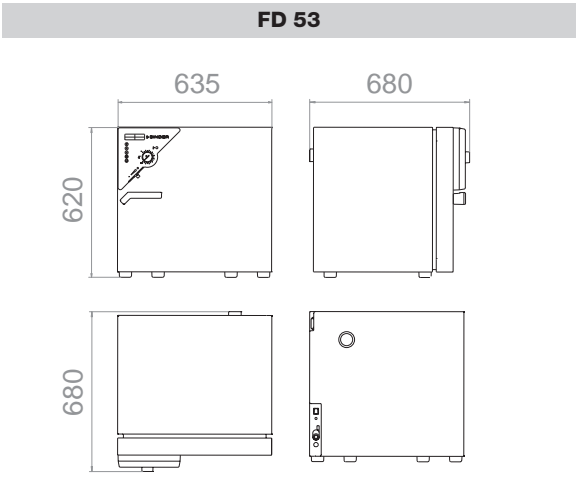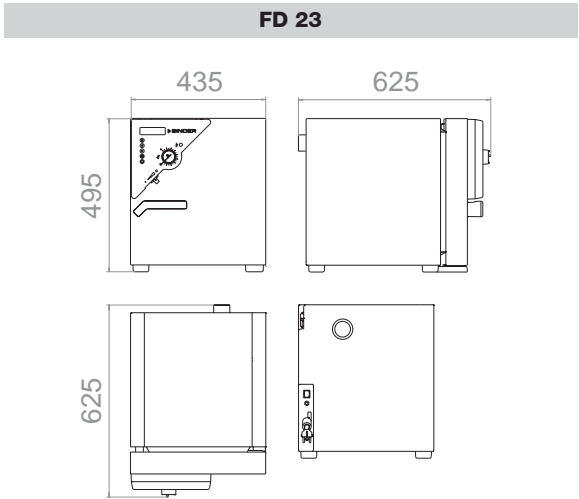

► INSTALLATION REQUIREMENTS


	FD 23	FD 53	FD 115	FD 240
Nominal voltage (±10 %) 50/60 Hz (V)	230 1 N ~	230 1 N ~	230 1 N ~	230 1 N ~
Nominal power (kW)	0.8	1.2	1.6	2.7

# Heating chambers with mechanical convection FED series

Our multi-talented series: almost unlimited capacity and particularly adaptable to the individual requirements of many different tests. With advanced timing functions and a digitally controllable fan, temperature and convection conditions are easily controlled.



► FED 240 model

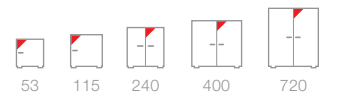

Available sizes (liters)

## ► EQUIPMENT


- Temperature range from 5 °C above ambient temperature to 300 °C
- MS controller with several timer functions
- Controller timer functions: delayed ON, delayed OFF, temperature dependent delayed OFF
- Digital temperature setting with an accuracy of one degree
- Adjustable fan speed
- Adjustable ventilation by means of front ventilation flap slide and rear exhaust Ø 50 mm
- Independent adjustable temperature safety device class 2 (DIN 12880), with visual temperature alarm
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Units up to 115 liters are stackable
- Two chrome-plated racks included

## ► FED SERIES | BEST DRYING RESULTS:



### Uniform drying conditions

- APT.line™ preheating chamber
  - Homogeneous temperature control with mechanical (forced) convection
  - Identical test conditions throughout the chamber interior independent of sample size and quantity

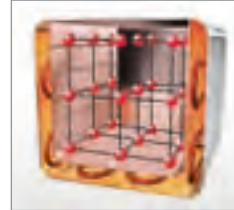

### Best quality and precision guaranteed

- High standard according to DIN 12880 (27-point measurement)
- Short delivery times
- Minimal maintenance and operating costs



### User-friendly controller

- Digital multifunctional controller (MS)
  - Advanced timing functions
  - In the user's field of vision
  - Easy to operate



### Convenient work environment

- User-friendly microprocessor control
- Hermetic door closure with 2-points door closure
- Low heat dissipation due to 60 mm insulation
- Rack with tilt protection for easy loading and unloading
- Complete stainless steel inner chamber without permanent fixtures
- Easy cleaning

## ► OPTIONS

- Access ports of various diameters with silicone plugs
- Racks, chrome-plated or stainless steel
- Perforated shelf, stainless steel
- Independent temperature safety device class 3.1 according to DIN 12880
- Door with window and interior lighting
- Door lock
- Door gasket made of FKM (Viton)
- Switchable audio alarm for overheating
- Analog output for temperature 4 – 20 mA with 6-pin DIN socket
- Temperature measurement according to DIN 12880
- Calibration certificate
- Extension to factory calibration certificate (additional measuring point)
- Data Logger Kits and Logger Software



Data Logger Kits



Reinforced shelves



Door with window and interior lighting




Heating chamber with special extension for loading without opening the door

► **BINDER  
INDIVIDUAL**  
More information  
on page 133

# Technical data for your planning and installation

## FED series

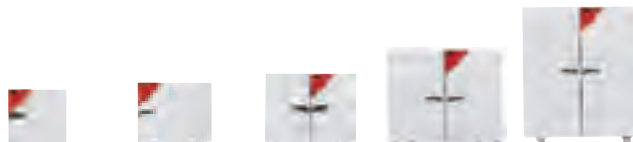

	FED 53	FED 115	FED 240	FED 400	FED 720
<b>Exterior dimensions</b>					
Width (mm/inch)	635 / 25.0	835 / 32.87	1035 / 40.75	1235 / 48.62	1235 / 48.62
Height (including feet/castors) (mm/inch)	620 / 24.41	705 / 27.76	825 / 32.48	1025 / 40.35	1530 / 60.24
Depth (mm/inch)	575 / 22.64	645 / 25.39	745 / 29.33	765 / 30.12	865 / 34.06
plus door handle, I-panel, exhaust duct (mm/inch)	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13
Number of doors	1	1	2	2	2
<b>Interior dimensions</b>					
Width (mm/inch)	400 / 15.75	600 / 23.62	800 / 31.50	1000 / 39.37	1000 / 39.37
Height (mm/inch)	400 / 15.75	480 / 18.90	600 / 23.62	800 / 31.50	1200 / 47.24
Depth (mm/inch)	340 / 13.39	410 / 16.14	510 / 20.08	510 / 20.08	610 / 24.02
Interior volume (l/cu.ft.)	53 / 1.9	115 / 4.1	240 / 8.6	400 / 14.3	720 / 25.7
Number of racks (standard/max.)	2 / 5	2 / 6	2 / 7	2 / 10	2 / 15
Load per rack (kg/lbs.)	15 / 33	20 / 44	30 / 66	35 / 77	45 / 99
Permitted total load (kg/lbs.)	40 / 88	50 / 110	70 / 155	90 / 199	120 / 265
Weight of the unit (empty) (kg/lbs.)	44 / 97	62 / 137	96 / 212	145 / 320	195 / 430
<b>Temperature data</b>					
Temperature range, 5 °C/9 °F above ambient up to °C/°F	300 / 572	300 / 572	300 / 572	300 / 572	300 / 572
Temperature uniformity <sup>1)</sup> at 150 °C/302 °F (±K)	2	1.8	2	2.5	2.0
Temperature fluctuation at 150 °C/302 °F (±K)	0.3	0.3	0.3	0.3	0.3
Heating-up time <sup>2)</sup> to 150 °C/302 °F (min)	24	30	27	35	39
Recov. time after door was opened for 30 sec. <sup>3)</sup> at 150 °C/302 °F (min)	5	8	10	17	20
Air change at 150 °C/302 °F (x/h)	43	32	20	18	12
<b>Electrical data</b>					
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230 1 N ~	230 1 N ~	230 1 N ~	400 3 N ~	400 3 N ~
Nominal power (kW)	1.2	1.6	2.7	3.4	5
Energy consumption at 150 °C/302 °F (Wh/h)	397	544	850	1200	1320
<b>Model no.</b>					
	9010-0210	9010-0212	9010-0214	9010-0216	9010-0218

<sup>1)</sup> value without window // <sup>2)</sup> to 98 % of the set value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %.

The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber.






All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



## ► DIMENSIONS

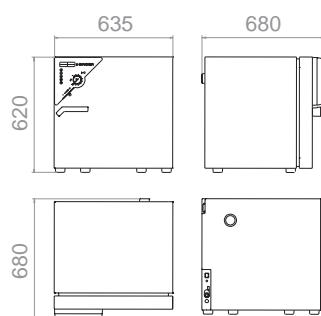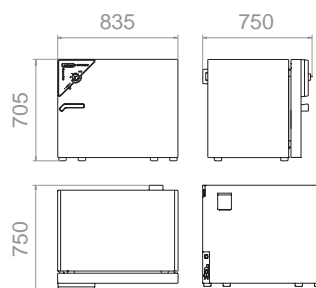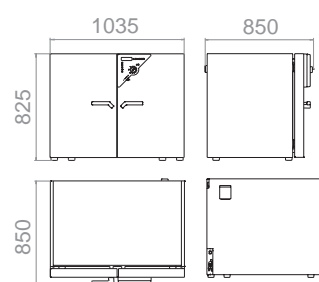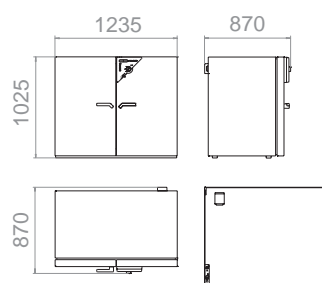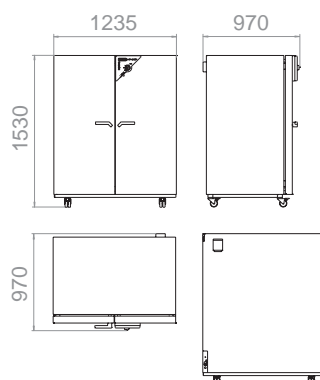**FED 53****FED 115****FED 240****FED 400****FED 720**

## ► INSTALLATION REQUIREMENTS

	FED 53	FED 115	FED 240	FED 400	FED 720
Nominal voltage ( $\pm 10\%$ ) 50/60 Hz (V)	230 1 N ~	230 1 N ~	230 1 N ~	400 3 N ~	400 3 N ~
Nominal power (kW)	1.2	1.6	2.7	3.4	5.0





# Material testing

## FP | M series

### Proven experts: BINDER Material test chambers

- ▶ Proven level of performance
  - ▶ Fan with increased airflow rate
  - ▶ Adjustable rapid air exchange rate
- ▶ Proven precision
  - ▶ High standard according to DIN 12880 (27-point measurement)
  - ▶ The specialists for demanding heating profiles
  - ▶ Uniform test conditions throughout the chamber interior
- ▶ Proven versatility
  - ▶ Communication interface
  - ▶ Digital multi-program controller
  - ▶ Adjustable ventilation, program-controlled (M series)

For demanding heating profiles, these chambers show what they're made of: Best APT.line™ preheating chamber for maximum precision, wide temperature range and comprehensive programming options, with which you can customize ramps, profiles and processes.



Tempering of seals



Drying of electronic components




Plastics testing

# Material test chambers with mechanical convection FP series

The FP series carries out the most demanding tests and scores particularly well with its comprehensive programming options. The mechanical (forced) convection provides reliably short drying and extremely fast heating times – even and especially for chambers under full loads.




► FP 115 model

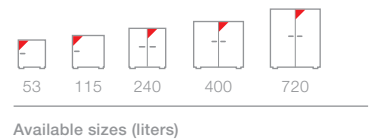

## ► EQUIPMENT

- Temperature range from 5 °C above ambient temperature to 300 °C
- MP controller with 2 programs with 10 sections each, alternatively switchable to program with 20 segments
- The time of an individual program step can be set to max. 999 hours and 59 minutes.
- Adjustable ramp function via program editor
- Integrated weekly program timer with real-time function
- Digital temperature setting accurate to tenths of a degree or to one degree
- Adjustable fan speed
- Adjustable ventilation by means of front ventilation flap slide and rear exhaust Ø 50 mm
- Elapsed time indicator
- Independent adjustable temperature safety device class 2 (DIN 12880), with visual temperature alarm
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Units up to 115 liters are stackable
- Two chrome-plated racks included

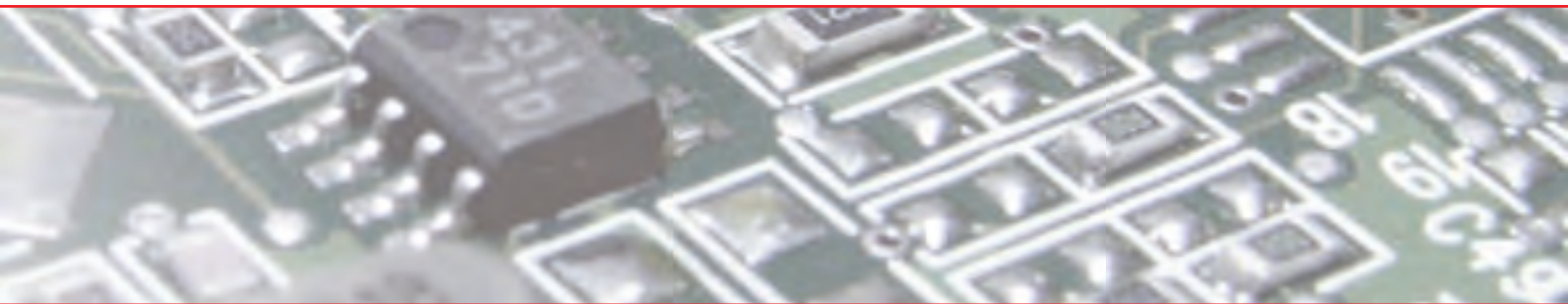


## ► FP SERIES | BEST TEST RESULTS:



### Uniform test conditions

- APT.line™ preheating chamber
  - Homogeneous temperature distribution
  - Identical test conditions throughout the chamber interior independent of sample size and quantity

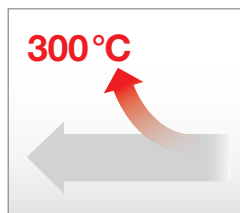

### Broad range of applications

- Rapid air exchange rate and large capacity reserves
- Adjustable fan speed
- Short heating up times



### Best quality and precision guaranteed

- High standard according to DIN 12880 (27-point measurement)
- Short delivery times
- Minimal maintenance and operating costs



### Convenient work environment

- Hermetic door closure with 2-points door closure
- Low heat dissipation due to 60 mm insulation
- Rack with tilt protection for easy loading and unloading
- Complete stainless steel inner chamber
- No permanent fixtures

## ► OPTIONS

- Access ports with silicone plugs
- Racks, chrome-plated or stainless steel
- Perforated shelf, stainless steel
- Reinforced rack, stainless steel
- Reinforced inner chamber with 2 reinforced racks
- Independent temperature safety device class 3.1 according to DIN 12880
- Door with window and interior lighting
- Door lock
- Analog output for temperature 4 - 20 mA with 6-pin DIN socket including DIN connector
- Additional measuring channel for display of specimen temperature (PT 100 sensor)
- Temperature measurement according to DIN 12880
- HEPA fresh-air filter, class EU 14
- Increased air exchange rate through high-performance fan
- Measurement of air exchange rate according to ASTM D5374
- Isolated switching outputs (6-pin DIN socket)
- Calibration certificate
- Extension to calibration certificate
- Data Logger Kits and Logger software



Door with heated viewing window with interior lighting



Access ports with silicone plugs



Calibrations and validations



Test chamber with special racks and modified airflow

► **BINDER  
INDIVIDUAL**  
More information  
on page 133

# Technical data for your planning and installation

## FP series



	FP 53	FP 115	FP 240	FP 400	FP 720
<b>Exterior dimensions</b>					
Width (mm/inch)	635 / 25.0	835 / 32.87	1035 / 40.75	1235 / 48.62	1235 / 48.62
Height (including feet/castors) (mm/inch)	620 / 24.41	705 / 27.76	825 / 32.48	1025 / 40.35	1530 / 60.24
Depth (mm/inch)	575 / 22.64	645 / 25.39	745 / 29.33	765 / 30.12	865 / 34.06
plus door handle, I-panel and exhaust duct (mm/inch)	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13
Number of doors	1	1	2	2	2
<b>Interior dimensions</b>					
Width (mm/inch)	400 / 15.75	600 / 23.62	800 / 31.50	1000 / 39.37	1000 / 39.37
Height (mm/inch)	400 / 15.75	480 / 18.90	600 / 23.62	800 / 31.50	1200 / 47.24
Depth (mm/inch)	340 / 13.39	410 / 16.14	510 / 20.08	510 / 20.08	610 / 24.02
Interior volume (l/cu.ft.)	53 / 1.9	115 / 4.1	240 / 8.6	400 / 14.3	720 / 25.7
Number of racks (standard/max.)	2 / 5	2 / 6	2 / 7	2 / 10	2 / 15
Load per rack (kg/lbs.)	15 / 33	20 / 44	30 / 66	35 / 77	45 / 99
Permitted total load (kg/lbs.)	40 / 88	50 / 110	70 / 155	90 / 199	120 / 265
Weight of the unit (empty) (kg/lbs.)	45 / 99	62 / 137	98 / 216	145 / 320	184 / 406
<b>Temperature data</b>					
Temperature range, 5 °C/9 °F above ambient up to °C/°F	300 / 572	300 / 572	300 / 572	300 / 572	300 / 572
Temperature uniformity <sup>1)</sup> at 150 °C/302 °F (±K)	2.0	1.8	2.0	2.5	2.0
Temperature fluctuation (±K)	0.3	0.3	0.3	0.3	0.3
Heating-up time <sup>2)</sup> to 150 °C/302 °F (min)	24	30	27	35	39
Recov. time after door was opened for 30 sec. <sup>2)</sup> at 150 °C/302 °F (min)	5	8	10	17	20
Air change at 150 °C/302 °F (x/h)	64	32	20	18	12
<b>Electrical data</b>					
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230 1 N ~	230 1 N ~	230 1 N ~	400 3 N ~	400 3 N ~
Nominal power (kW)	1.2	1.6	2.7	3.4	5.0
Energy consumption at 150 °C/302 °F (Wh/h)	300	544	850	1200	1320
<b>Model no.</b>					
	9010-0153	9010-0255	9010-0263	9010-0265	9010-0267

<sup>1)</sup> value without window // <sup>2)</sup> to 98 % of the set value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %.

The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber.


All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



## ► DIMENSIONS

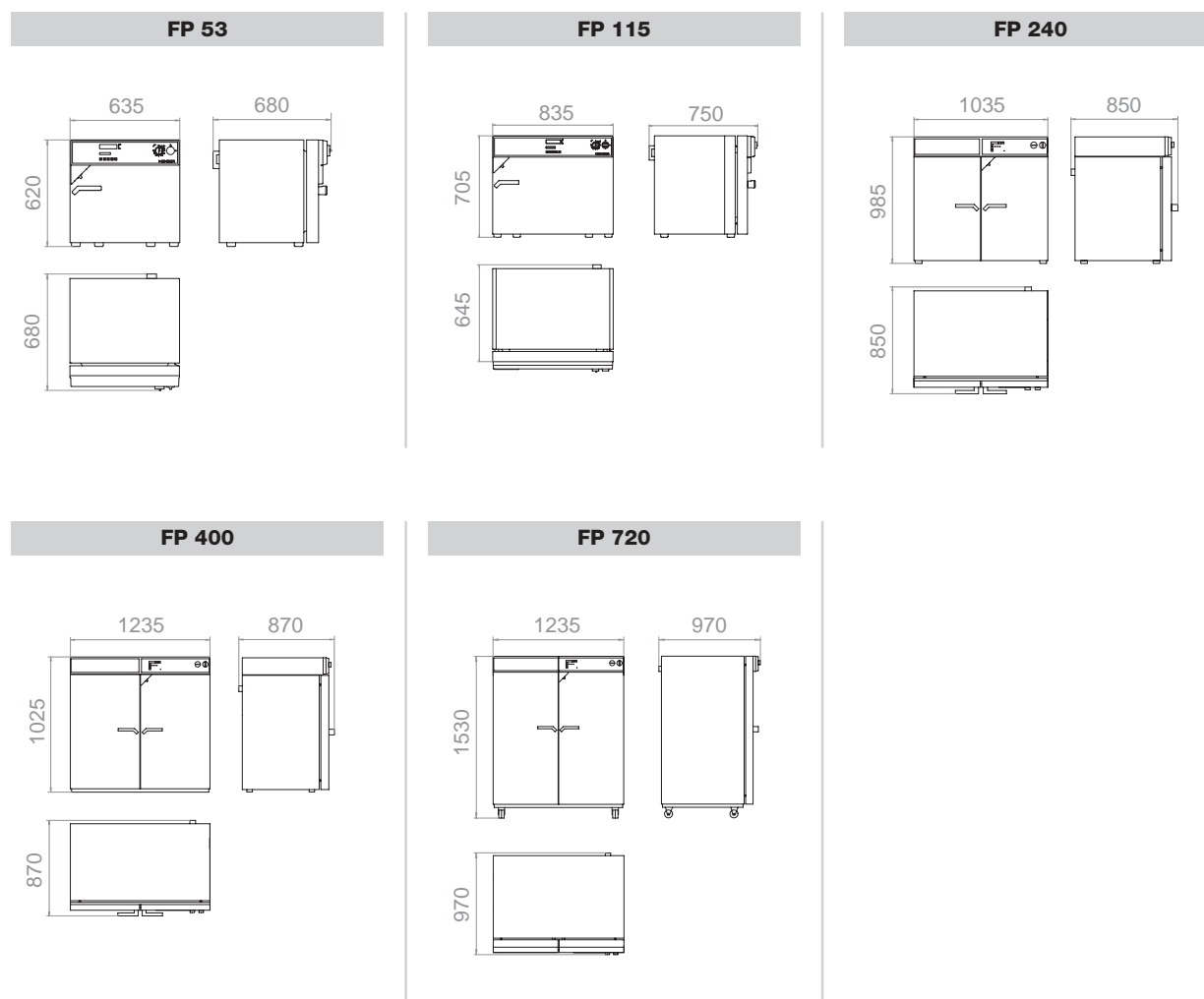

## ► INSTALLATION REQUIREMENTS


	FP 53	FP 115	FP 240	FP 400	FP 720
Nominal voltage ( $\pm 10\%$ 50/60 Hz (V))	230 1N~	230 1N~	230 1N~	400 3N~	400 3N~
Nominal power (kW)	1.2	1.6	2.7	3.4	5.0



# Temperature test chambers with individual programming

## M series

With a maximum temperature of 300 °C and comprehensive programming options, the M series is ideally suited for material and accelerated aging testing. The particularly powerful airflow rate and programmable exhaust air valve ensure fast heating, and test temperatures are maintained at a level with minimal spatial fluctuations more accurately than ever before.

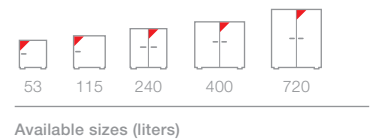

### ► EQUIPMENT

- Temperature range from 5 °C above ambient temperature to 300 °C
- MCS screen controller which can store 25 programs of 100 sections each for a maximum of 500 program segments:
  - User-friendly LCD screen
  - Easy-to-read menu guidance
  - Integrated electronic chart recorder
  - Variety of options for the graphic display of process parameters
  - Real-time clock
- Adjustable ramp function via program editor
- Program-controlled ventilation flap
- High air exchange rate through high-performance fan
- Adjustable fan speed
- Exhaust duct Ø 50 mm
- Independent adjustable temperature safety device class 2 (DIN 12880), with visual temperature alarm
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Units up to 115 liters are stackable
- Two chrome-plated racks included

## ► M SERIES | BEST TEST RESULTS:




### Uniform test conditions

- APT.line™ preheating chamber
  - Homogeneous temperature distribution
  - Same test conditions throughout the chamber interior independent of sample size and quantity



### Best quality and precision guaranteed

- High standard according to DIN 12880 (27-point measurement)
- Short delivery times
- Minimal maintenance and operating costs

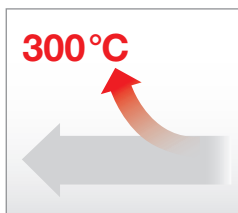


### Broad range of applications

- High air exchange rate and large power reserves
- Adjustable fan speed
- Short heating up times



### Convenient work environment

- Hermetic door closure with 2-points door closure
- Low heat dissipation due to 60 mm insulation
- Rack with tilt protection for easy loading and unloading
- Complete stainless steel inner chamber
- No permanent fixtures

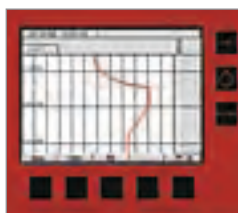

### Custom programming


- Controller with color display which can store 25 programs
- Integrated digital continuous-line recorder for monitoring limits and alarm function
- Convenient documentation
- All displayed values read at a glance

## ► OPTIONS

- Access ports with silicone plugs
- Racks, chrome-plated or stainless steel
- Perforated shelf, stainless steel
- Reinforced rack, stainless steel
- Reinforced inner chamber with 2 reinforced racks
- Door with window and interior lighting
- Door lock
- Analog output for temperature 4 - 20 mA with 6-pin DIN socket
- Additional measuring channel for display of specimen temperature (PT 100 sensor)
- Inert gas connection (gas inlet and outlet)
- Temperature measurement according to DIN 12880
- HEPA fresh-air filter, class EU 14
- Ventilation measurement according to ASTM D5374 with definition and protocol
- Calibration certificate
- Extension to calibration certificate
- Data Logger Kits and Logger software



Reinforced shelves

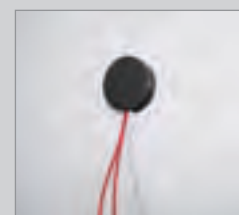

Various access ports



Data Logger Kits



Test chamber with electrical door lock

**BINDER  
INDIVIDUAL**  
More information  
on page 133

# Technical data for your planning and installation

## M series



	M 53	M 115	M 240	M 400	M 720
<b>Exterior dimensions</b>					
Width (mm/inch)	635 / 25.0	835 / 32.87	1035 / 40.75	1235 / 48.62	1235 / 48.62
Height (including feet/castors) (mm/inch)	780 / 30.71	865 / 34.06	985 / 38.78	1185 / 46.65	1695 / 66.73
Depth (mm/inch)	575 / 22.64	645 / 25.39	745 / 29.33	765 / 30.12	865 / 34.06
plus door handle, I-panel and exhaust duct (mm/inch)	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13	105 / 4.13
Quantity of doors	1	1	2	2	2
<b>Interior dimensions</b>					
Width (mm/inch)	400 / 15.75	600 / 23.62	800 / 31.50	1000 / 39.37	1000 / 39.37
Height (mm/inch)	400 / 15.75	480 / 18.90	600 / 23.62	800 / 31.50	1200 / 47.24
Depth (mm/inch)	340 / 13.39	410 / 16.14	510 / 20.08	510 / 20.08	610 / 24.02
Interior volume (mm/inch)	53 / 1.9	115 / 4.1	240 / 8.6	400 / 14.3	720 / 25.7
Quantity of racks (standard/max.)	2 / 5	2 / 6	2 / 7	2 / 10	2 / 15
Load per rack (kg/lbs.)	15 / 33	20 / 44	30 / 66	35 / 77	45 / 99
Permitted total load (kg/lbs.)	40 / 88	50 / 110	70 / 155	90 / 199	120 / 265
Weight of the unit (empty) (kg/lbs.)	61 / 135	89 / 196	131 / 289	173 / 382	203 / 448
<b>Temperature data</b>					
Temperature range, 5 °C/9 °F above ambient up to °C/°F	300 / 572	300 / 572	300 / 572	300 / 572	300 / 572
Temperature uniformity <sup>1)</sup> at 150 °C/302 °F (±K)	1.3	1.5	1.5	1.5	1.9
Temperature fluctuation (±K)	0.3	0.3	0.3	0.3	0.3
Heating-up time <sup>2)</sup> to 150 °C/302 °F (min)	15	16	19	18	21
Recov. time after door was opened for 30 sec. <sup>1)2)</sup> at 150 °C/302 °F (min)	3	3	3	3	3
Air change at 150 °C/302 °F (x/h)	192	96	60	54	36
<b>Electrical data</b>					
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	230 1 N ~	230 1 N ~	230 1 N ~	400 3 N ~	400 3 N ~
Nominal power (kW)	1.2	1.6	2.7	3.4	5.0
Energy consumption at 150 °C/302 °F (Wh/h)	300	544	850	1200	1320
<b>Model no.</b>					
	9010-0201	9010-0202	9010-0203	9010-0204	9010-0205

<sup>1)</sup> without glass door // <sup>2)</sup> to 98 % of the set value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %.

The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber.

All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.








Current information and values are available at:

[www.binder-world.com](http://www.binder-world.com)



## ► DIMENSIONS

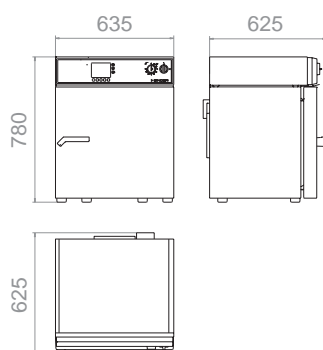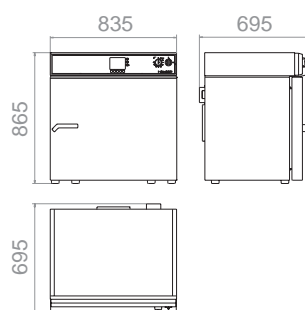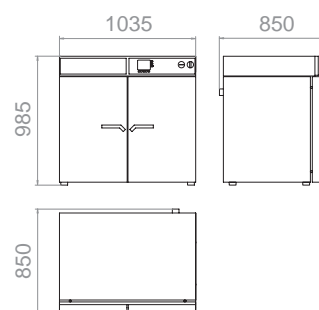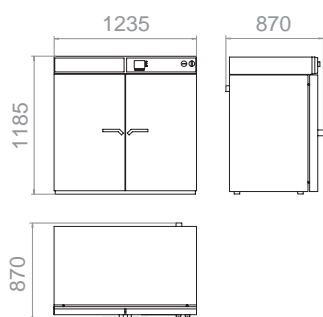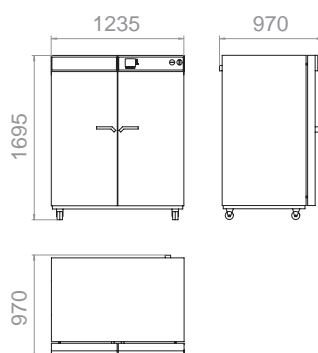**M 53**

**M 115**

**M 240**

**M 400**

**M 720**


## ► INSTALLATION REQUIREMENTS

	FP 53	FP 115	FP 240	FP 400	FP 720
Nominal voltage ( $\pm 10\%$ 50/60 Hz (V))	230 1N~	230 1N~	230 1N~	400 3N~	400 3N~
Nominal power (kW)	1.2	1.6	2.7	3.4	5.0



# Safety drying

## FDL | MDL series

### Unconditionally good: BINDER Safety drying ovens

- ▶ Unconditionally accurate
  - ▶ Fast, uniform drying
  - ▶ Wide temperature range up to 350 °C
  - ▶ Low heat dissipation due to 60 mm insulation
  - ▶ Specimen temperature display with temperature sensor
- ▶ Unconditionally safe
  - ▶ Meets EN 1539: Fresh air monitoring with automatic switchoff
  - ▶ Safety class IP 33
  - ▶ Hermetic door with 2-points door closure
- ▶ Unconditionally versatile
  - ▶ Extension for coil coating test available for MDL series
  - ▶ Digital multi-program controller
  - ▶ Communication interface

For the perfect finish to your solvent-based paints and coating materials, we offer you two series. Absolute temperature accuracy with a high air exchange rate and thus the basis for the best results qualitatively and reproducible tests. You can work safely at all times thanks to replaceable fresh air cartridges and electronic monitoring.



Surface technology / Coating materials



Drying adhesives



# Safety drying ovens

## FDL series

This series provides the perfect environment for all specimens containing solvents: the high efficiency filter cartridge and symmetric airflow provide a silicone free and dust-free inner chamber. The FDL meets all EN 1539 requirements and also ensures maximum work safety with the intelligent fresh air monitoring program.



► FDL 115 model



115

Available size (liters)

### ► EQUIPMENT

- All safety features compliant with EN 1539
- Temperature range from 5 °C above ambient temperature to 300 °C
- MP controller with 2 programs with 10 sections each, or alternatively one program with 20 sections
- The time of an individual program step can be set to max. 999 hours and 59 minutes
- Integrated weekly program timer with real-time function
- Digital temperature setting with an accuracy of one or a tenth of a degree
- Elapsed time indicator
- Door gasket made of FKM (up to max. 200 °C)
- Replaceable fresh-air filter cartridge, class F6 (EU6 fine particle filter for particle sizes between 1 µm and 10 µm)
- Independent adjustable temperature safety device class 2 (DIN 12880), with audible and visual alarm
- Fresh-air monitoring with audible and visual alarm and automatic deactivation of heating
- Rear exhaust connector Ø 100 mm
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Two chrome-plated racks included

## ► FDL SERIES | BEST TEST RESULTS:



### Uniform test conditions

- APT.line™ preheating chamber
  - Homogeneous temperature distribution
  - Identical test conditions throughout the chamber interior independent of sample size and quantity



### Convenient work environment

- Hermetic door closure with 2-points door closure
- Low heat dissipation due to 60 mm insulation
- Rack with tilt protection
- Complete stainless steel inner chamber
- No permanent fixtures
- Silicone free




### Maximum work safety


- Defined solvent quantity meets drying safety standards
- EN 1539 compliant: Fresh air monitoring with automatic switchoff
- Safety class IP 33
- Easy to replace fresh air filter

## ► OPTIONS


- Access ports, Ø 10 mm, Ø 30 mm, right, left
- Racks, chrome-plated or stainless steel
- Perforated shelf, stainless steel
- Specimen temperature display with temperature sensor and analog output 4 – 20 mA
- Replacement air filter (class F6/EU6 for particle sizes between 1 µm and 10 µm)
- Door lock
- Door gasket made of silicone for applications > 200 °C
- Calibration certificate, measurement in center
- Extension to factory calibration certificate (additional measuring point)
- Temperature measurement according to DIN 12880

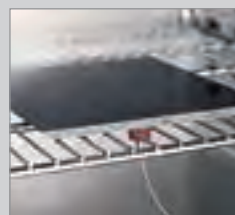

Temperature measurement of the specimen

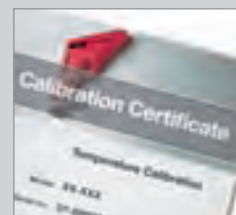

Calibrations and validations

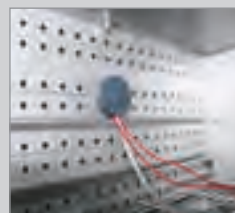

Various access ports



Safety drying oven with UV lamp for testing UV resistance

► **BINDER  
INDIVIDUAL**  
More information  
on page 133

# Technical data for your planning and installation

## FDL series



	FDL 115
<b>Exterior dimensions</b>	
Width (mm/inch)	835 / 32.87
Height (including feet) (mm/inch)	800 / 31.50
Depth (mm/inch)	685 / 26.97
plus door handle (mm/inch)	50 / 1.97
<b>Interior dimensions</b>	
Width (mm/inch)	600 / 23.62
Height (mm/inch)	435 / 17.13
Depth (mm/inch)	435 / 17.13
Interior volume (l/cu.ft.)	115 / 4.1
Quantity of racks (standard/max.)	2 / 5
Load per rack (kg/lbs.)	20 / 44
Permitted total load (kg/lbs.)	50 / 110
Weight of the unit (empty) (kg/lbs.)	90 / 199
<b>Temperature data</b>	
Temperature range, 5 °C/9 °F above ambient up to °C/°F	300 / 572
Temperature uniformity at 150 °C/302 °F (±K)	2.5
Temperature fluctuation (±K)	0.3
Heating-up time <sup>2)</sup> at 150 °C/302 °F (min.)	17
Recov. time after door was opened for 30 sec. <sup>2)</sup> at 150 °C/302 °F (min.)	3
Air change (approx. x/min.)	3
Air circulation (approx. x/min.)	40
Exhaust air volume flow (approx. L/Min. m³/h)	400 (24.0)
Air flow velocity (m/sec)	0.8 – 1.2
Highest permitted solvent quantity (g) (at T-180 °C, M-100 g/mol, U-40 g/m3, K=0,5)	6.65
<b>Electrical data</b>	
Housing protection acc. to EN 60529	IP 33
Nominal voltage (+10 %) 50/60 Hz (V)	230
Nominal power (kW)	2.9
Energy consumption at 150 °C/302 °F (Wh/h)	1098
<b>Model no.</b>	
	9010-0269

<sup>1)</sup> value without window // <sup>2)</sup> to 98 % of the set-point value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.




Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



► DIMENSIONS

**FDL 115**

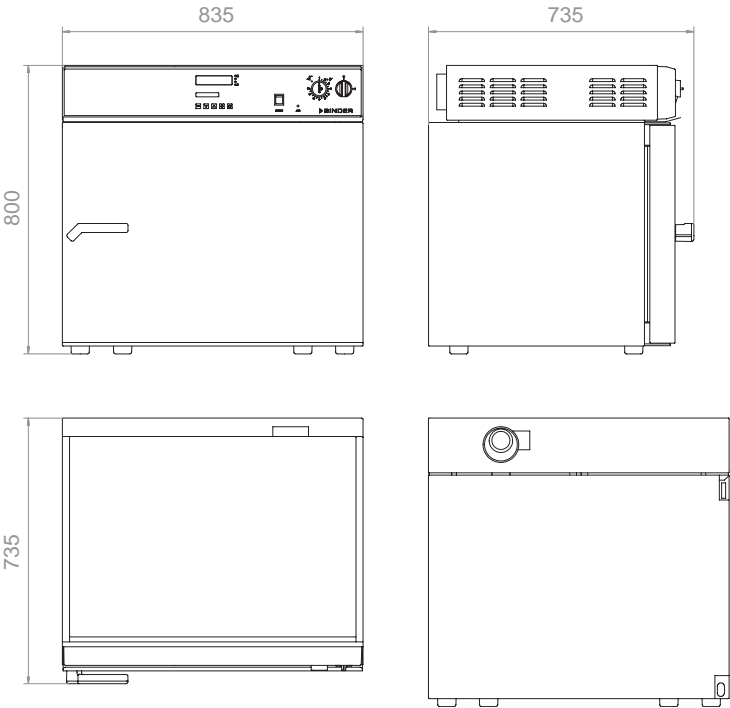

► INSTALLATION REQUIREMENTS

	FDL 115
Nominal voltage (+10 %) 50/60 Hz (V)	230
Nominal power (kW)	2.9

1
2
3
4
5
6

# Safety drying ovens with expanded temperature range

## MDL series

The MDL series works at temperatures up to 350 °C and at an extremely high airflow rate. It is perfectly suited for high-temperature testing such as in the coil coating testing sector. The preheating chamber technology with special airflow design allows evenly distributed burn-in over a very short period – and all that with maximum work safety due to electronic fresh air monitoring. The customized programming also gives you all the freedom you need for your work.



► MDL 115 model



115

Available size (liters)

### ► EQUIPMENT

- Compliant with all EN 1539 safety requirements
- Temperature range from 5 °C above ambient temperature to 350 °C
- MCS controller which can store 25 programs of 100 sections each for a maximum of 500 program segments
  - User-friendly LCD screen
  - Easy-to-read menu guidance
  - Integrated electronic chart recorder
  - Variety of options for the graphic display of process parameters
  - Real-time clock
- Heating output 9.0 kW
- Door gasket made of high temperature resistant silicone
- Rear exhaust connector Ø 100 mm
- Replaceable fresh-air filter cartridge, class F6 (EU6 fine particle filter for particle sizes between 1 µm and 10 µm)
- Independent adjustable temperature safety device class 2 (DIN 12880), with audible and visual alarm
- Fresh-air monitoring with audible and visual alarm and automatic heating feature deactivation
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Two chrome-plated racks included

## ► MDL SERIES | BEST TEST RESULTS:



### Uniform test conditions

- APT.line™ preheating chamber
  - Homogeneous temperature distribution
  - Identical test conditions throughout the chamber interior independent of sample size and quantity



### Convenient work environment

- Hermetic door closure with 2-points door closure
- Low heat dissipation due to 60 mm insulation
- Rack with tilt protection
- Complete stainless steel inner chamber
- No permanent fixtures



### Maximum work safety

- Defined solvent quantity meets drying safety standards
- EN 1539 compliant: Fresh air monitoring with automatic switchoff
- Safety class IP 33
- Easy to replace fresh air filter



### Custom programming

- Controller with color display which can store 25 programs
- Integrated digital continuous-line recorder for monitoring limits and alarm function
- Convenient documentation
- All measured values read at a glance

## ► OPTIONS

- Racks, chrome-plated or stainless steel
- Perforated shelf, stainless steel
- Extension in the door for coil coating tests
- Replacement air filter (class F6/EU6 for particle sizes between 1 µm and 10 µm)
- Additional measuring channel for digital display of specimen temperature with clip temperature sensor
- Door lock
- Lockable controller keyboard
- Calibration certificate, measurement in center
- Extension to factory calibration certificate (additional measuring point)
- Temperature measurement according to DIN 12880



Temperature measurement of the specimen



Extension for coil coating applications



Calibrations and validations



Safety drying oven with HEPA fresh-air filter EU 14

► **BINDER  
INDIVIDUAL**  
More information  
on page 133



# Technical data for your planning and installation

## MDL series



	MDL 115
<b>Exterior dimensions</b>	
Width (mm/inch)	835 / 32.87
Height (including feet) (mm/inch)	800 / 31.50
Depth (mm/inch)	685 / 26.97
plus door handle (mm/inch)	50 / 1.97
<b>Interior dimensions</b>	
Width (mm/inch)	600 / 23.62
Height (mm/inch)	435 / 17.13
Depth (mm/inch)	435 / 17.13
Interior volume (l/cu.ft.)	115 / 4.1
Quantity of racks (standard/max.)	2 / 5
Load per rack (kg/lbs.)	20 / 44
Permitted total load (kg/lbs.)	50 / 110
Weight of the unit (empty) (kg/lbs.)	90 / 199
<b>Temperature data</b>	
Temperature range, 5 °C/9 °F above ambient up to °C/°F	350 / 662
Temperature uniformity at 150 °C/302 °F (±K)	3.4
Temperature uniformity with door flap at 150 °C/302 °F (±K)	3
Temperature fluctuation (±K)	0.5
Heating-up time <sup>1)</sup> at 150 °C/302 °F (minutes)	6
Recov. time after door was opened for 30 sec. <sup>2)</sup> at 150 °C/302 °F (minutes)	2
Recov. time after door was opened for 30 sec. <sup>2)</sup> with door flap, at 150 °C/302 °F (minutes)	1
Air change (approx. x/min.)	3
Air circulation (approx. x/min.)	40
Exhaust air volume flow (approx. L/Min. m³/h)	400 (24.0)
Air flow velocity (m/sec)	0.8 – 1.2
Highest permitted solvent quantity (g) (at T-180 °C, M-100 g/mol, U-40 g/m³, K=0,5)	6.65
<b>Electrical data</b>	
Housing protection acc. to EN 60529	IP 33
Nominal voltage (+10 %) 50/60 Hz (V)	400 V / 3N~
Nominal power (kW)	9
Energy consumption at 150 °C/302 °F (Wh/h)	1130
<b>Model no.</b>	<b>9010-0200</b>

<sup>1)</sup> value without window // <sup>2)</sup> to 98 % of the set-point value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.




Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



► DIMENSIONS

MDL 115

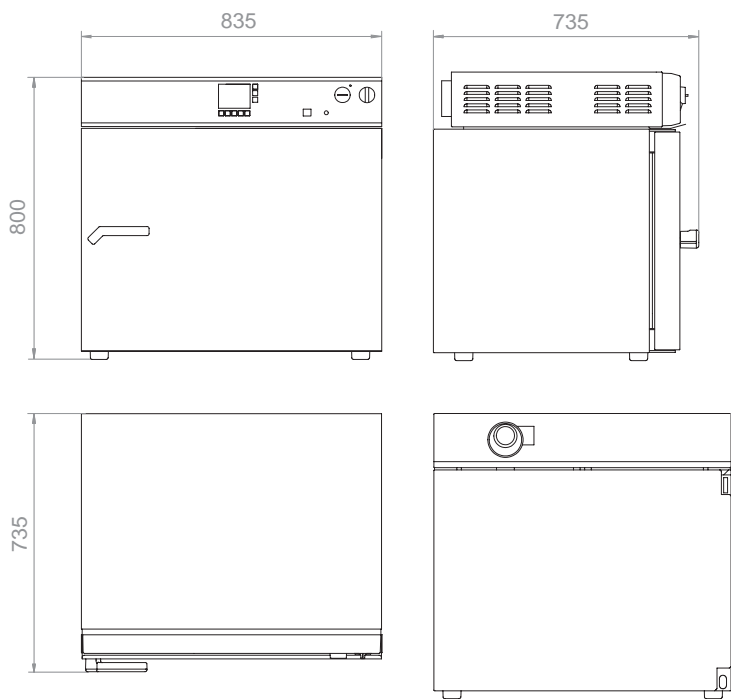

► INSTALLATION REQUIREMENTS

MDL 115	
Nominal voltage (±10 %) 50/60 Hz (V)	400 V / 3N~
Nominal power (kW)	9



# Vacuum drying

## VD | VDL series

### Extremely superior: BINDER Vacuum drying ovens

- ▶ Extremely fast
  - ▶ Direct heat transfer to the sample material through thermal conducting plates
  - ▶ Fast, condensation-free drying processes
  - ▶ Homogeneous sample drying
- ▶ Extremely safe
  - ▶ VDL series with explosion-proof inner chamber (ATEX compliant)
  - ▶ Maximum specimen protection with adjusted overshooting
  - ▶ Ensures optimal work safety through its one-of-a-kind safety concept
- ▶ Extremely durable
  - ▶ Inner chamber made of highly corrosion resistant stainless steel V4A (1.4571)
  - ▶ Coordinated, modular system
  - ▶ Application-specific Vacuum chemical membrane pumps

Both VD and VDL ovens can dry samples completely without residues, scaling or oxidation, and this is achieved in overdrive. Depending on your individual safety requirements, our unique safety concept sets a new standard and is combined with a first-rate performance and quality.



Semiconductors / Electronics



Chemistry / Plastic



Precision engineering

# Vacuum drying ovens for non-flammable solvents VD series

The VD series makes a strong impression with its drying abilities which provide a homogeneous temperature distribution ensured by its electronically controlled APT.line™ preheating chamber. The patented shelf expansion technology guarantees optimal heat transfer. The shelves are easy to clean and provide a flexible positioning.



► VD 53 model



23 53 115


Available sizes (liters)

## ► EQUIPMENT

- Temperature range from 15 °C above ambient temperature to 200 °C
- MP controller with two programs with 10 sections each or switchable to one program with 20 sections
- Integrated weekly program timer with real-time function
- Digital temperature setting with an accuracy of one degree
- Precision-adjustable ventilation valve (for VD 23, the inert gas connection is also used as the ventilation valve)
- Precision-adjustable inert gas valve with Cross-Flow Technology
- All electrical components are decoupled from the inner chamber
- Spring-mounted safety glass panel with shatter protection
- Independent adjustable temperature safety device class 2 (DIN 12880), with visual temperature alarm
- Analog pressure gauge (displays pressure difference between the inner chamber and ambient pressure)
- Electro polished inner chamber, suction and ventilation tubes, pressure container, expansion racks and ball valve are made of highly corrosion resistant stainless steel V4A (1.4571)
- Door gasket made of tempered silicone
- Two x 24 V DC (max 0.4 A) switching outputs, switched via two control contacts in the program editor
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Two patented, flexible aluminum expansion racks
- Also available as complete system with module and vacuum pump



## ► VD SERIES | FAST AND CONDENSATION-FREE DRYING PROCESS:

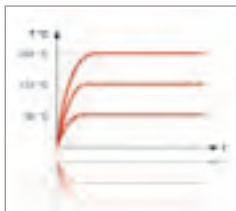

### Process stability for perfect results

- Innovative control concept for regulation of the entire temperature range without overshooting
  - Short heating up times
  - Easy operation
- APT.line™ preheating chamber
  - Gentle drying throughout the chamber interior
  - Direct heat transfer through large thermal conducting plates
  - Patented, flexible positioning of the expansion racks



### Accelerated drying process

- BINDER Cross-Flow Principle
- Even flow throughout the inner chamber from bottom to top
- Finely adjustable inert gas valve without turbulence for lighter samples
- Individually controlled



### Convenient, safe work environment

- BINDER Complete system
- Coordinated, modular system consisting of vacuum drying oven, application-specific vacuum pumps and vacuum module.
  - Connection kit for easy assembly
  - Optimal working height
  - Low noise
  - Pressure and temperature profiles are depicted simultaneously



### Tested security

- BINDER safety concept
- Spring-mounted safety glass panel with shatter protection
- Standard inert gas connection for interior flushing
- Electronic components are decoupled from the inner chamber



### Simple, time-saving cleaning

- Smooth inner chamber with rounded corners
- Inner chamber and all connections made of highly corrosion resistant stainless steel V4A (1.4571)
  - Fixtures are fully removable

## ► OPTIONS

- Expansion racks, stainless steel
- Connection kit with various small-flange components
- Measuring port for vacuum-tight access port of measuring lines into the device (9-pin)
- Temperature measurement of the specimen with flexible PT 100 sensor and digital specimen temperature display
- Digital pressure display, measuring range from 1 mbar to atm. pressure, display accuracy 1 mbar
- Calibration certificates
- Extension to calibration certificate (additional measuring points)
- Door gasket, FKM (Viton)
- Vacuum module with chemical membrane pump VP 1.1
- Vacuum module with chemical membrane pump VP 2.1
- Vacuum module with speed-controlled chemical membrane pump VP 3.1
- Vacuum module for installation of vacuum pumps
- Measuring port
- Specimen temperature display with PT 100 sensor
- Validations and calibrations
- Vacuum drying oven with special racks for large numbers of particularly flat samples



Measuring port



Specimen temperature display with PT 100 sensor



Validations and calibrations



Vacuum drying oven with special racks for large numbers of particularly flat samples

► BINDER  
INDIVIDUAL  
More information  
on page 133

# Technical data for your planning and installation

## VD series



	VD 23	VD 53	VD 115
<b>Exterior dimensions</b>			
Width VD (mm/inch)	515 / 20.28	635 / 25.0	740 / 29.13
Height VD (including feet) (mm/inch)	655 / 25.79	775 / 30.51	900 / 35.43
Total height VD + option "vacuum module" (mm/inch)	1280 / 50.39	1400 / 55.12	1525 / 60.04
Depth incl. door handle, connection (mm/inch)	600 / 23.62	650 / 25.59	770 / 30.31
<b>Interior dimensions</b>			
Width (mm/inch)	285 / 11.22	400 / 15.75	506 / 19.92
Height (mm/inch)	285 / 11.22	400 / 15.75	506 / 19.92
Depth (mm/inch)	295 / 11.61	340 / 13.39	460 / 18.11
Interior volume (l/cu.ft.)	23 / 0.81	53 / 1.87	115 / 4.06
Quantity of expansion racks (aluminum) (standard/max.)	2 / 4	2 / 5	2 / 6
Load per rack (kg/lbs.)	20 / 44	20 / 44	20 / 44
Permitted total load (kg/lbs.)	35 / 77	45 / 99	65 / 143
Weight of the unit (empty) (kg/lbs.)	63 / 139	95 / 210	153 / 338
<b>Temperature data</b>			
Temperature range, 15 °C/27 °F above ambient up to °C/°F	200 / 392	200 / 392	200 / 392
Temperature uniformity <sup>1)</sup> at 100 °C/212 °F (±K)	1.5	2	3.5
Temperature fluctuation <sup>1)</sup> (±K)	0.1	0.1	0.1
Heating up time <sup>1) 2)</sup> to 100 °C (min.) Position POWER	65	80	95
Permitted end vacuum (mbar / inchHg)	0,01 / 0.000295	0,01 / 0.000295	0,01 / 0.000295
Leak rate max. (bar/h / inHg/h)	0,01 / 2.95	0,01 / 2.95	0,01 / 2.95
<b>Electrical data VD</b>			
Housing protection acc. to EN 60529	IP 20	IP 20	P 20
Nominal voltage (±10 %) 50/60 Hz (V)	230	230	230
Nominal power (kW)	0.8	1.2	1.9
Energy consumption at 100 °C (Wh/h)	105	150	250
<b>Model no.</b>	<b>9030-0029</b>	<b>9030-0030</b>	<b>9030-0031</b>
<b>Electrical data VD-UL</b>			
Nominal voltage (±10 %) 60 Hz (V)	115	115	115
Nominal current (A)	10.5	14.8	7.0
Nominal power (kW)	1.2	1.7	0.8
<b>Model no. UL unit</b>	<b>9030-0035</b>	<b>9030-0036</b>	<b>9030-0037</b>




<sup>1)</sup> values measured with aluminum racks // <sup>2)</sup> to 98 % of the set-point value /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



► DIMENSIONS

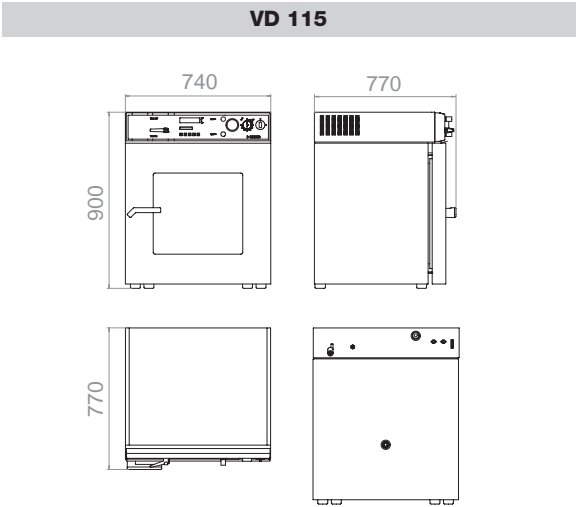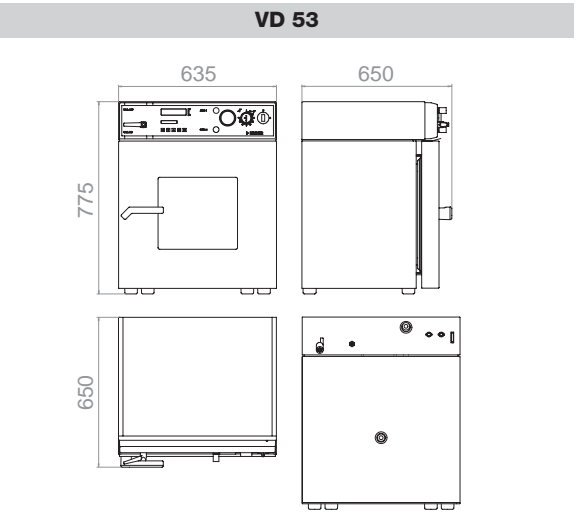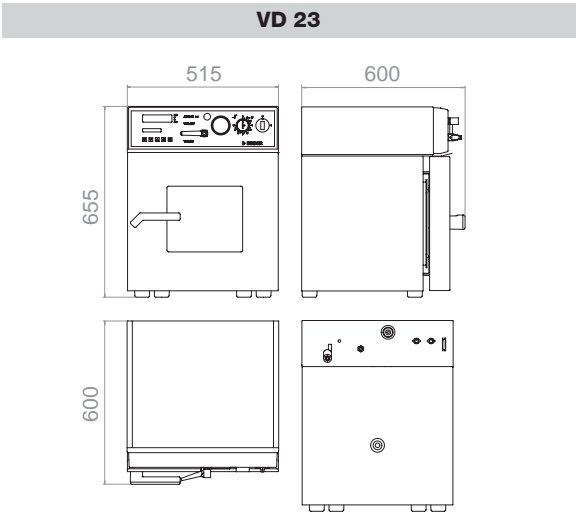

► INSTALLATION REQUIREMENTS

	VD 23	VD 53	VD 115
Vacuum connection with small flange (DN mm/inch)	16 / 0.63	16 / 0.63	16 / 0.63
Measuring access port with small flange (DN mm/inch)	16 / 0.63	16 / 0.63	16 / 0.63
Inert gas connection with flow limiter (RP")	3 / 8	3 / 8	3 / 8

# Safety vacuum drying ovens for flammable solvents

## VDL series

The safety package of the VDL series ensures maximum safety when drying organic solvents standard with TÜV/GS. The inner chamber is designed according to ATEX directive 94/9/EC for Zone EX II 3G.



► VDL 53 model




23 53 115

Available sizes (liters)

### ► EQUIPMENT


- Temperature range from 15 °C above ambient temperature to 200 °C
- MP controller with 2 programs with 10 sections each or switchable to one program with 20 segments
- Digital temperature setting with an accuracy of one degree
- Spring-mounted safety glass panel with shatter protection
- Pressure control device for heating activated < 125 mbar
- Over pressure capsuled instrument panel with compressed air connection and maintenance unit
- Flame protection gasket
- Precision-adjustable ventilation valve
- Precision-adjustable inert gas valve with Cross-Flow Technology
- Analog pressure gauge (displays pressure difference between the inner chamber and ambient pressure)
- Electro polished inner chamber, suction and ventilation tubes, pressure container, expansion racks and ball valve are made of highly corrosion resistant stainless steel V4A (1.4571)
- Door gasket made of tempered silicone
- Independent adjustable temperature safety device class 2 (DIN 12880), with visual temperature alarm
- RS 422 interface for APT-COM™ DataControlSystem communication software
- Two patented, flexible aluminum expansion racks
- Also available as complete system with module and vacuum pump

## ► VDL SERIES | FAST AND CONDENSATION-FREE DRYING PROCESS:

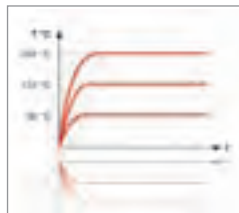

### Process stability for perfect results

- Innovative control concept for regulation of the entire temperature range without overshooting
  - Short heating up times
  - Easy operation
- APT.line™ preheating chamber
  - Gentle drying throughout the chamber interior
  - Direct heat transfer through large thermal conducting plates
  - Patented, flexible positioning of the expansion racks

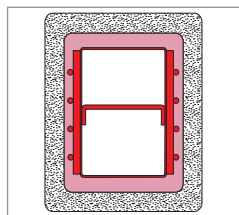

### Accelerated drying process

- BINDER Cross-Flow Principle
- Even flow throughout the inner chamber from bottom to top
- Finely adjustable inert gas valve without turbulence for lighter samples
- Individually controlled



### Standards compliant according to ATEX II 3G

- BINDER safety concept
- Explosion-protected inner chamber in accordance with EX II 3G
- Spring-mounted safety glass panel with shatter protection
- Standard inert glass connection for interior flushing
- Electronic components decoupled from the inner chamber
- Overpressure encapsulated instrument field
- Automatic heating activated < 125 mbar
- Patented flame protection gasket



### Simple, time-saving cleaning

Smooth inner chamber with rounded corners

- Inner chamber and all connections made of highly corrosion resistant stainless steel V4A (1.4571)
- Fixtures are fully removable

## ► OPTIONS

- Expansion racks, stainless steel
- Connection kit with various small-flange components
- Measuring port for vacuum-tight access port of measuring lines into the device (9-pin)
- Temperature measurement of the specimen with flexible PT 100 sensor and digital specimen temperature display
- Calibration certificates
- Extension to calibration certificate (additional measuring points)
- Door gasket, FKM (Viton)
- Vacuum module with ATEX chemical membrane pump VP 4
- Vacuum module with ATEX chemical membrane pump VP 5
- Vacuum module for installation of vacuum pumps



BINDER one-stop solution vacuum module with pump



ATEX chemical membrane pump VP 4



► BINDER  
INDIVIDUAL  
More information  
on page 133

Vacuum drying oven with custom-made front panel for additional protection against gas leakage

# Technical data for your planning and installation

## VDL series



	VDL 23	VDL 53	VDL 115
<b>Exterior dimensions</b>			
Width VDL (mm/inch)	515 / 20.28	635 / 25.0	740 / 29.13
Height VDL (including feet) (mm/inch)	655 / 25.79	775 / 30.51	900 / 35.43
Total height VDL + option "vacuum module" (mm/inch)	1280 / 50.39	1400 / 55.12	1525 / 60.04
Depth incl. door handle, connection (mm/inch)	600 / 23.62	650 / 25.59	770 / 30.31
<b>Interior dimensions</b>			
Width (mm/inch)	285 / 11.22	400 / 15.75	506 / 19.92
Height (mm/inch)	285 / 11.22	400 / 15.75	506 / 19.92
Depth (mm/inch)	295 / 11.61	340 / 13.39	460 / 18.11
Interior volume (l/cu.ft.)	23 / 0.81	53 / 1.87	115 / 4.06
Quantity of expansion racks (aluminum) (standard/max.)	2 / 4	2 / 5	2 / 6
Load per rack (kg/lbs.)	20 / 44	20 / 44	20 / 44
Permitted total load (kg/lbs.)	35 / 77	45 / 99	65 / 143
Weight of the unit (empty) (kg/lbs.)	63 / 139	95 / 210	153 / 338
<b>Temperature data</b>			
Temperature range, 15 °C/27 °F above ambient up to °C/°F	200 / 392	200 / 392	200 / 392
Temperature uniformity <sup>1)</sup> at 100 °C/212 °F (±K)	1.5	2	3.5
Temperature fluctuation <sup>1)</sup> (±K)	0.1	0.1	0.1
Heating up time <sup>1) 2)</sup> to 100 °C (min.) Position POWER	65	80	95
Permitted end vacuum (mbar/inchHg)	0.01 / 0.000295	0.01 / 0.000295	0.01 / 0.000295
Leak rate (mbar/h / inchHg/h)	10 / 0.295	10 / 0.295	10 / 0.295
Compressed air connection for pressure-encapsulation (Ø mm)	8	8	8
<b>Electrical data</b>			
Housing protection acc. to EN 60529	IP 54	IP 54	IP 54
Nominal voltage (±10 %) 50/60 Hz (V)	230	230	230
Nominal power (kW)	0.8	1.2	1.9
Energy consumption at 100 °C (Wh/h)	105	150	250
<b>Model no.</b>			
	9030-0038	9030-0039	9030-0040

<sup>1)</sup> values measured with aluminum racks // <sup>2)</sup> up to 98 % of the set-point value /// All technical data are specified for units with standard equipment at an ambient temperature of +25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance to factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.






Current information and values are available at  
[www.binder-world.com](http://www.binder-world.com)





► DIMENSIONS

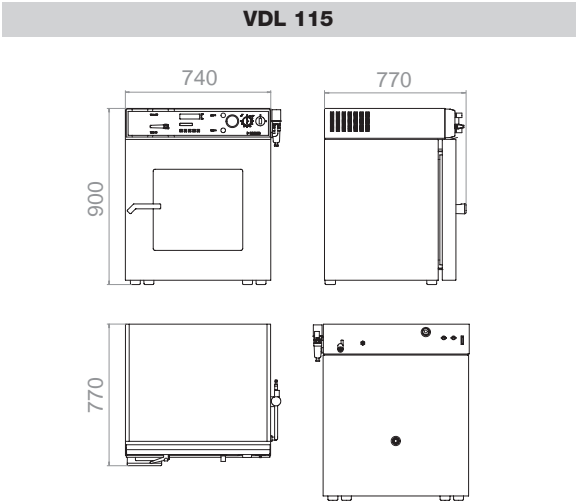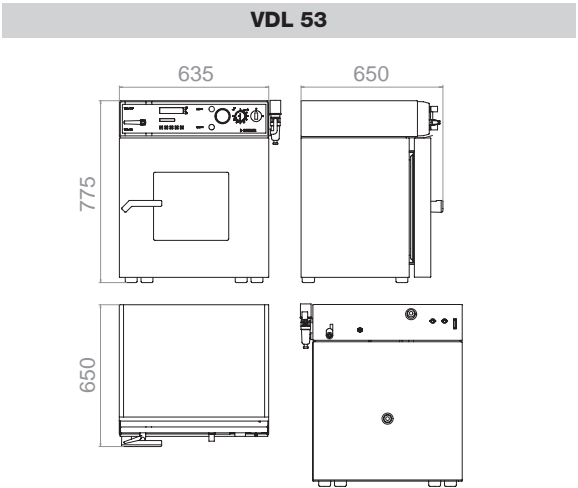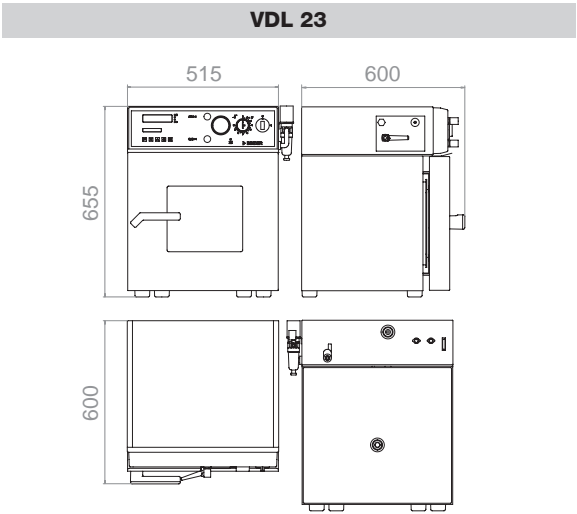

► INSTALLATION REQUIREMENTS

	VDL 23	VDL 53	VDL 115
Vacuum connection with small flange (DN mm/inch)	16 / 0.63	16 / 0.63	16 / 0.63
Measuring access port with small flange (DN mm/inch)	16 / 0.63	16 / 0.63	16 / 0.63
Inert gas connection with flow limiter (RP")	3 / 8	3 / 8	3 / 8

# Vacuum module with pump

## VD series



For the VD series, you get an optional separate vacuum module for an ergonomic working height and to **reduce noise**.

The associated chemical membrane pump is available in three different designs - depending on individual needs. The speed-controlled vacuum pump VP 3.1, for example, adapts to the relevant process cycle and offers up to 30 % time savings.

Additional benefits:

- Excellent pumping speed even at low pressure
- Long lifespan due to its oil-free operation
- Special fluoroplastics provide high chemical resistance



Vacuum module with integrated chemical membrane pump	Vacuum module with chemical membrane pump	Vacuum module with chemical membrane pump	Vacuum module with speed-controlled chemical membrane pump
Type	VP 1.1	VP 2.1	VP 3.1
Nominal airflow	2.0 m³/h	3.4 m³/h	4.6 m³/h
Final pressure	7 mbar	1.5 mbar	1.5 mbar
Electrical connection (50 – 60 Hz)	230 V / 115 V	230 V / 115 V	230 V / 115 V
VD 23 series	•	•	•
VD 53 series	•	•	•
VD 115 series	•	•	•

# Vacuum module with pump

## VDL series



In the vacuum module of the VDL series, there are two different chemical membrane pumps to choose from with different suction capacity and final pressure. Both models are ATEX approved and distinguish themselves by the pressure-capsulated motor with integrated, self-locking overload and overheating protection.

In potentially explosive areas, only ATEX-compliant operating materials may be used since July 1, 2003. All BINDER vacuum pumps for the VDL series are ATEX-compliant.



Vacuum module with integrated chemical membrane pump	Vacuum module with chemical membrane pump, explosion protected	Vacuum module with chemical membrane pump, explosion protected
Typ	VP 4	VP 5
ATEX approval according to ATEX 94/9/EC	yes	yes
Nominal airflow	1.9 m³/h	3.7 m³/h
Final pressure	12 mbar	3 mbar
Electrical connection (50 Hz)	230 V	230 V
VDL 23 series	•	–
VDL 53 series	•	–
VDL 115 series	•	•
Application profile	Approved ATEX-compliant Membrane pump for VDL 23 VDL 53 series	Approved ATEX-compliant Membrane pump with low final vacuum for VDL 115 series

• Option

– not available

We reserve the right to alter technical specifications at any time.







# Environmental simulation

## | Always the right climate

Research and quality assurance is solely focused on the optimization and monitoring of the performance standards, of functions and service life of materials, products and systems. It is expected from a climate chamber that it simulates the effects of temperature, humidity and light on the chemical, physical and biological behavior of the test material in a reproducible manner.

BINDER offers a range of 10 series of constant climate chambers specifically designed for these demanding requirements. They provide innovative technical solutions and comprehensive equipment options, with which specimens can be safely and reliably tested, whether actively or passively. For special technical requirements, users can also rely on the individually tailored solutions from BINDER INDIVIDUAL.

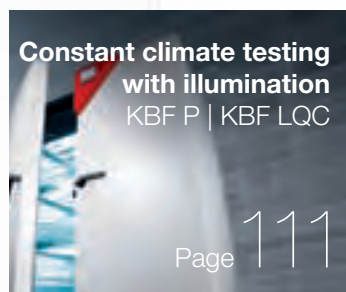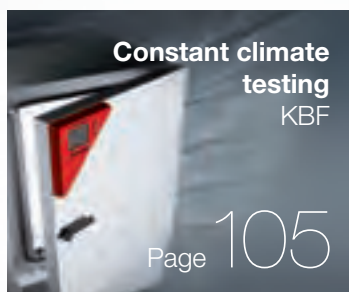





# Constant climate testing

## KBF series

### Proven superior: Constant climate chambers

- ▶ Proven precision
  - ▶ Double-sided horizontal air flow for maximum parameter uniformity
  - ▶ Automatic water and wastewater management
  - ▶ Responsive vapor pressure humidification
  - ▶ Powerful cooling system for safe operation up to 32 °C ambient temperature
  - ▶ Stability tests according to ICH guideline Q1A(R2)
- ▶ Proven documentation
  - ▶ Documentation in support of GMP/GLP compliance
  - ▶ FDA-compliant software APT-COM™
  - ▶ IQ, OQ, PQ validation documentation

The long-term constant temperature and humidity values of BINDER constant climate chambers are impressive. The KBF series is ideal for norm-compliant work according to ICH guidelines. As an inexpensive alternative to using multiple types of climate chambers, the BINDER constant climate chambers can be used for a number of applications.



Basic research



Pharmaceutical industry



Cosmetic industry

# Constant climate chambers

## KBF Series

The KBF is the specialist for reliable stability testing and precise maintenance of constant climatic conditions. From programming to documentation, it meets all important guidelines such as ICH, FDA, GMP and GLP. With its high capacity reserves, automatic water and wastewater management and many options, the BINDER KBF is well prepared for the future.




► KBF 240 model



Available sizes (liters)

### ► EQUIPMENT

- Temperature range from 0 °C to 70 °C (without humidity)
- Humidity range: 10 % to 80 % r. H.
- MCS controller which can store 25 programs of 100 sections each for a maximum of 500 program segments
- User-friendly LCD color screen
- Easy-to-read menu guidance
- Integrated electronic chart recorder
- Variety of options for the graphic display of process parameters
- Real-time clock
- Electronically controlled humidification and dehumidification system with capacitive humidity sensor
- Suitable for stability tests according to ICH guideline Q1A(R2)
- Inner glass door with seal
- Independent temperature safety device class 3.1 (DIN 12880) with visual and audible alarm
- Access port with silicone plug Ø 30 mm, left side
- Complete safety connection kit for water supply and drainage, up to 1 m high
- Ethernet interface for communication software APT-COM™ DataControlSystem
- Two shelves, stainless steel

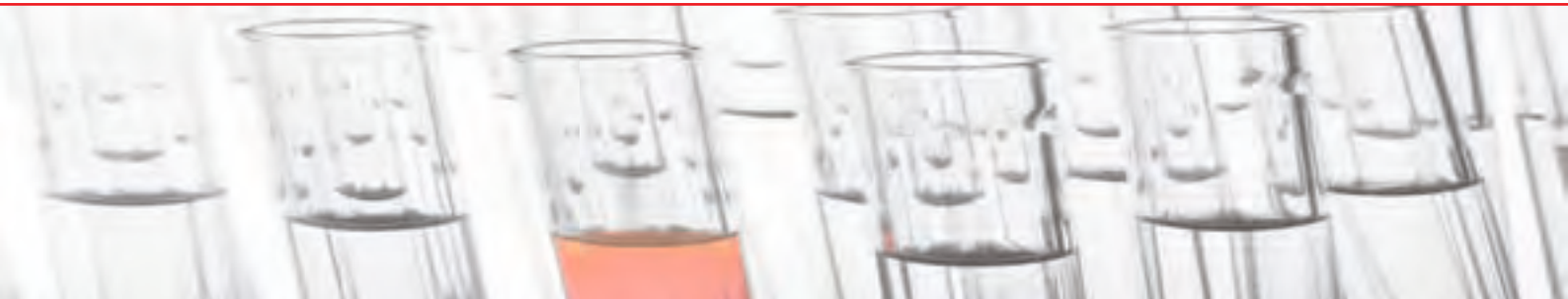

## ► KBF SERIES | BEST TEST RESULTS:



### Precise climate conditions

- APT.line™ preheating chamber
  - Constant and gentle circulation of air through large-surface side walls even under a full load
  - Homogeneous climate conditions throughout test specimens



### Flexible water management

- Sewage pump for discharges up to 1 m high
- Solutions independent of installation sight
- Water treatment with BINDER PURE AQUA SERVICE
- External water supply




### Fast and precise humidification

- Vapor pressure humidification with fast response times
- Drift-free, capacitive humidity sensor
- Short recovery times after door opening
- Finely adjustable humidity control



### Comprehensive additional services

- Data Logger Kits
- Years of experience with recognized validation materials and documents
- Customer-specific modifications

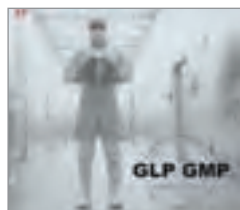

### Norm-compliant work

- Meets all ICH-relevant requirements in accordance with ICH Q1A(R2)
- Documentation and software FDA compliant

## ► OPTIONS

- Access ports of various diameters with silicone plug
- Shelves, stainless steel
- Perforated shelf, stainless steel
- Reinforced rack, stainless steel
- Lockable controller keyboard
- Interior lighting
- Waterproof interior socket 230 V AC
- Additional PT 100 temperature sensor
- RS 422 interface
- External water supply set
- BINDER PURE AQUA SERVICE
- Calibration certificate and extension to calibration certificate
- Measurement of temperature accuracy according to DIN 12880
- Data Logger Kit and software
- Independent temperature safety device class 3.3 (DIN 12880)
- Analog outputs 4 – 20 mA for temperature and humidity measurements
- Isolated alarm outputs for temperature and humidity
- Door lock



Data Logger Kits



BINDER PURE AQUA SERVICE water treatment system



External water supply



Climate chamber with full-view glass door and manual access ports

**BINDER  
INDIVIDUAL**  
More information  
on page 133


# Technical data for your planning and installation

## KBF series



	KBF 115	KBF 240	KBF 720
<b>Exterior dimensions</b>			
Width (mm/inch)	885 / 34.84	930 / 36.61	1255 / 49.41
Height (including feet/casters) (mm/inch)	1050 / 41.34	1460 / 57.48	1925 / 75.79
Depth (mm/inch)	650 / 25.59	800 / 31.50	890 / 35.04
including door handle, I-triangle, connection (mm/inch)	730 / 28.74	880 / 34.65	970 / 38.19
Quantity of doors / Quantity of inner glass doors	1 / 1	1 / 1	2 / 2
<b>Interior dimensions</b>			
Width (mm/inch)	600 / 23.62	650 / 25.59	973 / 38.31
Height (mm/inch)	483 / 19.02	785 / 30.91	1250 / 49.21
Depth (mm/inch)	351 / 13.82	485 / 19.09	576 / 22.68
Interior volume (l/cu.ft.)	102 / 3.60	247 / 8.73	700 / 24.73
Quantity of racks (standard/max.)	2 / 5	2 / 9	2 / 15
Load per rack (kg/lbs.)	30 / 66	30 / 66	45 / 99
Permitted total load (kg/lbs.)	100 / 221	100 / 221	150 / 331
Weight (empty) (kg/lbs.)	129 / 285	184 / 406	309 / 682
<b>Temperature data</b>			
Temperature range <sup>1)</sup> (°C/°F)	0 – 70 / 32 – 158	0 – 70 / 32 – 158	0 – 70 / 32 – 158
Temperature uniformity at 25 °C / 77 °F (±K)	0.2	0.2	0.2
Temperature fluctuation (±K)	0.1	0.1	0.1
Max. heat compensation up to 40 °C / 104 °F (W)	200	300	600
<b>Climatic data</b>			
Temperature range <sup>1)</sup> (°C/°F)	10 – 70 / 50 – 158	10 – 70 / 50 – 158	10 – 70 / 50 – 158
Temperature uniformity at 25 °C / 77 °F and 60 % r. H. (±K)	0.2	0.3	0.2
Temperature fluctuation at 25 °C / 77 °F and 60 % r. H. (±K)	0.1	0.1	0.1
Humidity range (% r. H.)	10 – 80	10 – 80	10 – 80
Humidity fluctuation at 25 °C / 77 °F and 60 % r. H. (±% r. H.)	≤ 2	1.5	1.5
Recovery time after doors were open for 30 sec <sup>2)</sup> at 25 °C / 77 °F and 60 % r. H.	15	4	2
<b>Electrical data</b>			
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~	200 – 240 1N~
Nominal power at 240 V (kW)	2.0	2.1	3.1
Energy consumption <sup>3)</sup> at 40 °C / 104 °F / 75 % r. H. (Wh/h)	470	650	620
Model no.	9020-0139	9020-0150	9020-0108

### ► TEMPERATURE-HUMIDITY CHART

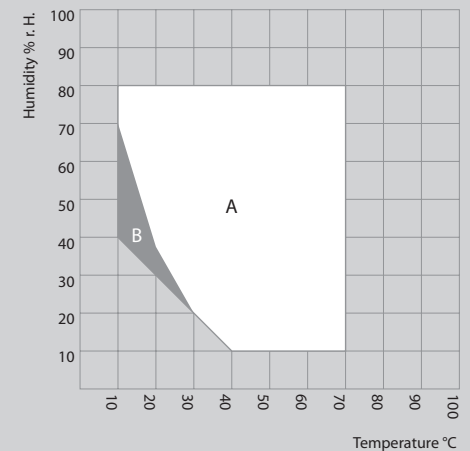

A: Guaranteed condensation-free range  
B: Time-limited operation (max. 24 hours)




<sup>1)</sup> Lower values are valid up to an ambient temperature of max. 25 °C / 77 °F // <sup>2)</sup> to 98 % of the set value // <sup>3)</sup> Use this value for dimensioning air conditioning systems. /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



## ► DIMENSIONS

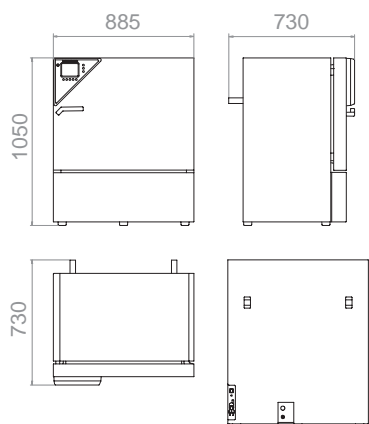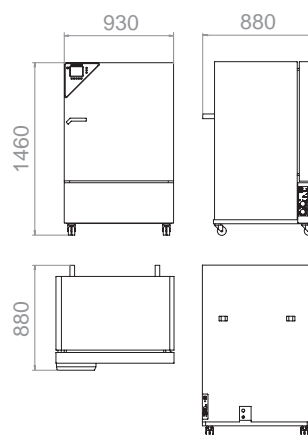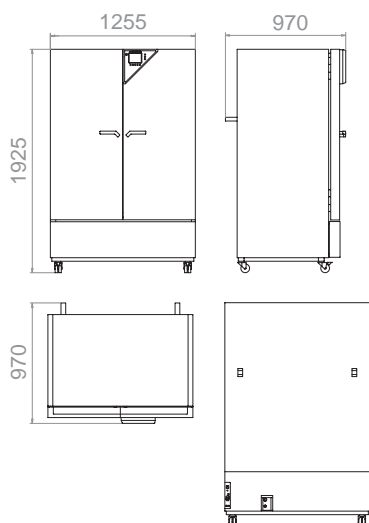**KBF 115**

**KBF 240**

**KBF 720**


## ► INSTALLATION REQUIREMENTS

	KBF 115	KBF 240	KBF 720
Nominal voltage ( $\pm 10\%$ ) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~	200 – 240 1N~
Nominal power at 240 V (kW)	2.0	2.1	3.1





# Constant climate testing with illumination

## KBF P | KBF LQC series

### Consistently better: Constant climate chambers with illumination

- ▶ Consistently better with illumination
  - ▶ Compliance with ICH guideline Q1B, Option 2.
  - ▶ Height-adjustable illumination cassettes
  - ▶ Homogeneous light distribution with the patented Synergy Light™
- ▶ Consistently better with Light Quantum Control
  - ▶ Spherical ball-type sensors for omnidirectional measurement of light intensity
  - ▶ Accurate light valuation directly at the sample location
  - ▶ Automatic shutdown of the light source upon reaching the programmed light dosage

BINDER's constant climate chambers with ICH-compliant illumination provide impressive long-term constant climate and illumination values. With the patented Light Quantum Control (LQC), they offer complete reliable solution for photostability testing in one chamber.



Pharmaceutical industry




Food / Beverage

# Constant climate chambers with ICH-compliant illumination KBF P series

Our expert for photostability testing. In terms of standards compliance, performance and functionality, the KBF P is a true KBF and is also equipped with ICH-compliant illumination. This provides clear and meaningful test results according to ICH guideline Q1B, Option 2. The variable position illumination cassettes with the unique BINDER Q1B Synergy Light™ ensure homogeneous lighting conditions over the entire usable area.



► KBF P 720 model

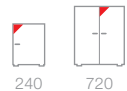

Available sizes (liters)

## ► EQUIPMENT

- Temperature range 0 °C to 70 °C (without illumination and humidity)
- Temperature range 10 °C/20 °C with 60 °C (with illumination and humidity)
- Humidity range 10 % to 75% r. H. (with illumination)
- Humidity range 10 % to 80% r. H. (without illumination)
- Variable position illumination cassettes with BINDER Q1B Synergy Light™
- MCS controller which can store 25 programs of 100 sections each for a maximum of 500 program segments
- User-friendly LCD color screen
- Easy-to-read menu guide
- Integrated electronic chart recorder
- Variety of options for the graphic display of process parameters
- Real-time clock
- Electronically controlled humidification and dehumidification system with capacitive humidity sensor
- Suitable for stability tests according to ICH guideline Q1A(R2) and Q1B Option 2
- Inner glass door with seal
- Independent temperature safety device class 3.1 (DIN 12880) with visual and audible alarm
- Access port with silicone plug Ø 30 mm, left side
- Complete safety connection kit for water supply and drainage, up to 1 m high
- Ethernet interface for communication software APT-COM™ DataControlSystem
- Shelves, stainless steel

## ► KBF P SERIES | BEST TEST RESULTS:



### Homogeneous light distribution

- Unique lighting concept
  - ICH-Q1 B Option 2 compliant light spectrum with BINDER Synergy Light™
  - Variable position and removable illumination cassettes
  - Homogeneous light over the entire usable space



### Precise climate conditions

- BINDER APT.line™ preheating chamber
  - Constant and gentle circulation of air through large-surface side walls even under a full load
  - Homogeneous climate conditions throughout test specimens



### Fast and precise humidification

- Vapor pressure humidification with fast response times
- Drift-free, capacitive humidity sensor
- Short recovery times after door opening
- Finely adjustable humidity control



### Flexible water management

- Sewage pump for discharges up to 1m in height
- Solutions independent of installation sight
- Water treatment with BINDER PURE AQUA SERVICE
- External water supply



### Norm-compliant work

- Meets all ICH-relevant requirements in accordance with ICH Q1A(R2), ICH Q1B, Option 2
- Documentation and software ensure FDA and GMP/GLP compliance

## ► OPTIONS

- Access ports of various diameters with silicone plug
- Shelves, stainless steel
- Perforated shelf, stainless steel
- Reinforced rack, stainless steel
- Lockable controller keyboard
- Additional PT 100 temperature sensor
- RS 422 interface
- Replacement fluorescent tube set for one illumination cassette
- External water supply set
- BINDER PURE AQUA SERVICE
- Calibration certificate and extension to calibration certificate
- Measurement of temperature accuracy according to DIN 12880
- Data Logger Kit and software
- Independent temperature safety device class 3.3 (DIN 12880)
- Analog outputs 4 – 20 mA for temperature and humidity measurements
- Isolated alarm outputs for temperature and humidity
- Door lock



APT-COM™ DataControlSystem  
GLP Edition



Data Logger Kits



BINDER PURE AQUA SERVICE  
water treatment system



Climate chamber with dimmable  
illumination cassettes

► BINDER  
INDIVIDUAL  
More information  
on page 133


# Technical data for your planning and installation

## KBF P series




	KBF P 240	KBF P 720
<b>Exterior dimensions</b>		
Width	930 / 36.61	1255 / 49.41
Height (including feet/casters) (mm/inch)	1460 / 57.48	1925 / 75.79
Depth (mm/inch) including door handle, I-triangle, connection (mm/inch)	880 / 34.65	970 / 38.19
<b>Interior dimensions</b>		
Width / Height / Depth (mm)	650 / 785 / 485	973 / 1250 / 576
Width / Height / Depth (inch)	25.59 / 30.91 / 19.09	38.31 / 49.21 / 22.68
Interior volume (l/cu.ft.)	247 / 8.73	700 / 24.73
Number of light cassettes	2	3
Number of racks (standard/max.)	2 / 9	3 / 12
Weight (empty) (kg/lbs.)	213 / 470	374 / 826
<b>Temperature data</b>		
Temperature range <sup>1)</sup> (°C/°F)	0 – 70 / 32 – 158	0 – 70 / 32 – 158
Temperature range without light cassettes <sup>1)</sup> (°C/°F)	0 – 70 / 32 – 158	0 – 70 / 32 – 158
Temperature range with light cassettes, with illumination <sup>1)</sup> (°C/°F)	10 – 60 / 50 – 140	10 – 60 / 50 – 140
Temperature uniformity at 25 °C / 77 °F (±K)	0.2	0.2
Temperature fluctuation (±K)	0.1	0.1
Max. heat compensation up to 40 °C / 104 °F (W)	400	1000
<b>Climatic data</b>		
Temperature range with light cassettes, with illumination <sup>1)</sup> (°C/°F)	10 – 60 / 50 – 140	20 – 60 / 68 – 140
Temperature uniformity with illumination at 25 °C / 77 °F and 60 % r. H. (±K)	0.6	1.2
Temperature fluctuation with illumination at 25 °C / 77 °F and 60 % r. H. (±K)	0.2	0.2
Humidity range with light cassettes, with illumination <sup>1)</sup> (% r. H.)	10 – 75	10 – 75
Humidity fluctuation at 25 °C / 77 °F and 60 % r. H. with illumination (±% r. H.)	1.5	2
Recovery time after doors were open for 30 sec <sup>2)</sup> at 25 °C / 77 °F and 60 % r. H. with illumination (min)	3	1
<b>Illumination data per light cassette</b>		
ICH compliant illumination for photostability testing <sup>3)</sup> (Lux / UVA W/m <sup>2</sup> )	7500 / 1.4	9900 / 2.1
<b>Electrical data</b>		
Housing protection acc. to EN 60529	IP 20	IP 20
Energy consumption <sup>4)</sup> at 40 °C / 104 °F / 75 % r. H. (Wh/h)	600	2350
<b>Model no.</b>	<b>9020-0160</b>	<b>9020-0143</b>

► **TEMPERATURE-HUMIDITY CHART WITH ILLUMINATION KBF P 240**

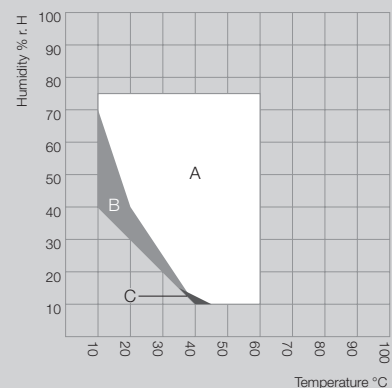

► **TEMPERATURE-HUMIDITY CHART WITH ILLUMINATION KBF P 720**

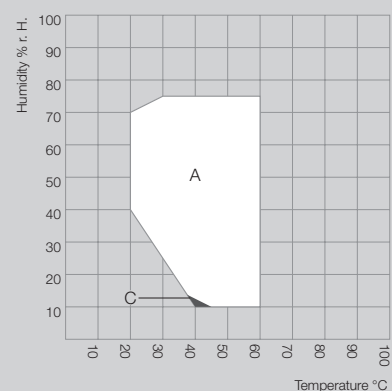

A: Guaranteed condensation-free range  
B: Time-limited operation (max. 24 hours)  
C: Deviations of technical data may be possible



<sup>1)</sup> Lower values are valid up to an ambient temperature of max. 25 °C / 77 °F. // <sup>2)</sup> to 98 % of the set value // <sup>3)</sup> Average value, measured with a spherical sensor (±10 %) by 12 cm / 4.7 in below the light cassette // <sup>4)</sup> Use this value for dimensioning air conditioning systems. /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)



► DIMENSIONS

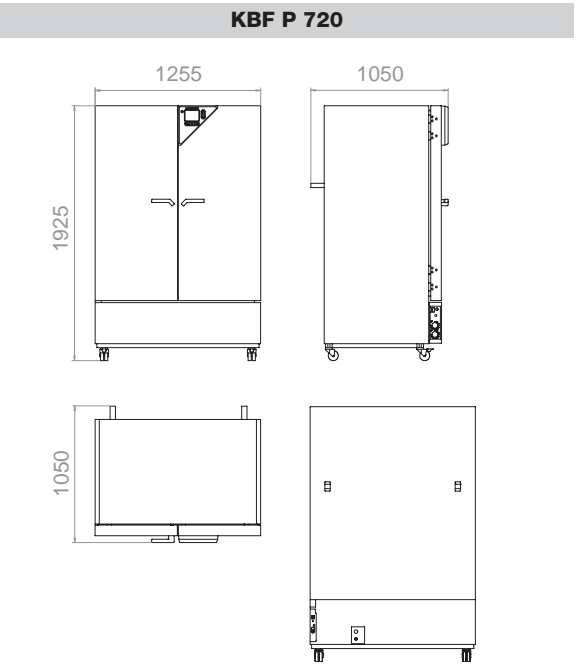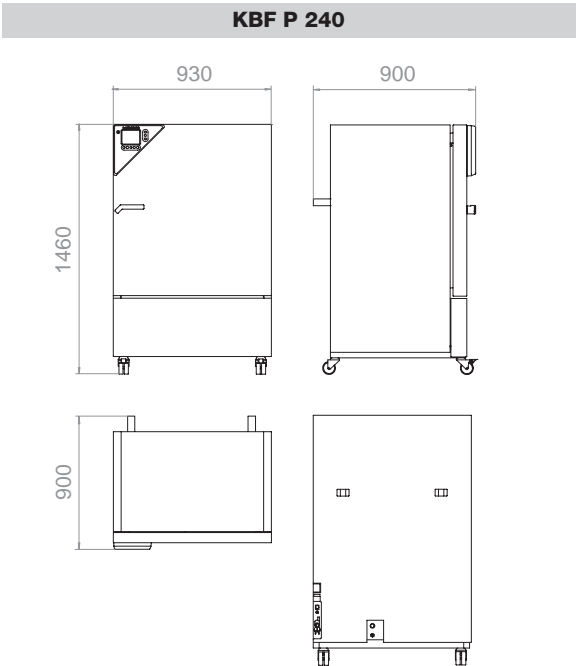

► INSTALLATION REQUIREMENTS

	KB F P 240	KB F P 720
Wall clearance rear (mm/inch)	100 / 3.94	100 / 3.94
Wall clearance side (mm/inch)	100 / 3.94	200 / 7.87
Nominal voltage (±10 %) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~
Nominal power at 240 V (kW)	2.4	3.5


# Constant climate chambers with Light Quantum Control

## KBF LQC series

The special advantage of the KBF LQC series is the photometry, one of our patented enhancements for photostability testing. Two spherical 3D ball-type sensors capture the visible light intensity and UV light intensity at the sample more realistically than any other systems. If the desired light dosage is achieved, the light sources switch off automatically. This convenient solution provides absolute standards compliance.



► KBF LQC 240 model

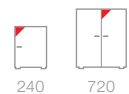

240

720

Available sizes (liters)

### ► EQUIPMENT

- Temperature range 0 °C to 70 °C (without illumination and humidity)
- Temperature range 10 °C/20 °C with 60 °C (with illumination and humidity)
- Humidity range 10 % to 75 % r. H. (with illumination)
- Humidity range 10 % to 80 % r. H. (without illumination)
- Light Quantum Control (LQC), realistic measurement and integration of the light intensity and the UV light intensity at the sample location
- Variable position illumination cassettes with BINDER Q1B Synergy Light™
- MCS controller which can store 25 programs of 100 sections each for a maximum of 500 program segments (User-friendly LCD color screen, easy-to-read menu guide, integrated electronic chart recorder, variety of options for the graphic display of process parameters, real-time clock)
- Electronically controlled humidification and dehumidification system with capacitive humidity sensor Suitable for stability tests according to ICH guideline Q1A(R2) and Q1B Option 2
- Inner glass door with seal
- Independent temperature safety device class 3.1 (DIN 12880) with visual and audible alarm
- Access port with silicone plug Ø 30 mm, left side
- Complete safety connection kit for water supply and drainage, up to 1 m high
- Ethernet interface for communication software APT-COM™ DataControlSystem
- Shelves, stainless steel



## ► KBF LQC SERIES | BEST TEST RESULTS:



### Precise climate conditions

- BINDER APT.line™ preheating chamber
- Constant and gentle circulation of air through large-surface side walls even under a full load
- Homogeneous climate conditions throughout test specimens



### Precise photometry

- LQC - Light Quantum Control
- 3D ball-type sensors capture light true light intensity, can be flexibly placed on the sample material
- Automatically switched off after required light intensity



### Homogeneous light distribution

- Unique lighting concept
- ICH-Q1 B Option 2 compliant light spectrum with BINDER Synergy Light™
- Variable position and removable illumination cassettes
- Homogeneous light over the entire usable space



### Flexible water management

- Sewage pump for discharges up to 1m in height
- Solutions independent of installation sight
- Water treatment with BINDER PURE AQUA SERVICE
- External water supply



### Fast and precise humidification

- Vapor pressure humidification with fast response times
- Drift-free, capacitive humidity sensor
- Short recovery times after door opening
- Finely adjustable humidity control



### Norm-compliant work

- Meets all ICH-relevant requirements in accordance with ICH Q1A(R2), ICH Q1B, Option 2
- Documentation and software ensure FDA and GMP/GLP compliance

## ► OPTIONS

- Access ports of various diameters with silicone plug
- Shelves, stainless steel
- Perforated shelf, stainless steel
- Reinforced rack, stainless steel
- Lockable controller keyboard
- Additional PT 100 temperature sensor
- RS 422 interface
- Replacement fluorescent tube set for one illumination cassette
- External water supply set
- BINDER PURE AQUA SERVICE
- Calibration certificate and extension to calibration certificate
- Measurement of temperature accuracy according to DIN 12880
- Data Logger Kit and software
- Independent temperature safety device class 3.3 (DIN 12880)
- Analog outputs 4 – 20 mA for temperature and humidity measurements
- Isolated alarm outputs for temperature and humidity
- Door lock



Data Logger Kits



Validations and calibrations



External water supply set



Climate chamber with various access ports

► BINDER  
INDIVIDUAL  
More information  
on page 133


# Technical data for your planning and installation

## KBF LQC series




	KBF LQC 240	KBF LQC 720
<b>Exterior dimensions</b>		
Width	930 / 36.61	1255 / 49.41
Height (including feet/casters) (mm/inch)	1460 / 57.48	1925 / 75.79
Depth (mm/inch) including door handle, I-triangle, connection (mm/inch)	880 / 34.65	970 / 38.19
<b>Interior dimensions</b>		
Width / Height / Depth (mm)	650 / 785 / 485	973 / 1250 / 576
Width / Height / Depth (inch)	25.59 / 30.91 / 19.09	38.31 / 49.21 / 22.68
Interior volume (l/cu.ft.)	247 / 8.73	700 / 24.73
Number of light cassettes	2	3
Number of racks (standard/max.)	2 / 9	3 / 12
Weight (empty) (kg/lbs.)	213 / 470	374 / 826
<b>Temperature data</b>		
Temperature range <sup>1)</sup> (°C/°F)	0 – 70 / 32 – 158	0 – 70 / 32 – 158
Temperature range without light cassettes <sup>1)</sup> (°C/°F)	0 – 70 / 32 – 158	0 – 70 / 32 – 158
Temperature range with light cassettes, with illumination <sup>1)</sup> (°C/°F)	10 – 60 / 50 – 140	10 – 60 / 50 – 140
Temperature uniformity at 25 °C / 77 °F (±K)	0.2	0.2
Temperature fluctuation (±K)	0.1	0.1
Max. heat compensation up to 40 °C / 104 °F (W)	400	1000
<b>Climatic data</b>		
Temperature range with light cassettes, with illumination <sup>1)</sup> (°C/°F)	10 – 60 / 50 – 140	20 – 60 / 68 – 140
Temperature uniformity with illumination at 25 °C / 77 °F and 60 % r. H. (±K)	0.6	1.2
Temperature fluctuation with illumination at 25 °C / 77 °F and 60 % r. H. (±K)	0.2	0.2
Humidity range with light cassettes, with illumination <sup>1)</sup> (% r. H.)	10 – 75	10 – 75
Humidity fluctuation at 25 °C / 77 °F and 60 % r. H. with illumination (±% r. H.)	1.5	2
Recovery time after doors were open for 30 sec <sup>2)</sup> at 25 °C / 77 °F and 60 % r. H. with illumination (min)	3	1
<b>Illumination data per light cassette</b>		
ICH compliant illumination for photostability testing <sup>3)</sup> (Lux / UVA W/m <sup>2</sup> )	7500 / 1.4	9900 / 2.1
<b>Electrical data</b>		
Housing protection acc. to EN 60529	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V) / Nominal power at 240 V (kW)	200 – 240 1N~ / 2.4	200 – 240 1N~ / 3.5
Energy consumption <sup>4)</sup> at 40 °C / 104 °F / 75 % r. H. (Wh/h)	600	2350
Model no.	9020-0177	9020-0171

### ► TEMPERATURE-HUMIDITY CHART WITH ILLUMINATION KBF LQC 240

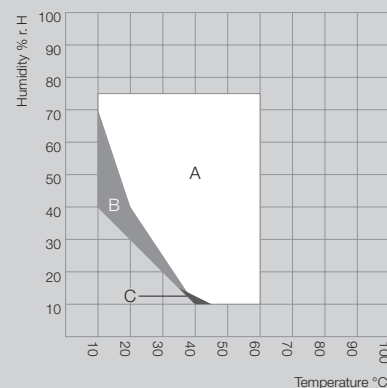

### ► TEMPERATURE-HUMIDITY CHART WITH ILLUMINATION KBF LQC 720

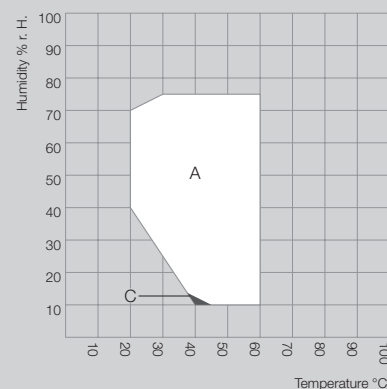

A: Guaranteed condensation-free range  
B: Time-limited operation (max. 24 hours)  
C: Deviations of technical data may be possible


<sup>1)</sup> Lower values are valid up to an ambient temperature of max. 25 °C / 77 °F. // <sup>2)</sup> to 98 % of the set value // <sup>3)</sup> Average value, measured with a spherical sensor (±10 %) by 12 cm / 4.7 in below the light cassette // <sup>4)</sup> Use this value for dimensioning air conditioning systems. /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.



Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)




► DIMENSIONS

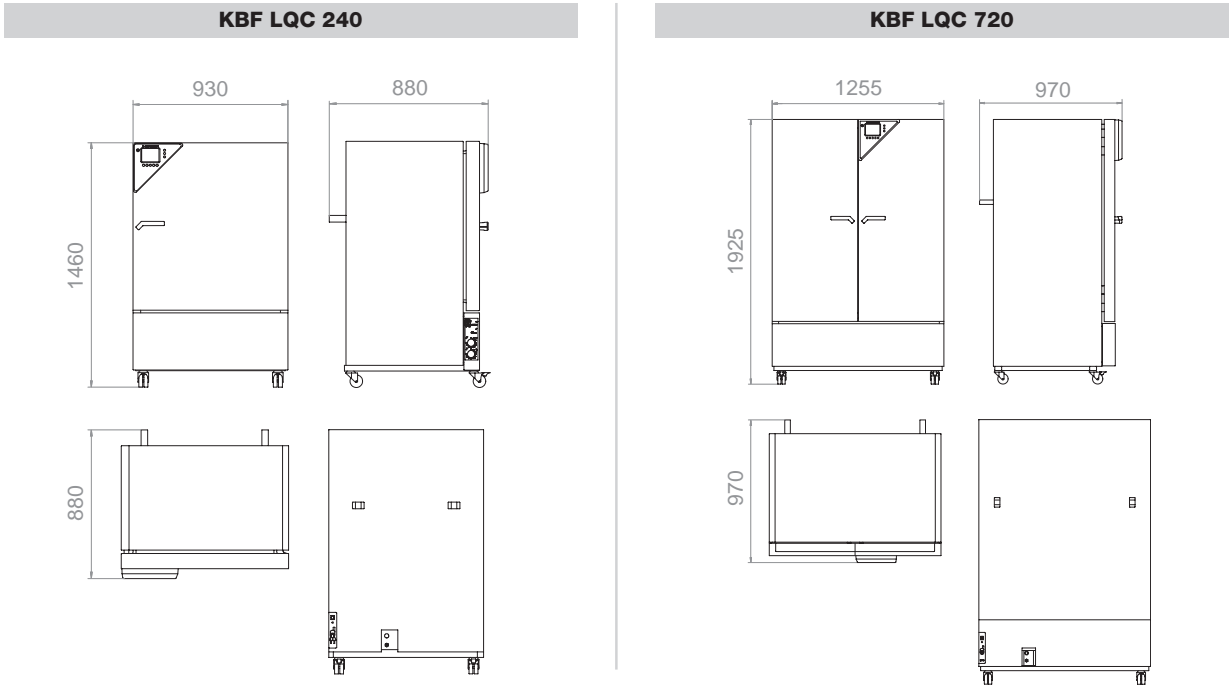

► INSTALLATION REQUIREMENTS

	KBF LQC 240	KBF LQC 720
Wall clearance rear (mm/inch)	100 / 3.94	100 / 3.94
Wall clearance side (mm/inch)	100 / 3.94	200 / 7.87
Nominal voltage (±10 %) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~
Nominal power at 240 V (kW)	2.4	3.5

 **BINDER**

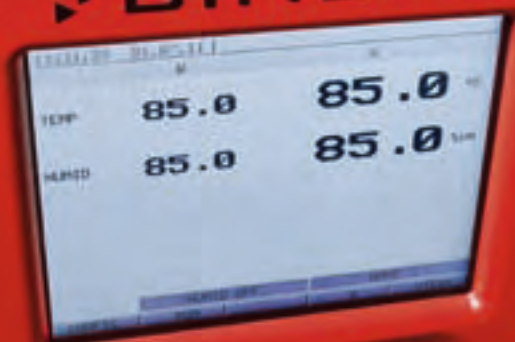

EXIT

ON

ENTER



# Dynamic constant climate testing

## KMF series

### More options: Dynamic constant climate chambers

- ▶ More flexibility
  - ▶ Extended temperature range from -10 °C to 100 °C
  - ▶ Broader climate range up to 90 % r. H.
  - ▶ Defined temperature change rates
  - ▶ Automatic water and wastewater management
- ▶ More reliability
  - ▶ Responsive vapor pressure humidification
  - ▶ Powerful cooling system for safe operation up to 32 °C ambient temperature
  - ▶ Long-term stress testing, e. g. at 85 °C / 85 % r. H.

The dynamic constant climate chambers from BINDER ensure absolutely constant test conditions throughout the testing area. Their greatest advantage is the low space requirement and flexibility regarding the water supply. The wide temperature and humidity ranges make them ideally suited for stress testing series.



Automotive



Plastics industry



Packaging industry

# Constant climate chambers for stress testing

## KMF series

The KMF is the specialist for unconditionally reliable stress testing and precise maintenance of constant test conditions. It has particularly large power reserves and an extremely broad climate range: from -10 °C to 100 °C. It works condensation-free up to 90% r. H. These features make the KMF unique in its class.



► KMF 240 model



Available sizes (liters)

### ► EQUIPMENT

- Temperature range from -10 °C to 100 °C
- Humidity range 10 % to 90 % r. H.
- MCS controller which can store 25 programs of 100 sections each for a maximum of 500 program segments
- User-friendly LCD color screen
- Easy-to-read menu guide
- Integrated electronic chart recorder
- Variety of options for the graphic display of process parameters
- Real-time clock
- Electronically controlled humidification and dehumidification system with capacitive humidity sensor
- Inner glass door
- Independent temperature safety device class 3.1 (DIN 12880) with visual and audible alarm
- Access port with silicone plug Ø 30 mm, left side
- Complete safety connection kit for water supply and drainage, up to 1m in height
- Ethernet interface for communication software APT-COM™ DataControlSystem
- Shelf, stainless steel
- Four casters, two with brakes (KMF 240/720)



## ► KMF SERIES | BEST TEST CONDITIONS:



### Precise climate conditions

- BINDER APT.line™ preheating chamber
  - Constant and gentle circulation of air through large-surface side walls even under a full load
  - Homogeneous climate conditions throughout test specimens



### Fast and precise humidification

- Vapor pressure humidification with fast response times
- Drift-free, capacitive humidity sensor
- Short recovery times after door opening
- Finely adjustable humidity control



### Convenient assembly and operation

- Large access area
- Control elements accessible from the front
- Optimal ratio of usable space and footprint



### Flexible water management

- Sewage pump for discharges up to 1m in height
- Solutions independent of installation sight
- Water treatment with BINDER PURE AQUA SERVICE
- External water supply



### Comprehensive additional services

- Data Logger Kits
- Years of proven validation and documentation materials
- Customer-specific modifications

## ► OPTIONS

- Access ports of various diameters with silicone plug
- Shelf, stainless steel
- Perforated shelf, stainless steel
- Reinforced rack, stainless steel
- Lockable controller keyboard
- Interior lighting
- Additional PT 100 temperature sensor
- RS 422 interface
- External water supply set
- BINDER PURE AQUA SERVICE
- Calibration certificate and extension to calibration certificate
- Measurement of temperature accuracy according to DIN 12880
- Data Logger Kit and software
- Independent temperature safety device class 3.3 (DIN 12880)
- Analog outputs 4 – 20 mA for temperature and humidity measurements
- Door lock



**BINDER PURE AQUA SERVICE**



**Numerous access ports**



**Data Logger Kits**



**Climate chamber with windows, doors and access ports for connecting several measuring instruments**

**BINDER INDIVIDUAL**  
More information on page 133


# Technical data for your planning and installation

## KMF series



	KMF 115	KMF 240	KMF 720
<b>Exterior dimensions</b>			
Width (including access port) (mm/inch)	885 / 34.84	930 / 36.61	1255 / 49.41
Height (including feet/casters) (mm/inch)	1050 / 41.34	1460 / 57.48	1925 / 75.79
Depth (mm/inch)	650 / 25.59	800 / 31.50	890 / 35.04
Depth including door handle, I-triangle, connection, cable (mm/inch)	730 / 28.74	880 / 34.65	970 / 38.19
<b>Interior dimensions</b>			
Width (mm/inch)	600 / 23.62	650 / 25.59	973 / 38.31
Height (mm/inch)	483 / 19.02	785 / 30.91	1250 / 49.21
Depth (mm/inch)	351 / 13.82	485 / 19.09	576 / 22.68
Interior volume (l/cu.ft.)	102 / 3.6	247 / 8.6	700 / 24.6
Number of racks (standard/max.)	1 / 5	1 / 9	1 / 15
Load per rack (kg/lbs.)	30 / 66	30 / 66	45 / 99
Permitted total load (kg/lbs.)	100 / 221	100 / 221	150 / 331
Weight (empty) (kg/lbs.)	127 / 280	185 / 408	309 / 682
<b>Temperature data</b>			
Temperature range <sup>1)</sup> (°C/°F)	-10 – 100 / 14 – 212	-10 – 100 / 14 – 212	-10 – 100 / 14 – 212
Average heating up time acc. to IEC 60068-3-5 (K/min.)	1.3	1.1	1.0
Average cooling down time acc. to IEC 60068-3-5 (K/min.)	0.5	0.6	0.4
Max. heat compensation up to 25 °C / 77 °F (W)	150	350	400
<b>Climatic data</b>			
Temperature range <sup>1)</sup> (°C/°F)	+10 – +90 / 50 – 194	+10 – +90 / 50 – 194	+10 – +90 / 50 – 194
Temperature uniformity <sup>3)</sup> (±K)	0.3 – 1.0	0.3 – 1.5	0.2 – 1.0
Temperature fluctuation <sup>3)</sup> (±K)	0.1 – 0.2	0.1 – 0.5	0.1 – 0.5
Humidity range (% r. H.)	10 – 90	10 – 90	10 – 90
Humidity fluctuation <sup>3)</sup> (± % r. H.)	≤ 2.5	≤ 2	≤ 2
<b>Electrical data</b>			
Housing protection acc. to EN 60529	IP 20	IP 20	IP 20
Nominal voltage (±10 %) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~	200 – 240 1N~
Nominal power at 240 V (kW)	2.0	2.1	3.1
Energy consumption <sup>4)</sup> at 85 °C / 185 °F / 85 % r. H. (Wh/h)	570	500	1050
<b>Model no.</b>			
	9020-0187	9020-0145	9020-0185

### ► TEMPERATURE-HUMIDITY CHART

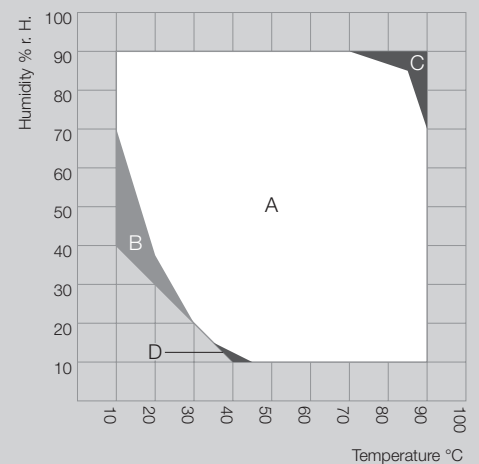




- A: Guaranteed condensation-free range
- B: Time-limited operation (max. 24 hours)
- C: Condensation in the inner chamber may be possible
- D: Deviations of technical data may be possible

<sup>1)</sup> Lower values are valid up to an ambient temperature of max. 25 °C / 77 °F // <sup>2)</sup> to 98 % of the set value // <sup>3)</sup> depending on the set point // <sup>4)</sup> Use this value for dimensioning air conditioning systems. /// All technical data are specified for units with standard equipment at an ambient temperature of 25 °C and a voltage fluctuation of ±10 %. The temperature data are determined in accordance with factory standard following DIN 12880 respecting the recommended wall clearances of 10 % of the height, width and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.






Current information and values are available at:  
[www.binder-world.com](http://www.binder-world.com)

## ► HEATING UP AND COOLING DOWN RATE

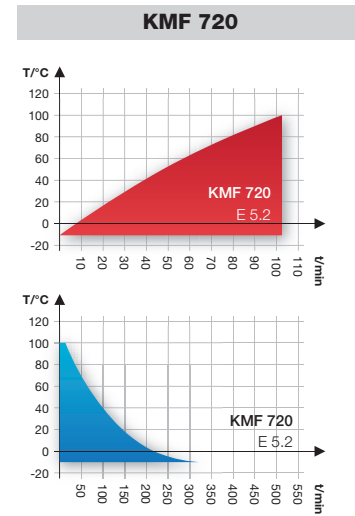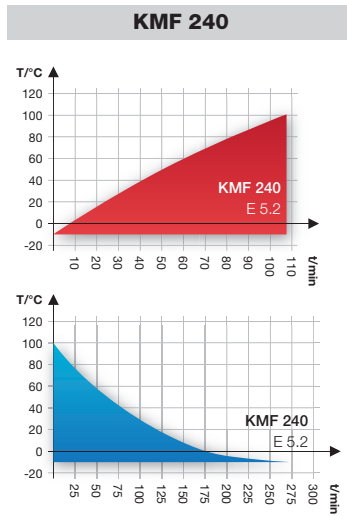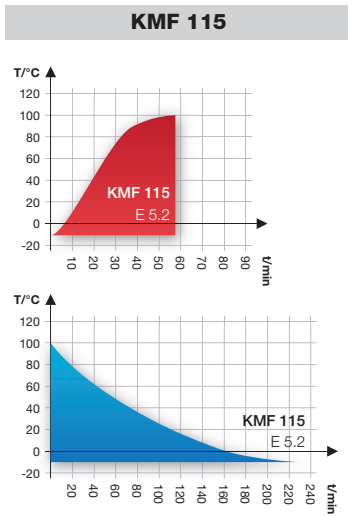

## ► DIMENSIONS

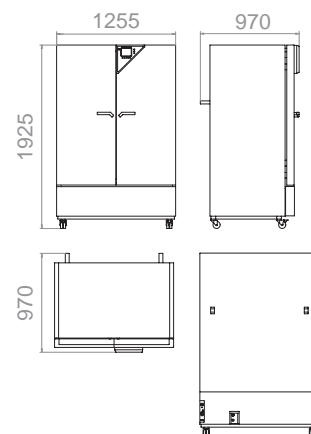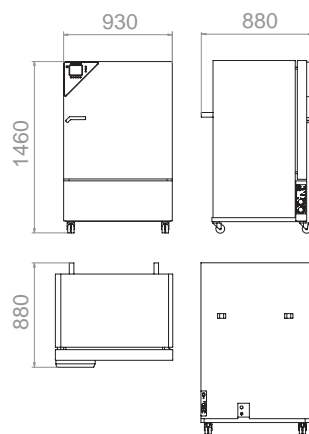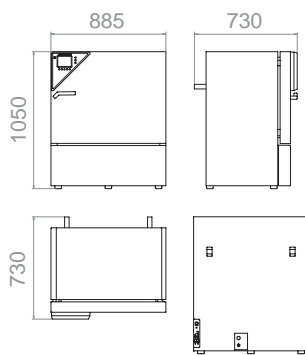

## ► INSTALLATION REQUIREMENTS

	KMF 115	KMF 240	KMF 720
Wall clearance rear (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94
Wall clearance side (mm/inch)	100 / 3.94	100 / 3.94	100 / 3.94
Nominal voltage ( $\pm 10\%$ ) 50/60 Hz (V)	200 – 240 1N~	200 – 240 1N~	200 – 240 1N~
Nominal power at 240 V (kW)	2	2.1	3.1



# Process documentation APT-COM™

## | APT-COM™ DataControlSystem

Working in compliance with standards always requires excellent documentation. APT-COM™ resulted in standard software with a uniquely universal capability: Users can easily achieve process safety and data security, as well as validation of the entire system using standard features. Easy. Cutting edge.

The system, which consists of software in three different editions and the connected equipment, offers features that are needed for tasks ranging from the simplest measurements to guideline-compliant work: seamless monitoring of processes and documentation of process data.

Documentation is automatically generated in electronic format and as a hard copy. This produces guideline-compliant documentation without extra effort, just like having a tailor-made suit for every PC user. Easy to use for a broad range of applications.



# APT-COM™ DataControlSystem

## ► BINDER control and documentation system

### ► Performance potential in 3 classes

The motto here is not “as much as possible”, but rather “as much as necessary”. This has less to do with the ever-present pressure to reduce costs, and much more to do with the fact that processes today have to be as efficient as possible to achieve the best results. This includes a software system that can meet the individual requirements of a multitude of different tests and users, all while maintaining optimal adaptability. This is the reason why we developed three different comprehensive versions of the APT-COM™ software:

#### BASIC


permits remote adjustment of test parameters for the connected equipment, graphic interface programming, and manual documentation of your data.

#### STANDARD

links up several units within a network and provides automatic documentation if required.

#### GLP-Edition


For the highest demands we provide maximum functionality. Most regulatory requirements are met in this area in no time at all. Another strength of this version is remote alerts for limit violations or communication problems by means of independent monitoring and alert functions.

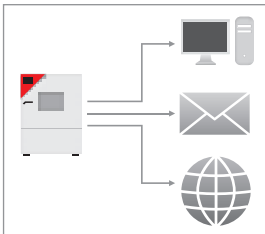

### ► Clear presentation of process data

**Always up-to-date:** Process data can be constantly accessed everywhere: either locally, by email, or through the Internet.


**Important information always available at a glance:** Making a decision on whether or not a process is running well is easy, thanks to the control console function.

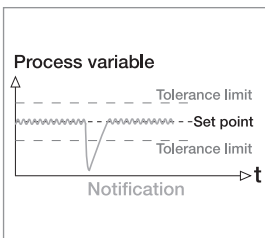

### ► Unsurpassed process safety and security

**Tolerance limits for each monitored parameter:** No parameter can exceed the specified tolerance limits without setting off an automatic alert. The alert is sent via an intranet, the Internet, as email, or as a phone call to the person responsible.

**Access restriction** User IDs and passwords control access to sensitive processes. Different levels of authorization for system changes ensure proper system administration.

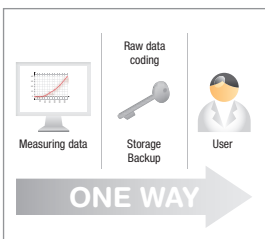


### ► Guideline-compliant data security

**Storage of measured data:** Protected against manipulation by an encrypted format, access restricted to the author or the administrator.

**Backup of measured data:** Automatic backup at user-defined time intervals for storage on all available storage media. Backed up information is easy to find thanks to automatic naming of the backup with a time-stamp.

**Documentation of operator interventions:** Complete documentation of any operator interventions, with user ID, time-stamp, archived protected against manipulation, and automatic backup of measured data.



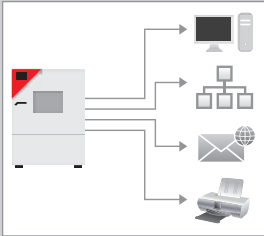


#### ► Time-saving documentation and presentation of results

**Generation of measured data on the monitor:** Measured data is constantly regenerated from protected raw data and protected against manipulation.

**Display of measured data on the internet:** Users can access the process sequence with a standard browser on any PC connected to an intranet or the Internet, even without the APT-COM™ software.

**Coordinated file archiving:** Enables quick and easy display and printing of any past test runs.

**Clear printouts:** The measured data can be printed out automatically at adjustable time intervals. Form fields for comments with respect to the measurements and for authentication ensure correct data assignment and coordination. Signature fields and page numbering provide an easy way to meet documentation requirements with minimum effort.

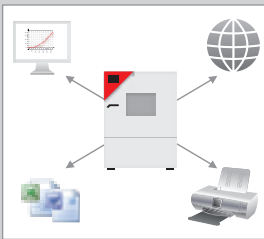

#### ► Control and programming

**Remote monitoring of setpoints over great distances:** Equipment isn't always within close proximity of the workstation. This is why we have provided the option of transmitting process variables to the equipment via PC and reviewing equipment settings.


**Graphic program editor:** Facilitates the easy generation of extensive programs, which can be reviewed and transferred to different units. This saves time and increases transparency.



#### ► System qualification

**Qualification folders with records for IQ and OQ:** Customized for the actual equipment and software configuration. Facilitates system validation, enjoys an excellent reputation among auditors, and saves time when implementing systems. Together with equipment qualification, it is a complete solution that takes full advantage of our experience.

**System qualification:** Our highly experienced BINDER service technicians supervise the proper commissioning and functionality of the system on site, and document these inspections in IQ/OQ protocols. The documentation of these important qualification steps is a comprehensive, time-saving service to ensure successful qualification.

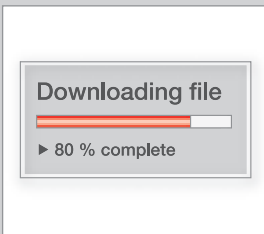

#### ► Always state-of-the-art with updates

**Software updates without added costs:** It goes without saying that our software is constantly being improved and updated. In addition to adding new equipment models, we also incorporate new guidelines and customer suggestions for improvements. Updates are available for free download from the BINDER website. Qualification documents are available for every version of the software.

1

2

3

4

5

6





# Services

## | Always the right support

The professional customer in the scientific lab expects more from a chamber than just the technical features it has when it leaves the factory. In fact, the customer expects a comprehensive solution offering of competent contacts, competent consultation and above all, customer solutions that allow him to perform his daily responsibilities efficiently, effectively and successfully.

BINDER is aware of these customer requirements and, with its extensive service offering, provides the right answer to the individual requirements and expectations of the customer. This extensive market service clearly distinguishes BINDER from the competition and guarantees the additional value that the customer needs.

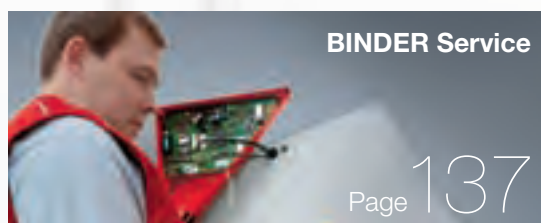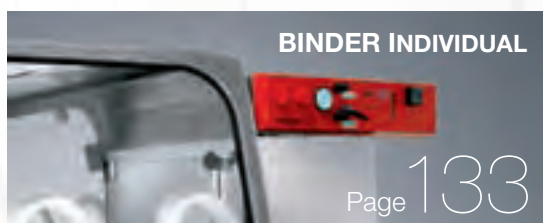





## Customized solutions for customized requirements

Special applications sometimes require a special solution. Precisely for this reason, BINDER has created the BINDER INDIVIDUAL division. The already implemented solutions in this division are very diverse: stainless steel housings and individual access ports are just the tip of the iceberg.

In many cases, special projects require more technical know-how: for example, some customers have particularly heavy specimens that can't be stored in a conventional chamber, while others need a unit with a particulate filter. We have already successfully mastered countless tasks for a wide range of requirements. This is the same for what particular features should distinguish your future BINDER chamber: BINDER INDIVIDUAL will implement them.

## CONTACT

Phone: +49 7462 2005 -0  
Fax: +49 7462 2005 -100  
Email: [BINDER-INDIVIDUAL@binder-world.com](mailto:BINDER-INDIVIDUAL@binder-world.com)

1

2

3

4

5

6

# Reasons for BINDER INDIVIDUAL

## ► Know-how

BINDER INDIVIDUAL is a new name for a proven service. As a manufacturer present in many markets, sectors and applications, BINDER offers you a broad knowledge base gained through years of experience. Some examples:

- Customized optimization of heat, refrigeration, humidity, light, air, CO<sub>2</sub>, or O<sub>2</sub> supply
- Customized measurement, management, control and switching
- Customized connections and outlets
- Customized design of parameters
- Customized integration of accessories (e. g. rollers)

## ► Reasons

Some of the reasons to discuss a customized solution with BINDER:

- To take advantage of know-how transfer through BINDER's extensive experience in a wide range of markets
- To create innovative solutions through individual technical consultation
- To implement simple and cost-effective customized solutions
- To optimize company-specific functionality
- To increase ease-of-use through customized designs
- Because an integrated solution is better than an in-house solution produced after the fact

## ► Service

Our technical support team is made up of application specialists, technical consultants, and engineers. This has the distinct advantage of being able to provide comprehensive services:

- Individual consultation
- Professional planning
- Full application support
- Certified calibration and validation (factory calibration certificate)
- State-of-the-art production

## ► Our Guarantee

All of our components complement each other since they come from a single source. All work meets the requirements of ISO 9001 standards. We also guarantee our customized solutions, and provide an operating manual that covers the additional modifications. In addition we guarantee a supply of spare parts for 10 years; all drawings and components for customized solutions have individual identification codes which are archived, in order to ensure that any spare parts for your customized application are provided correctly and promptly.



► Customized solutions



Drying oven with subdivided interior and hangers in guide rails for tubing



Integrated drying oven in a conveyor operation



Vacuum drying oven glove box



CO<sub>2</sub> incubator with side extension for a rotation mechanism



Safety drying oven with an ultraviolet lamp for testing UV resistance



Climatic simulation chamber with additional windows, doors and access ports to accommodate the various instrument connection leads



CO<sub>2</sub> incubator with integrated rotation mechanism



Vacuum drying oven with specific shelves for the placement of large numbers of very flat specimens



Climatic chamber with full-view glass door and manual access openings



Environmental simulation chamber with reinforced inner chamber and solidly mounted perforated shelves for very heavy test specimens

1

2

3

4

5

6



# BINDER Service

## Best service for your success

Not only does a BINDER chamber provide the “Best conditions for your success” even years later, but it also provides the “Best service for your success”. BINDER Service offers competent consultation and complete services that clearly distinguish BINDER from the competition and guarantee the additional value that the customer needs.

1

2

3

4

5

6

# BINDER Service offering



## ► Validation service

Reduce your equipment qualification and validation costs with a BINDER qualification package as part of your equipment order. BINDER's professional validation service guarantees confidence in meeting validation requirements.



## ► Calibration service

By letting BINDER calibrate your equipment, you can be sure that it will meet all of the requirements for maximum process safety. Our support for you: quick, reliable on-site execution, qualified calibration certificate, manufacturer's inspection plate on the unit and maintenance recommendations.



## ► Replacement parts service

Use only BINDER original replacement parts to ensure full and satisfactory equipment performance and complete, unrestricted manufacturer warranty services. We provide a standard 24-hour delivery service to ensure that you receive the optimum solution for your problems as quickly as possible.



## ► Repair and maintenance service

The BINDER manufacturer service is particularly important when it comes to repairs or maintenance. They know the technology, the little changes and updates the best, invest most in training, diagnostic software, documentation and service products for your success.

## ► Service contracts – Prevention is prudent

BINDER provides you with first-class customized services, precisely tailored to meet your specific requirements. These contracts also cover comprehensive consulting services, current updates and exclusive Internet service support. Customers have the invaluable advantage of extended warranty coverage, including the option of lifetime coverage as part of BINDER's manufacturer service.

- Optimal functionality through preventative maintenance
- Security of constant results through calibration, certificates, etc.
- Discount on replacement parts
- Software updates (APT-COM™)
- Response times based on your requirements
- Intensive service support
- Pool contracts
- Warranty extension



► BINDER INTERNATIONAL SALES- AND SERVICE ORGANISATION




► BINDER GmbH  
Im Mittleren Ösch 5  
D-78532 Tuttlingen



► BINDER Inc.  
Bohemia, NY | USA  
[www.binder-world.us](http://www.binder-world.us)



► BINDER Russia & CIS  
Moscow | Russia  
[www.binder-world.ru](http://www.binder-world.ru)

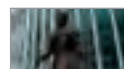

► BINDER Asia Pacific Ltd  
Kwun Tong Kowloon,  
Hong Kong | China



► Shanghai Office  
Shanghai | China

► SERVICE CONTACT

**International Service**





Hotline: +49 7462 2005 555  
Fax: +49 7462 2005 93555  
Email: [service@binder-world.com](mailto:service@binder-world.com)

**USA**

Phone: Tollfree +1 866 816 8191  
Email: [service@binder-world.us](mailto:service@binder-world.us)

# Product selection chart

## ► INNER CHAMBER VOLUME AND TEMPERATURE RANGE

			23 L / 0.7 cuft.	53 L / 1.9 cuft.	115 L / 4.1 cuft.	150 L / 5.4 cuft.	210 L / 7.5 cuft.	240 L / 8.5 cuft.	300 L / 10.6 cuft.	400 L / 14.3 cuft.	500 L / 17.7 cuft.	700 L / 24.7 cuft.	720 L / 25.4 cuft.	
Application	Product	Series												Temperature range
 Growth Page 11	CO <sub>2</sub> incubator	C				●								ambient 7 °C – 50 °C / 13 °F – 122 °F
	CO <sub>2</sub> incubator	CB		●		●	●							ambient 7 °C – 60 °C / 13 °F – 140 °F
	Incubator	BD	●	●	●			●		●			●	ambient 5 °C – 100 °C / 9 °F – 212 °F
	Incubator	BF		●	●			●		●			●	ambient 5 °C – 100 °C / 9 °F – 212 °F
	Refrigerated incubator	KB	●	●	●			●		●			●	-5 °C – 100 °C / 23 °F – 212 °F
	Growth chambers	KBW						●		●			●	0 °C – 70 °C / 32 °F – 158 °F
	Growth chambers	KBWF								●			●	0 °C – 70 °C / 32 °F – 158 °F
 Storage Page 47	Ultra low temperature freezer	UF V							●		●	●		-40 °C – -86 °C / -40 °F – -123 °F
 Drying and tempering Page 55	Drying oven with gravity convection	ED	●	●	●			●		●			●	ambient 5 °C – 300 °C / 9 °F – 572 °F
	Drying oven w. mechanical convection	FD	●	●	●			●						ambient 5 °C – 300 °C / 9 °F – 572 °F
	Heating chamber w. mechanical convection	FED		●	●			●		●			●	ambient 5 °C – 300 °C / 9 °F – 572 °F
	Temperature test chamber	FP		●	●			●		●			●	ambient 5 °C – 300 °C / 9 °F – 572 °F
	Temperature test chamber	M		●	●			●		●			●	ambient 5 °C – 300 °C / 9 °F – 572 °F
	Safety drying oven	FDL			●									ambient 5 °C – 300 °C / 9 °F – 572 °F
	Safety drying oven	MDL			●									ambient 5 °C – 350 °C / 9 °F – 662 °F
	Vacuum drying oven	VD	●	●	●									ambient 15 °C – 200 °C / 27 °F – 392 °F
	Safety vacuum drying oven	VDL	●	●	●									ambient 15 °C – 200 °C / 27 °F – 392 °F
 Environmental simulation Page 103	Constant climate chamber	KBF			●			●					●	0 °C – 70 °C / 32 °F – 158 °F
	Constant climate chamber with light	KBF P						●					●	0 °C – 70 °C / 32 °F – 158 °F
	Constant climate chamber with LQC	KBF LQC						●					●	0 °C – 70 °C / 32 °F – 158 °F
	Dynamic constant climate chamber	KMF			●			●					●	-10 °C – 100 °C / 14 °F – 212 °F



► TEMPERATURE-TIME FUNCTIONS

Product	Series	LCD screen control with extended program functions	Constant temperature	Ramp function	Delayed OFF	Delayed ON	Temperature dependent delayed OFF	Program mode	Week program functions	Fan	Interface	Temperature safety device
CO <sub>2</sub> incubator	C	●	●									3.1
CO <sub>2</sub> incubator	CB		●	●	●	●	●	●	●		RS422	3.1
Incubator	BD		●	●	●						RS422	3.1
Incubator	BF		●	●	●	●	●			🌀	RS422	3.1
Refrigerated incubator	KB		●	●	●	●	●	●	●	🌀	RS422	3.1
Growth chambers	KBW		●	●	●	●	●	●	●	🌀	RS422	3.1
Growth chambers	KBWF	●	●	●	●	●	●	●	●	🌀	RS422	3.1

Ultra low temperature freezer	UF V		●								RS422	
-------------------------------	------	--	---	--	--	--	--	--	--	--	-------	--

Drying oven with gravity convection	ED		●	●	●						RS422*	2
Drying oven w.mechanical convection	FD		●	●	●					●		2
Heating chamber w. mechanical convection	FED		●	●	●	●	●			🌀	RS422	2
Temperature test chamber	FP		●	●	●	●	●	●	●	🌀	RS422	2
Temperature test chamber	M	●	●	●	●	●	●	●	●	🌀	RS422	2
Safety drying oven	FDL		●	●	●	●	●	●	●	●	RS422	2
Safety drying oven	MDL	●	●	●	●	●	●	●	●	●	RS422	2
Vacuum drying oven	VD		●	●	●	●	●	●	●		RS422	2
Safety vacuum drying oven	VDL		●	●	●	●	●	●	●		RS422	2

Constant climate chamber	KBF	●	●	●	●	●	●	●	●	🌀	Ethernet	3.1
Constant climate chamber with light	KBF P	●	●	●	●	●	●	●	●	🌀	Ethernet	3.1
Constant climate chamber with LQC	KBF LQC	●	●	●	●	●	●	●	●	🌀	Ethernet	3.1
Dynamic constant climate chamber	KMF	●	●	●	●	●	●	●	●	🌀	Ethernet	3.1

● available / 🌀 speed control fan / ★ option

## BINDER Product Range

### ► Scientific Products

Drying ovens	ED   FD
Heating chambers	FED
Vacuum drying ovens	VD   VDL
Incubators	BD   BF   KB
CO <sub>2</sub> -incubators	C   CB
Ultra low temperature freezers	UF V
Constant climate chambers	KBF   KBF P/LQC   KMF
Growth chambers	KBW   KBWF
Temperature test chambers	FP   M
Safety drying ovens	FDL   MDL

### ► Industrial Products

Heating chambers	FED
Temperature test chambers	FP   M
Environmental test chambers	MK
Climatic test chamber	MKF
Low temperature test chambers	MKT   MKFT
Constant climate chambers	KMF
Vacuum drying ovens	VD   VDL
Safety drying ovens	FDL   MDL

## Polyscientific Enterprise Sdn Bhd

百利企业有限公司 <sup>152064 D</sup>

**Melaka (HQ)**  
272, Taman Asean,  
Jalan Malim,  
75250 Melaka

T +606 3350690 F +606 3351631

**Shah Alam**  
14, Jalan Serendah 26/39,  
Kawasan Perindustrian HICOM,  
40400 Shah Alam, Selangor

T +603 51036920 F +603 51036980

**Penang**  
16, Jalan Sungai Tiram 7,  
11900 Bayan Lepas,  
Penang

T +604 6371500 F +604 6371600

**Johor**  
No. 365 Jalan Ekoperniagaan 11,  
Taman Ekoperniagaan 2,  
Senai Airport City, 81400 Senai,  
Johor

T +607 5955243 F +607 5955745

[info@polyscientific.com.my](mailto:info@polyscientific.com.my)

[www.poly.my](http://www.poly.my)